



# QUANTAX EDS

Energy-dispersive X-ray spectrometer  
for electron microscopy

- **User Manual**

Bruker Nano GmbH  
Am Studio 2D  
12489 Berlin

Tel. +49 (30) 670 990-0  
Fax +49 (30) 670 990-30  
E-mail [support.bna.emea@bruker.com](mailto:support.bna.emea@bruker.com)  
Internet [www.bruker.com](http://www.bruker.com)

The reproduction, transmission or use of this document or its contents is not permitted without express written authority. Offenders will be liable for damages. All rights reserved.

We have checked the contents of this manual for agreement with the hardware and software described. Since deviations cannot be precluded entirely, we cannot guarantee full agreement. However, the data in this manual are reviewed regularly and any necessary corrections are included in subsequent editions. Suggestions for improvement are welcome.

Order no. DOC-M82-EXS066  
© 2015 Bruker Nano GmbH, Berlin, Germany.

All trademarks and registered trademarks are the sole property of their respective owners.  
Printed in the Federal Republic of Germany.

# Contents

<b>1 Introduction .....</b>	<b>5</b>
<b>2 Safety Information .....</b>	<b>7</b>
2.1 Radiation Safety .....	9
2.2 Beryllium .....	10
2.3 Electrical Safety.....	10
2.4 Electromagnetic Compatibility .....	11
<b>3 The QUANTAX EDS System .....</b>	<b>13</b>
3.1 Hardware System Components.....	15
3.1.1 XFlash® Energy Dispersive X-ray Detectors .....	15
3.1.2 Signal Processing Unit.....	15
3.1.3 Spectrometry Data Interface and EM Control.....	16
3.1.4 Computer Workstations .....	17
3.2 ESPRIT 2 Software.....	18
3.3 General Operating Instructions .....	19
3.3.1 Detector Maintenance.....	19
3.3.2 Electronic Units .....	20
3.3.3 Software and Data Handling.....	21
<b>4 The ESPRIT 2 Software.....</b>	<b>23</b>
4.1 Start-up and Login .....	25
4.2 User Administration .....	26
4.3 Workspaces.....	28
4.4 Display Control .....	29
4.5 Configurator Bar .....	31
4.5.1 Sample Configuration .....	32
4.5.2 Standards Configuration .....	32
4.5.3 Microscope Configuration .....	33
4.5.4 Scan Configuration .....	34
4.5.5 EDS Configuration .....	36
4.5.6 Additional Configurators .....	37
4.5.7 Report.....	38
4.5.8 Project .....	38
4.6 Loading and Saving Data and Results .....	39
4.7 System Functions .....	41
4.7.1 System .....	41
4.7.2 Appearance.....	43
4.7.3 Spectrometer .....	44
4.7.4 Imaging.....	45
4.7.5 WDS .....	47
4.7.6 Microscope.....	47
4.7.7 Stage .....	47
4.8 The Help System.....	47
<b>5 Step-by-Step Guides.....</b>	<b>49</b>
5.1 Preparatory Steps.....	51
5.2 Energy-Channel Calibration .....	53

5.3	Image Capture .....	55
5.4	Image Extension .....	57
5.5	Drift Correction .....	59
5.6	Image Processing .....	61
5.7	Spectrum Acquisition .....	63
5.8	Using the Spectrum Chart .....	65
5.9	Sample Coating Correction .....	68
5.10	System Factor Calibration/ Beam Current Setting.....	70
5.11	Standards Library .....	73
5.11.1	Maintain Standards Library for Bulk Samples .....	73
5.11.2	Maintain Standards Library for Electron Transparent Samples – Cliff-Lorimer Factors .....	76
5.11.3	Maintain Standards Library for Electron Transparent Samples – Zeta Factors .....	78
5.12	Quantitative EDS Analysis of Bulk Samples .....	81
5.12.1	Using the Method Editor .....	81
5.12.2	Standardless Quantification .....	83
5.12.3	Standard-based Quantification with a Temporary Reference .....	86
5.12.4	Standard-based Quantification with Standards Library .....	88
5.13	Quantitative EDS of Electron Transparent Samples .....	90
5.13.1	Standardless Cliff-Lorimer Quantification .....	90
5.13.2	Standard-based Cliff-Lorimer Quantification .....	92
5.13.3	Standard-based Zeta Quantification .....	94
5.14	Object Analysis .....	96
5.15	Line Scan .....	98
5.16	Quantitative Line Scan (QLine) .....	100
5.17	Map Acquisition .....	102
5.18	Quantitative Mapping (QMap) .....	104
5.19	Map Processing .....	106
5.20	Phase Analysis .....	109
5.21	Creating a Report .....	112
5.22	Managing Projects .....	114
<b>Index</b> .....		<b>117</b>

# 1 Introduction

*QUANTAX is the ideal all-purpose energy dispersive microanalysis system for industry, research and education. Different system levels and various options are provided for scaling and tuning of QUANTAX to a broad range of analysis tasks and application environments. QUANTAX microanalysis systems are suitable for scanning electron microscopes, transmission electron microscopes, electron beam microprobes, dual beam devices, as well as other X-ray spectrometry applications.*

*All QUANTAX systems provide state-of-the-art qualitative and true standardless quantitative microanalysis for bulk specimens, polished samples, thin layers, particles, and rough surfaces. This is accomplished by optimized automatic or interactive spectrum analysis methods, fundamental parameter approaches (e.g. P/B-ZAF) and the most exact and comprehensive atomic database available. Different standard-based quantification packages are optionally available. The unique capability of combining standard-based and standardless analysis methods (hybrid quantification) completes the range of analytical possibilities.*

*An intuitive user interface, the flexible project management package, as well as various options for quick and comprehensive report generation complement the analysis toolboxes. The software tools are tailored to meet both the needs of the novice as well as the experienced user. All QUANTAX systems' ESPRIT software includes an online help system and supports remote diagnosis and assistance via the Internet.*

---

*This manual provides a general software overview and a practical step-by-step description of the most common measurement procedures that can be performed with the QUANTAX EDS system. The integrated assistants and the online help of the ESPRIT software provide additional support. Details on the individual hardware parts and additional technical data are contained in separate manuals. The individual device manuals also contain further references regarding operational safety. Please follow all safety instructions closely to avoid hazards to personal safety and equipment. In case of necessary maintenance, reinstallation, severe computer breakdown, hardware changes, etc. the Bruker customer support or your local supplier have to be contacted for further assistance and instructions.*

*According to the modular structure of QUANTAX not all parts of this manual may apply in detail to a given installation nor should be taken literally in all cases. For instance, where the general term "electron microscope" (EM) is used, this applies to the whole class of electron-beam or dual-beam systems; SEM includes also STEM, microprobes, or other scanning devices.*

*The ESPRIT software layout may vary according to the modules installed on your system and therefore may differ slightly from what is shown in the illustrations and figures.*



## 2 Safety Information



2.1	Radiation Safety .....	9
2.2	Beryllium .....	10
2.3	Electrical Safety.....	10
2.4	Electromagnetic Compatibility .....	11





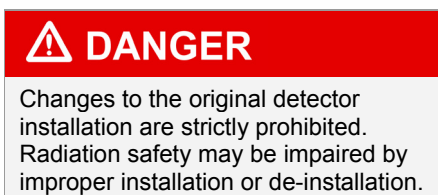


Operation of the QUANTAX EDS microanalysis system is restricted to trained personnel familiar with the system as well as the product documentation, general safety precautions, and laboratory rules. A briefing on safety issues is given at the time of installation or during user training.

Local, state, or federal regulations have to be taken into account additionally to the safety instructions given here. Additional instructions for specific parts of the QUANTAX system - e.g. the X-ray detectors - may be contained in the according manuals.

The QUANTAX EDS microanalysis system may only be used in combination with electron microscopes or similar devices. Any other use beyond that is considered non-intended usage. The operator, not the manufacturer, assumes sole liability for all personal injury and material damage arising from non-intended usage.

## 2.1 Radiation Safety



The QUANTAX microanalysis system uses the X-radiation that is generated by sample interaction with electrons from the electron source of the microscope or photons emitted by the optional X-ray source. Microscopes and attachments approved by the microscope vendors are generally designed to shield this type of radiation sufficiently.



The QUANTAX XFlash® detector is installed according to the applicable radiation safety regulations. Changes to the original installation and equipment, including flanges, vacuum sealing, support, etc. are strictly prohibited.

In case the XFlash® detector has to be uninstalled, make sure that the **original blind flange** of the electron microscope vendor is used to seal the sample chamber port. In case the original flange is not available contact the microscope vendor for support.

## 2.2 Beryllium

### DANGER

Beryllium is highly toxic!

If a beryllium window is destroyed despite all caution, absolutely all fragments must be carefully collected and disposed of according to regulations.



*Beryllium safety issues apply only to X-ray detectors with beryllium X-ray entrance windows. Low energy windows do not contain toxic substances.*

**Beryllium is toxic** if parts are inhaled or swallowed. With X-ray detectors, however, this is only possible in the rare case of a window breaking. Special care must be taken with XFlash® detectors, as they do not contain enough vacuum to suck in the remains of a broken window, as a UHV-Dewar normally does.

In case of an accidental destruction of a beryllium window, gather all fragments thoroughly and collect them using pieces of adhesive tape. The waste has to be labeled and disposed of according to the local safety hazard regulations. Wear a filter mask during cleaning up.

## 2.3 Electrical Safety

### DANGER

High voltage inside. Do not remove covers.



All parts have been designed according to the safety requirements for electrical equipment for measurement, control and laboratory use, or the European Low Voltage Directive, respectively. The system must be correctly installed and used only for the purposes it is designed for. Especially grounding of the system or system parts - as performed during installation - must not be changed for any reason. In case of any supplements or replacement parts being installed, supply voltage settings must be checked and adjusted to the local mains voltage.

Certain parts of the system may contain dangerous voltages. No covers need to be removed during regular operation; maintenance should only be carried out by trained and certified personnel.

## 2.4 Electromagnetic Compatibility



Due to its physical construction, the highly sensitive X-ray detector may not fully comply with common standards regarding electromagnetic immunity of general electronic equipment. This is not a disadvantage under normal circumstances because the QUANTAX system works under the same conditions as electron microscopes and the laboratory environment is designed to meet the requirements of these instruments.

Bruker devices fulfill all requirements regarding active electromagnetic compatibility (emission rules).



# 3 The QUANTAX EDS System



3.1	Hardware System Components.....	15
3.2	ESPRIT 2 Software.....	18
3.3	General Operating Instructions .....	19



## 3.1 Hardware System Components

### Typical QUANTAX EDS system parts

- XFlash® energy dispersive X-ray detector (s)
- Signal processing unit
- Interface and EM control
- QUANTAX server computer
- Optional client workstations
- ESPRIT 2 software package

### 3.1.1 XFlash® Energy Dispersive X-ray Detectors

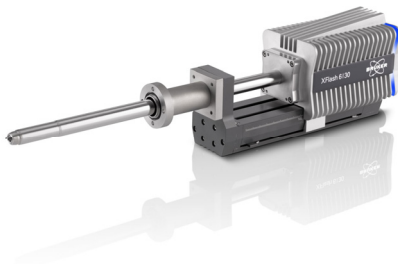


Fig. 3.1-1 Bruker XFlash® detector

### 3.1.2 Signal Processing Unit

System components of QUANTAX comprise the X-ray detector, interface hardware, one or more computers, and several software modules. All parts are selected for performance and trouble free interaction. Exchanging any part of the QUANTAX system with non-approved items (e.g. user defined computers) may impair the functionality.

**Single photon counting.** Energy dispersive X-ray spectrometers take advantage of the photon nature of light. In the X-ray range the energy of a single photon is just sufficient to produce a measurable voltage pulse at the output of an ultra-low noise preamplifier connected to the semiconductor detector crystal. The individual pulse heights are a statistical measure of the corresponding quantum energy. By digitally recording and counting a great number of such pulses within a so-called multi-channel analyzer (MCA), a complete image of the X-ray spectrum is built up almost simultaneously. This digital quantum counting technique makes the energy dispersive spectrometry extremely reliable.

**Bruker XFlash® Detector.** The Bruker XFlash® detector is an energy dispersive X-ray detector that works according to the principle of the silicon drift detector (SDD). It uses a special arrangement of drift electrodes and active components on the detector crystal chip to raise the signal strength to a level that can be measured even at room temperature.

XFlash® detectors are maintenance-free, durable, lightweight, completely vibration free, and do not require any consumables.

The signal processing unit is the heart of the spectrometer electronics. Different auxiliary utilities, like signal inspectors and count rate meters are incorporated in the signal processing unit, as well as all the necessary supply and operational circuits. The data interface connects the spectrometer and the QUANTAX server.



Fig. 3.1-2 Signal processing unit SVE 6

### 3.1.3 Spectrometry Data Interface and EM Control

**Spectrometry signal processing.** Inevitable electronic noise requires the detector signal to be averaged over a certain time for the purpose of pulse height analysis. This averaging or noise filtering is performed by the so-called signal shaper. The necessary time per pulse (filter time) limits the maximum acquisition speed of a spectrometer. The applicable filter times are determined by the properties of the detector noise and the required energy resolution (peak separation).

**Pulse throughput.** With optimum signal shaper the QUANTAX the XFlash® can output 20 000 to 130 000 cps, depending on the type of detector used (refer to the detector's hardware manual for specifications). Faster than optimum shapers can be selected at the expense of energy resolution, which is acceptable for ultrahigh-speed applications like element mapping.

The QUANTAX spectrometer electronics is configured according to the chosen detector and QUANTAX system level.

**Data interface.** The proprietary Bruker MegaLink high-speed data interface transfers the spectrometer data in real time to the QUANTAX server. The PCI plug-in board Bruker IO-Scan can connect up to 4 spectrometers simultaneously. Bruker MegaLink is a serial interface protocol featuring full galvanic isolation for improved noise immunity at connection length of up to 30 m. For basic spectrometer functions that do not require real time spectrometry a standard computer interface (RS232) can be used as an alternative.



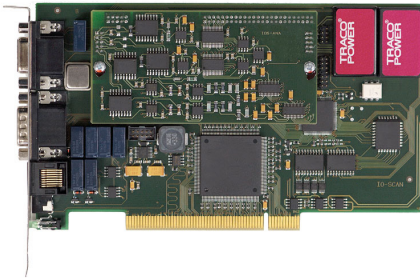


Fig. 3.1-3 PCI-plug in IO-Scan


**QUANTAX scan system.** The PCI plug-in board IO-Scan also contains the necessary hardware for controlling the x/y electron beam deflection as well as two separate digitizers for the image signals of the electron microscope (EM). Position tagged spectrometry data acquisition is provided for advanced line scan and element mapping. Digital signal processing supports the server computer in time critical tasks. Control signals for hardware scan switching are provided as well.

**Switch Box.** The optional switch box allows sharing the external scan interface of the EM between QUANTAX and third party applications. The switching is fully automatic.

**Microscope data transfer.** Microscope data transfer is based on different networking technologies according to the type of microscope. For older brands, serial connections (RS232) are also provided.

### 3.1.4 Computer Workstations

**Client-server architecture.** All QUANTAX systems feature client-server architecture. The server part controls the spectrometer and microscope hardware and performs basic data evaluation; the client system is what is visible to the user and provides the graphical user interface (GUI). Client and server can share one computer or run on different computers. Multiple client workstations can connect to a QUANTAX server with network access (LAN).

 In some cases client and/or server application may share computers with the SEM installation.

**Server workstation.** Based on the operating Windows® system, the QUANTAX server can be used as standalone unit or network component. The connections to the optional client workstations are accomplished either via a separate network or via the intranet of the company or lab. With appropriate upgrades it is also possible to connect to wide area networks. The QUANTAX server performs the measurements and primary spectrum mathematics. It also provides the public and private data areas of the QUANTAX user.

## 3.2 ESPRIT 2 Software

**System requirements.** QUANTAX requires Windows® 7 with a service pack level specified at time of installation. Operating system updates are to be approved by Bruker Nano. Only computers provided or approved by Bruker Nano may be used.


Windows® administrator rights are required for the installation of QUANTAX. Ask your system administrator for help.

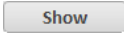
**Hardware driver.** During the first startup of the server PC the driver for the IO-Scan plug-in must be installed. The driver is available from the ESPRIT installation CD, which should always be at hand.

**License file.** A user and system specific license file is contained on a separate CD. During the initial start of ESPRIT the user is prompted to provide this CD for registration.

The license file (Licence.pdf) is stored to the server computer; after completion the license CD should be kept in a safe location.

**Software options.** Licenses control general access to the ESPRIT software as well as installed options. Selected options or the complete license can be time-limited. Updated license files obtained from Bruker Nano can be read in the **System** workspace (see section 4.7). Changes will be enabled after restarting ESPRIT.

 Do not save the license with a pdf-reader. This procedure destroys the license.

The license is also provided as hard copy. The contents can be checked against the file, which can be opened by any pdf-reader or from the **System / System** workspace by clicking the  button.

**Hardware key.** The system license is bound to a system specific hardware key coded into the QUANTAX server hardware. Client workstations can receive certification from a licensed server via the network. Standalone PCs (data stations) must be equipped with an optional USB dongle replacing the hardware coded key.

**USB Dongle.** The QUANTAX dongle for standalone systems (systems without access to a QUANTAX server) fits any USB-port on the regarding PC. A software driver for the dongle must be installed during system installation; the dongle must be present to use the data station.


Install the QUANTAX base system selecting option **Data Station (with dongle)**. Leave the ESPRIT install CD in the drive and plug in the dongle into a convenient USB-port. For details refer to the separate ESPRIT Installation Manual.

Windows® will detect the new hardware automatically and display an installation wizard. Select the options shown in the images below and finish the installation. Note that the dongle does not replace but only enables a valid QUANTAX license.


## 3.3 General Operating Instructions

### 3.3.1 Detector Maintenance

The detector is the most sensitive part of the spectrometer. According to the type of detector at hand, different operations regulations apply. These must be followed strictly.


 After complete shutdown of the XFlash® detector, an extended warmup period and/or energy channel recalibration (see section 5.2) may be necessary for high precision measurements.

**XFlash® detectors.** Bruker XFlash® detectors are maintenance-free. There is no additional wearout caused by permanently powering the detector; nonetheless detectors for light element analysis may exhibit slight parameter changes after very long periods (several weeks) of continuous operation.

 For windowless detectors the standby mode must be enabled before the microscope chamber is vented to avoid contamination on the detector crystal.

**Standby.** Standby mode for the thermoelectrically cooled detector can be chosen between work sessions to prevent unnecessary heating of the detector housing. From the standby mode, the detector can be quickly put back into a stable measuring mode. Standby mode can be entered via the Detector configuration dialog (see section 4.5.5) or by confirming the corresponding option when terminating a work session.

**Cooling mode.** A mode selection for the detector cooling is provided for XFlash® detectors. **Maximum** cooling will provide the optimum energy resolution. **Thermostat mode** will reduce possible changes due to varying environment temperature. Though these changes are normally negligible, thermostat mode is the proposed setting with most QUANTAX applications. Please check the XFlash® Reference manual to decide which of the above options is the most suitable for your purpose.


 A broken detector window caused by any user interaction impairs warranty and can result in a total damage of the X-ray detector. Additionally, a broken beryllium window represents a substantial health risk.

**Detector window.** The X-ray detectors are equipped with a beryllium window (detection range starting from sodium) or a light element window. Both kinds of windows are extremely thin, fragile, and contact sensitive. In general, cleaning of the windows is neither necessary nor permissible. If the detector window is contaminated, contact customer support or your local supplier for assistance.

### 3.3.2 Electronic Units

**Setup.** Unauthorized changes to the electronic and electric installation are not permitted. Besides safety regulations (see section 3.1.2) also electromagnetic immunity and noise pick up issues are addressed during installation.

Note that electronic parts may have to be adapted to the local mains voltages, if this should change for whatsoever reason. Also replacement parts have to be checked for proper settings before being connected to the mains supply. Details are described in the according reference manual.

 The spectrometers are highly sensitive measurement devices, which can be impaired or destroyed by inappropriate use.

**Electrostatic damage.** All electronic components and computers require the typical operating precautions for electronic products. Especially, all usual preventative measures against electrostatic discharge (ESD) must be carried out. All components of the QUANTAX system are designed for laboratory use only.

**Cleaning.** Covers of electronic compartments must never be opened. Only dry cleaning (dusting) of the outer compartments is permissible. The detector window must not be cleaned at all.

### 3.3.3 Software and Data Handling

**Backing up data.** Even state-of-art, complex and versatile software systems cannot be designed to completely exclude the possibility of damage or impairment by improper operation and/or misuse. Additionally no existing computer system is actually immune against potential hardware or software errors. Appropriate backing-up of all relevant data is strongly recommended. No warranty can be accepted in case of data loss for whatsoever reasons.

**Maintaining system integrity.** The QUANTAX system is installed on special selected and configured computers. Any change of the system configuration, user status, access rights, or else can impair system integrity. Also new applications, especially multi-media, internet, and game software can interfere with QUANTAX. The internal client-server architecture of QUANTAX depends on Windows® communication services that may be adversely modified by subsequent installations.

**Recovering from errors.** In cases of program lockup due to improper operation or unexpected circumstances a restart of ESPRIT is normally sufficient. In more severe cases a restart of the personal computer and/or the QUANTAX server may be required. More severe system breakdowns - if any - should be handled by trained service personnel.



## 4 The ESPRIT 2 Software



4.1	Start-up and Login .....	25
4.2	User Administration .....	26
4.3	Workspaces.....	28
4.4	Display Control .....	29
4.5	Configurator Bar .....	31
4.6	Loading and Saving Data and Results .....	39
4.7	System Functions .....	41
4.8	The Help System.....	47





## 4.1 Start-up and Login

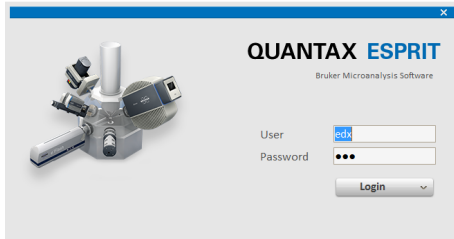






Fig. 4.1-1 ESPRIT Login screen

 System functions, especially concerning the measurement system (e.g. calibrations) will affect subsequent users. It is advised to exercise special caution here.

The **ESPRIT** software runs on the QUANTAX server and the client workstations. However the user only has to start the client application; the server software and communication drivers will start and log in automatically.

**Start-up.** To start ESPRIT click the program icon  on the Windows® desktop or in the Windows® Start menu. The ESPRIT login screen is displayed prompting user name and password.

**Login.** Enter user name and password and click the  button. The password is stored on the local workstation, if the corresponding option is checked. The name of the last user is always retained.

**Multi-user systems.** On multi-user systems several users can access, manage, and evaluate private and shared data simultaneously. The unique user name and password will give access to the private data and settings. Select the server to log in from the pull down menu below after clicking on the arrow symbol of the  button.

After logging in with a unique user name the current settings and data will be loaded from the private user profile volume on the QUANTAX server.

**Remote login.** With client-server installations (option LAN) it is possible to log in from any connected remote workstation. The first user to start a measurement will gain access to the spectrometer and imaging system. This user will be able to perform measurements from the remote workstation. The access to the acquisition hardware remains locked to others as long as this user is logged in.

## 4.2 User Administration



Fig. 4.2-1 QUANTAX communication server access

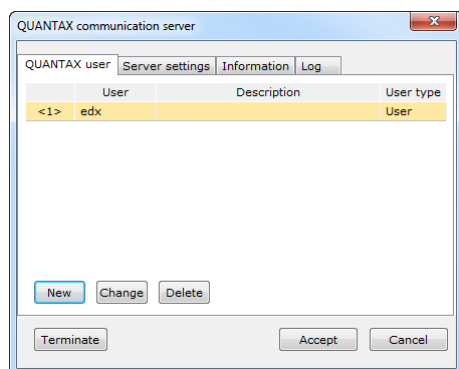


Fig. 4.2-2 QUANTAX user list

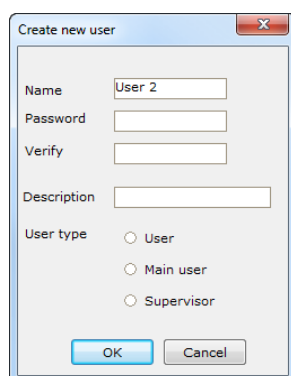




Fig. 4.2-3 Assigning a new user

The **QUANTAX communication server** controls the user assignments as well as the internal and external client-server communication. Service tools and information for troubleshooting are also provided. The communication server runs on the QUANTAX server computer. It is also present in single computer and standalone systems.

Clicking the Bruker icon  in the Windows® system tray of the server computer opens the console of the communication server. With exception of the actions described below no changes to the communication server settings must be made. Unauthorized altering of settings can render the system useless and lead to severe data loss.

**New users** are to be assigned from the QUANTAX communication server. Double clicking the Bruker icon  in the Windows® system tray opens the console of the communication server.


The tab **QUANTAX user** lists the current assigned users. A new user can be added after clicking the **New** button. A password can be entered (leaving the field empty is permitted) and an optional user description can be added. As user type **User**, **Main user**, and **Supervisor** can be selected.

**User** has limited access to the ESPRIT **System** workspace. This user level only allows access to the **Appearance** tab and can review the connected instruments and license.

The **Main user** has limited access to the ESPRIT **System** workspace. This user level has full access to the tabs **System**, **Appearance** and can calibrate the connected EDS detectors as well as the imaging system. The **Main user** has no access to the service settings of the WDS detector and the electron microscope driver.

The **Supervisor** has full access to all tabs in the ESPRIT **System** workspace.

Each new user is automatically assigned a folder structure in a private volume of the server. Default methods, a predefined user profile, and demo data are copied to the user folders.

**Altering user name and password** can be done at any time from the QUANTAX communication server (icon  in the Windows® task bar) by clicking **Change** in **QUANTAX user** tab.


The user data and profile are automatically copied to the new user data structures. Backups of important data, which are always recommended, are especially advisable when changing user names.

Normally there is no need to delete assigned users. If users have to be deleted, it has to be made sure that all user data to be retained are copied to public volumes.

**To delete a user** highlight the user name in the list provided in the **QUANTAX user** tab of the QUANTAX communication server. After clicking **Delete** the deletion of the user data has to be confirmed separately. Caution, once confirmed, the deletion cannot be revoked! If deletion of the data is not confirmed, the data are retained invisibly; a new user with the same name as a previously deleted will be assigned to the idle data.



Fig. 4.2-4 Access to QUANTAX server settings

 The communication server port in the top box must not be altered; it must always match the settings on the client workstations.

The number of assigned measure client ports determines how many users can log in simultaneously to the server computer. Per default only one port is assigned, so that only one user at a time can log in to the server computer. Even when multiple ports are present, the hardware can only be controlled by one user at a time.

Select a new port number in the field **Measure client ports** and click **Add >**. Usually, the number following the last assigned port in natural order is adequate. Ask your system administrator for help.

The **connection setting** for a client workstation is accessed from the ESPRIT login screen. Select the entry **Configuration** from the **Server** dropdown list. The **QUANTAX SERVER SETTINGS** dialog will open.

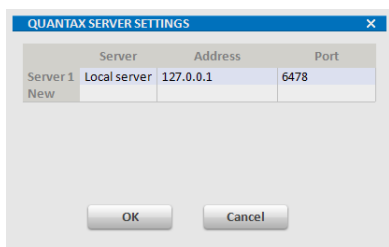


Fig. 4.2-5 Assigned measure client ports

## 4.3 Workspaces

A random name can be given to a connection. If the client workstation also runs a server, select "Local server" for a name and input network address 127.0.0.1. The port number must match the setting of the communication server (see section above); default is 6478. Ask your system administrator for help with network addresses and port numbers.

Workspaces are ESPRIT screen areas in which images and spectra are displayed, the analysis is performed and the results are managed. A workspace can be selected by clicking on the corresponding workspace icon. The active workspace icon is highlighted.

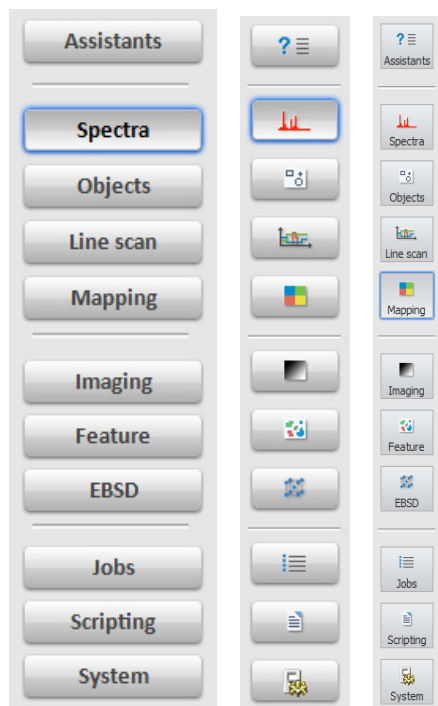


Fig. 4.3-1 ESPRIT workspace list in normal, fullscreen and large style buttons mode



The name of each workspace is displayed on the icon, if the **Large style buttons** checkbox under **System / Appearance** is activated.

### Assistants

gives access to a step-by-step guide for common analysis procedures.

### Spectra

allows spectrum acquisition and analysis of (saved) X-ray spectra.

### Objects

permits point, multipoint and area EDS analysis.

### Line scan

is used to perform qualitative and quantitative line scan EDS analysis.

### Mapping


allows the acquisition of maps: intensity maps and HyperMaps (Bruker's position-tagged spectrometry tool (spectral imaging)). Mapping also includes the chemical phase analysis tool **AutoPhase**.

### Imaging

allows capturing and processing saved electron microscope images.

### Feature

is used to perform particle analysis and chemical classification. Refer to the ESPRIT Feature User manual for details.

 When changing the workspace during an active measurement, a confirmation dialog box pops up and after clicking OK the current measurement will be terminated.

### DANGER

Take special care when changing any parameters in the **System** workspace. These changes affect all QUANTAX user profiles and may influence measurement results. Using the system with wrong **Microscope** and **Stage** parameters may damage sample, EDS detector and electron microscope.

## 4.4 Display Control

### EBSD

gives access to the fully integrated EBSD (electron backscattered diffraction) analysis package. Refer to the QUANTAX EBSD User manual and the EBSD Quick reference guide for details.

### Jobs

is the workspace for configuring ESPRIT's powerful analysis automation tool and running unattended fully automatic analysis tasks.






### Scripting

permits writing and running user-defined scripts.


### System


is used to adjust global settings and calibrate the X-ray spectrometer and image system. Additional functions: display settings (language, font size); license management; communication with electron microscope and motorized sample stage.

**Program window and title bar.** The title bar at the top of the ESPRIT program window contains the general control buttons common to the Windows® applications.

The complete ESPRIT window can be resized by dragging and pulling the handle on the bottom right-hand corner or on one of the border lines. To switch to full screen display, click the  icon in the title bar; switch back by clicking the  icon. The  icon minimizes the program window. The  icon activates the program help system. Please note that the  icon closes ESPRIT completely, not just the active workspace.

To scale individual parts of the program window the splitter (dark grey lines) between different screen areas on the ESPRIT interface can be dragged (a changing mouse cursor indicates availability of this option).

**Full screen / dual screen mode.** An image or a diagram chart can be decoupled from the ESPRIT window and switched to full screen by clicking the full screen symbol  on the upper right-hand corner of the corresponding screen area.

If a second monitor is attached, the decoupled window can be dragged and pulled to the second monitor and ESPRIT can be used in extended display mode. Revert from full screen mode by pressing the ESC key or clicking the full screen symbol  again.


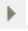
**Configurator bar.** The configurator bar compiles all attached hardware parts as well as information about the sample and standards, the scan settings and the report and project editor. All parts of the configurator bar are described in detail in section 4.5.

**Image capture chart.** The workspaces **Objects**, **Line scan**, **Mapping** and **Imaging** have a dedicated chart for controlling the electron image capture. Analysis spots, objects, scan lines, and mapping areas can be defined on the captured image. Image capture is controlled directly from the current workspace. Additionally, a special window for fast image preview is provided. For details please see section 5.3.

One or two of the image detectors (typically the SE and BSE detectors) can be connected to the QUANTAX hardware and selected for image capture.

**Spectrum and Result chart.** The workspaces **Spectra**, **Objects**, **Line scan**, and **Mapping** have a dedicated chart for displaying spectra and associated spectrum list or results. Details are described in section 5.8.

**Icons and buttons.** A number of icons and buttons associated to the different workspaces and display charts are provided. All icons can be displayed in normal or large style button mode which can be changed by activating/ deactivating the corresponding checkbox in the **System/Appearance** workspace.

Using the triangles  at the control buttons opens drop-down menus to enter acquisition and analysis options. Diagram axis labels and column titles can be selected with .

Data lists and thumbnail images will be highlighted by clicking onto them. Multiple items can be selected by using the SHIFT or CTRL key to perform batch processing.

## 4.5 Configurator Bar

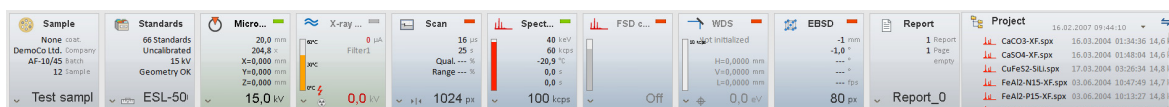


Fig. 4.5-1 Configurator bar (large)





Fig. 4.5-2 Configurator bar (minimized)

The configurator bar consists of a horizontal row of buttons (configurators) in the upper part of the ESPRIT window that allow the configuration of the hardware components available in the QUANTAX system. The number of configurators that actually appear on the screen depends on the license and may differ from the images shown in this manual.

The individual configurators display relevant adjustable parameters of the sample, standard libraries, X-ray excitation sources, scan unit, spectrometers and other detectors, as well as the report and project tools. They will be explained in detail in the next sections.

The configurator bar is resizable by moving the lower splitter (dark grey line). Pulling the splitter upwards minimizes the configurator bar. When minimized, the individual configurators show only one parameter.

The parameter to be displayed can be set by the user by clicking with the right mouse key on the displayed value in the configurator and selecting the desired parameter in the **Select entry** menu.

Use the downward-facing arrow in the left bottom corner of a configurator ( or  icon, depending on the display mode) to access the corresponding configuration dialog.

### 4.5.1 Sample Configuration

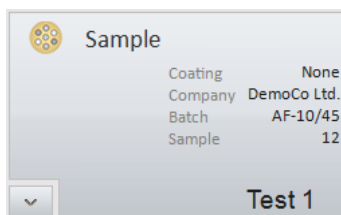


Fig. 4.5-3 Sample configurator

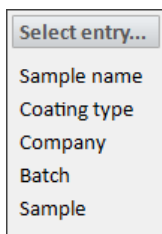


Fig. 4.5-4 Sample configurator select entry menu

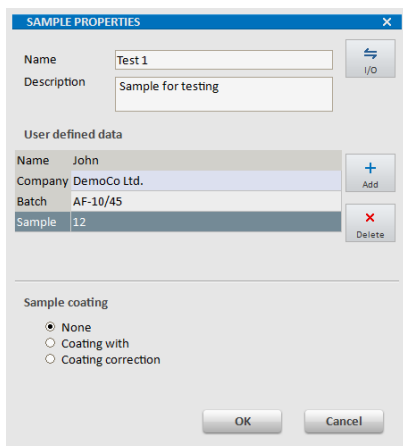

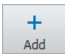


Fig. 4.5-5 Sample properties dialog

The **Sample configurator** is used to display and manage information about the sample. The information displayed is divided into the following parameters:

- Sample type
- Coating type
- Company
- Batch
- Sample

The parameter to be displayed can be set by the user by clicking with the right mouse key on the displayed value in the configurator and selecting the desired parameter in the **Select entry** menu.

When clicking on the  icon in the bottom left corner of the **Sample** configurator, the **SAMPLE PROPERTIES** dialog pops up. Before measurement, general properties of the sample can be entered in this dialog. All subsequently collected data (spectrum, image, line scan, map) will contain these parameters. Default input parameters are name and description. Additional user-defined data can be added clicking the  icon.

**Sample coating.** If the sample is coated to enhance electrical conductivity, this can be entered under **Sample coating**.

- **None.** Select this option, if the sample is not coated.
- **Coating with.** Select this option, if the sample is coated. Select the element of the coating material in the periodic table which pops up after selecting this option. The selected element will be considered for the peak deconvolution algorithm but will not be quantified. The display changes to **Coating with "Element name"**.
- **Coating calibration.** Select this option, if the sample is coated and a coating correction data file (.ccc) is available. Open the \*.ccc file in the popup dialog.

### 4.5.2 Standards Configuration

The **Standards configurator** is available, if a standard-based EDS quantification license is



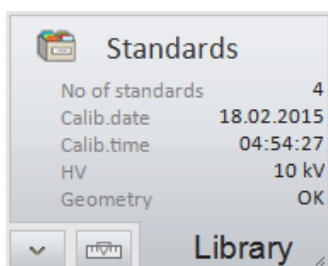




Fig. 4.5-6 Standards configurator

activated. The configurator displays the following information about the system and the standards library:

- Number of standards in the standards library
- Calibration date (of the system factor)
- Calibration time (of the system factor)
- HV
- Geometry

**Standards library.** When clicking on the  icon in the bottom left corner of the Standards configurator, the **ADMINISTRATION OF STANDARD LIBRARY** dialog pops up. For further details see section 5.12.4.

**System calibration.** When clicking on the  icon in the bottom left of the Standards configurator, the **SYSTEM FACTOR CALIBRATION** dialog pops up. For further details refer to section 5.10.

### 4.5.3 Microscope Configuration

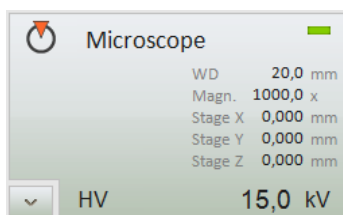


Fig. 4.5-7 Microscope configurator

The **Microscope configurator** displays the main parameters of the electron microscope.

- Magnification: Image magnification set on the electron microscope.
- High voltage: Acceleration voltage set on the electron microscope.
- WD (working distance): Sample – pole piece distance [mm].
- Stage X, Y, Z: Microscope stage position coordinates are displayed, if available

The parameter to be displayed can be set by the user by clicking with the right mouse key on the displayed value in the configurator and selecting the desired parameter in the **Select entry** menu.

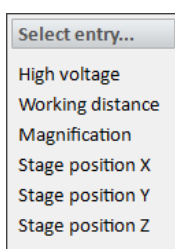



Fig. 4.5-8 Microscope configurator select entry menu

When clicking on the  icon in the bottom left corner of the **Microscope** configurator, the **MICROSCOPE CONFIGURATION** dialog pops up. If the data communication between microscope and ESPRIT is active, the microscope parameters **Magnification**, **High voltage** and **Working distance** will be transferred from the microscope to ESPRIT and displayed in this dialog.

**Communication active.** This option should be

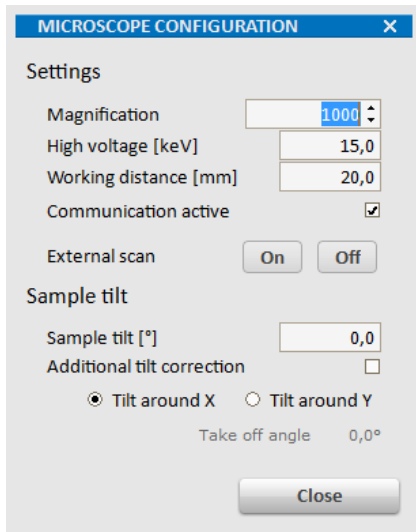


Fig. 4.5-9 Microscope configuration dialog

Because sample tilt affects the take-off angle and hence the Bremsstrahlung background, its value affects the quantification results.

#### 4.5.4 Scan Configuration

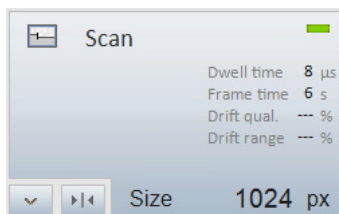


Fig. 4.5-10 Scan configurator

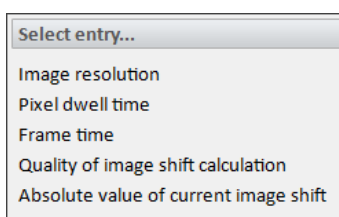


Fig. 4.5-11 Scan configurator select entry menu

checked to maintain communication between the microscope and the ESPRIT software, otherwise relevant microscope parameters for image scaling and quantitative analysis will not be transferred to ESPRIT. The communication is set up during installation and automatically enabled after starting ESPRIT.

If the communication is disturbed or not available with the current type of microscope, the microscope data has to be entered manually. To disable automatic data transfer intentionally, uncheck the control box **Communication active**.


**Sample tilt [°]**. If the sample is tilted, the value of the sample tilt (in degrees) has to be entered in this field manually.

**Additional tilt correction.** Check this option, if the sample is tilted and the microscope has no image tilt correction capability. This is recommended when performing EBSD analysis using large tilt angles to correct image distortion of microscope images, element or orientation maps.

The **Scan configurator** displays the scanning parameters:

- Image resolution: Image size in pixels
- Dwell time: Signal collection time of a single pixel during image acquisition
- Frame time
- Drift qual.: Quality of image shift calculation
- Drift range: Absolute value of current image shift

The parameter to be displayed can be set by the user by clicking with the right mouse key on the displayed value in the configurator and selecting the desired parameter in the **Select entry** menu.

When clicking on the  icon in the bottom left corner of the **Scan** configurator, the **SCAN CONFIGURATION** dialog pops up.

**Image resolution.** The entered value applies to the x-axis of the image. Since the aspect ratio of the image is defined by the microscope and set during installation, the value for the y-axis is set automatically and is not displayed.

**Image inputs.** If more than one image detector is connected to the QUANTAX scan system, the

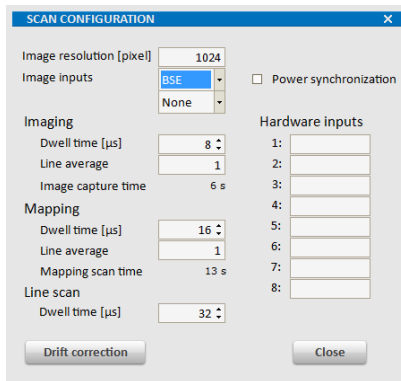



Fig. 4.5-12 Scan configuration dialog

The dwell time can only be set in increments of numbers which are a power of two. Direct input of arbitrary numbers is not possible. Use the arrows in the input field.

Very high image resolution will slow down operation.

For **Imaging**, **Mapping** and **Line scan** different timing settings are maintained. However, the electron image captured in the left image window of each workspace always follows the settings for **Imaging**.

The scan speed can also be set from the dialog that opens after clicking the  icon next to the image chart.

desired detector can be selected here. For each channel only one image detector can be selected at a time. If only Ch1 and Ch2 can be selected from the dropdown menu, note that the image channel names can be changed in the System workspace under Microscope to e.g. SE, BSE, HAADF,... (refer to section 4.7.6).

**Imaging.** The **Dwell time** is the time the electron beam stays on a pixel while capturing an image. Select expanded dwell times to allow filtering and averaging of the image signal and obtain less noisy images. The **Line average** factor controls the number of times a horizontal line is scanned in the slow scan direction and averaged before proceeding to the next line. Increasing this factor provides noise reduction and expanded image scan time while limiting the increased stress on the sample by maintaining a high scan speed. The line average factor will multiply the total image acquisition time. The total **Image capture time** is calculated and displayed.

**Mapping.** The **Dwell time** is the time the electron beam stays on a pixel while acquiring an EDS map. The **Line average** factor for mapping controls the number of times a horizontal line is scanned in the slow scan direction and averaged before proceeding to the next line. The capture time (**Mapping scan time**) for one mapping frame is calculated by multiplying the numbers of pixels and lines with the dwell time and the line average factor. The real scan time may be longer due to the time needed for data processing by the computer.

**Drift correction.** When clicking on the **Drift correction** button, the **IMAGE DRIFT CORRECTION SETTINGS** dialog pops up. It is described in detail in section 5.5.

**Power synchronization.** When the according control box is checked, the scan process is synchronized to the cycles of the AC mains voltage. Power synchronization substantially reduces blurring of vertical lines caused by electromagnetic interference (at the expense of the scan speed). It is applicable for capturing any image or map. However, it is normally only useful in connection with very high magnification of the microscope. A special AC/AC adapter to be connected to the QUANTAX server is provided.

**Hardware inputs.** The IO scan card provides 8

### 4.5.5 EDS Configuration

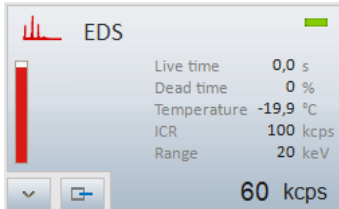


Fig. 4.5-13 EDS configurator

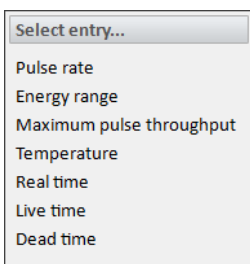


Fig. 4.5-14 EDS configurator select entry menu

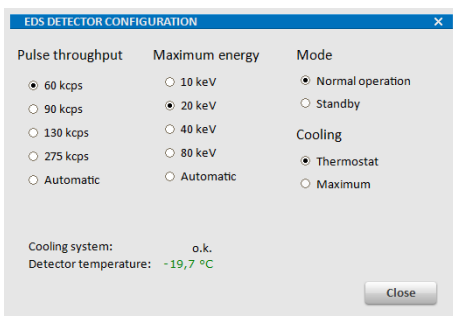


Fig. 4.5-15 EDS detector configuration dialog, pulse throughput choices may vary according to detector type installed

The best energy resolution (peak separation) is available at lower maximum pulse throughput settings.

The maximum energy setting affects the width of an energy channel.

counter inputs for TTL impulse signals. Detailed information is given in the IO scan reference manual.

The **EDS configurator** displays the measuring parameters of the available EDS detector(s).

- Pulse rate (input count rate)
- Energy range
- Maximum pulse throughput
- Temperature
- Real time
- Live time
- Dead time

The parameter to be displayed can be set by the user by clicking with the right mouse key on the displayed value in the configurator and selecting the desired parameter in the **Select entry** menu.

The rectangle in the upper right corner of the **EDS** configurator indicates the status of the instrument. A green rectangle means that the detector is activated and ready to use, a red rectangle indicates that the detector is idle. Toggling between activated and deactivated states can be achieved by clicking on the EDS configurator.

When clicking on the icon in the bottom left corner of the **EDS** configurator, the **EDS DETECTOR CONFIGURATION** dialog pops up. The following parameters can be set:

**Pulse throughput.** This control box lists the signal processor settings available for the spectrometer hardware. When **Automatic** is checked, QUANTAX selects the throughput setting according to the current pulse load at the beginning of a spectrum acquisition or measurement.

**Maximum energy.** The available energy ranges are listed. When **Automatic** is checked, QUANTAX sets the energy range according to the microscope high voltage (kV).

**Standby mode** can be activated from the device control box or when terminating the work session. At start-up, detectors in standby mode will automatically be switched to full operation. Detectors are **not** switched to standby mode by deselecting them in the configurator bar.

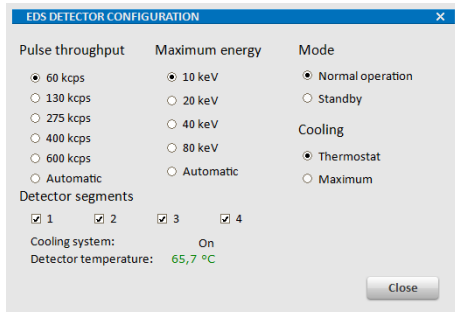



Fig. 4.5-16 Enabling detector segments


### 4.5.6 Additional Configurators


**Cooling mode.** The XFlash® detectors can be operated in **Thermostat** mode, where the Peltier cooling operates at constant temperature (default setting) or in **Maximum** cooling mode, where the Peltier cooling operates at full power.

**Detector segments.** If the detector has multiple segments (or if several detectors were installed as such), only one EDS configurator is shown in the configurator bar. In this case the detectors or detector segments are listed as **Detector segments**. The individual detectors can be selected by enabling the respective checkboxes in the detector segments list.

**Multiple detectors.** Multiple detectors can be registered with individual configurators. In this case, parameters of the spectrometers have to be set individually. When multiple detectors are used, activate the detectors by clicking on the corresponding configurators.

**Motorized detector axis.** If the EDS detector is equipped with a motorized axis, the  icon in the bottom left corner of the EDS configurator (see Fig. 4.5-17) displays and controls the position of the EDS detector. The icon shows one of the two states:

The EDS detector is inserted and is at measurement position, if the  icon is displayed. The system is ready for use. When clicking on the icon in this status, the message **Detector will be moved to park position** pops up and after confirming with OK, the detector will be retracted.

The EDS detector is retracted and in park position, if the  icon is displayed. When clicking on the icon in this status, the message **Detector will be moved to acquisition position** pops up and after clicking on OK, the detector moves to the measurement position.

If additional detectors or X-ray sources are attached, their configurators are displayed in the configurator bar.

**EDS detector(s).** If additional EDS detectors are used, the parameters of each individual detector can be set in the individual configurators. For details see also the XFlash® Reference Manual.

**WD spectrometer.** If a WD spectrometer is attached to the system, it can be configured under the WDS detector configurator. For details see the QUANTAX WDS User Manual and the XSense Reference Manual.

**EBSD detector.** If an EBSD detector is connected, it can be configured under the EBSD detector configurator. For details see the QUANTAX EBSD Quick Reference Guide and the eFlash Reference Manual.

**X-ray source.** If an X-ray tube is connected, it can be configured under the X-ray source configurator. For details see the QUANTAX Micro-XRF User Manual and the XTrace Reference Manual.

### 4.5.7 Report

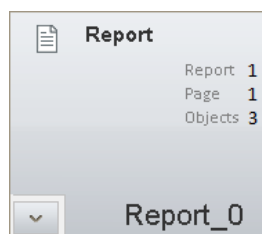



Fig. 4.5-18 Report configurator

The **Report configurator** shows the number of opened reports, the number of pages, objects and the name of the selected report.

Clicking on the  icon in the bottom left corner of the **Report** configurator opens the report editor. Further details on how to use the report function are described in section 5.21.

### 4.5.8 Project

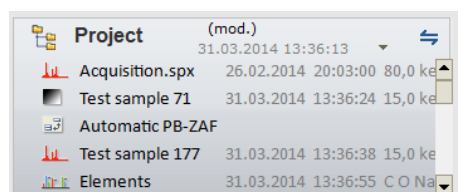
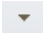
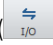


Fig. 4.5-19 Project overview

The **Project** tool serves as data clipboard for data transfer between ESPRIT workspaces or save all data added to **Project** as one file to disk (.rtx file format). Images, spectra, element selection, quantification methods, result tables, map/phase images, line scan data (including point spectra) can be added to and stored in a project.

The **Project** is either docked into the configurator bar on the top of the screen or docked onto the right side of the screen above the **Report/ Report preview**. The two positions of the **Project** (and **Report**) can be toggled by clicking on the  icon next to the project name. See section 5.22 for further details.

## 4.6 Loading and Saving Data and Results

**Loading and saving data or adding to project or report** can be done in two different ways: using the Import/ Export function (  icon) or with a right mouse click into the data field.

A local, workspace-relevant or chart-relevant menu pops up providing the options:

- Load
- Open
- Add to project
- Add to report
- Save
- Copy
- I/O
- Print
- Export
- Properties.

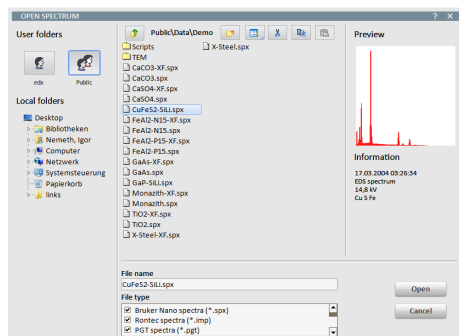




Fig. 4.6-1 File browser

**Loading data.** Data files can be loaded or imported into the current project, into an ESPRIT report or directly into the workspace. To open the file browser for loading data into the workspace, open the according Import/Export menu (  icon) and click **Load** (see also sections 5.21 and 5.22 for information on loading reports and projects, respectively).


Enter a filename or select a file in the list. A preview image and an info box will be displayed, if applicable.

The private data volume on the QUANTAX server is selected via the topmost icon labeled with the user name. To open the shared server folder or access the local or network drives, click the button labeled **Public** or a drive letter in the list below.

Any volume can contain an unrestricted number of folders and subfolders. To go one level up, click the  icon in the headline or use the Backspace button.

A list of file types selects multiple files for browsing. Only file types selected under **File type** are shown in the browser.

Spectra can be loaded to spectrum charts; images can be loaded to the processing area of the **Imaging**

 In client-server installations, local and network drives refer to the client computer. Access rights are defined by the local settings and network administration.

Note that different logical volumes are provided for general data, projects, reports, templates, and methods on the QUANTAX server.

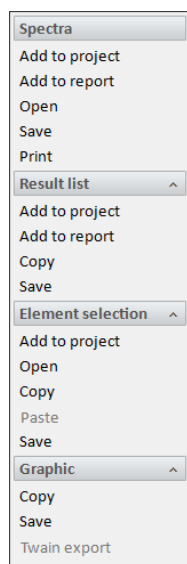

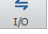


Fig. 4.6-2 Import/Export menu (e.g. Spectrum workspace)

 The file browsers for loading and saving files also allow the control of the folder structure, copying and moving files between private, public, and local volumes, and performing other common file actions.

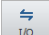
workspace or to the **Report**. Other data types including result tables and free texts can be imported to the **Report** or to the current **Project**. Multiple spectra can be selected for loading by holding the SHIFT or CTRL key when clicking onto the file list.

**Saving data to files.** To save ESPRIT items (spectra, images, etc.) open the Import/Export menu by clicking the  icon in the according spectrum chart or image chart. Decide whether a graphic representation or data file is appropriate and click **Save** below the corresponding headline. A browser for saving files will open. The same selection of volumes and folders is provided as for loading data.

In the file browser enter or edit the file name and check the desired data format.

**Graphical data output.** Graphic export is possible using common graphical file types or the Windows® clipboard. To export graphical data to files use the option **Save** under the subsection **Graphic (Image, Map image)**. To exchange graphics via the Windows® clipboard, click **Copy** in the according subsection of the Import/Export menu or in the local menu after a right mouse click into the corresponding window area.

**Projects.** Storing data in a project and saving the whole project to a project file is the recommended way of storing data. Working with projects is described in section 5.22.

**Result tables/lists.** Qualitative and quantitative results are automatically included in spectra files. To save result tables separately, open the spectrum chart Import/Export menu ( icon) and **Save** under Result list/tables or use the local popup menu after a right mouse click into the result table.

**Element selection.** Element selections are a special data type within QUANTAX that control the display of line markers and provide element lists for spectrum quantification, line scan and mapping. Element selections can be saved to file, transferred between workspaces using **Copy** and **Paste** or added into a project (see also section 5.22).

**Twain export.** Images or other data from ESPRIT can be exported via the standardized interface Twain export. It is enabled only when data is being requested by image processing applications.



## 4.7 System Functions



### Attention!

Any changes in the system settings will influence the global ESPRIT software. The user must proceed with caution when altering the system settings!

Clicking the **System** button in the main menu gives access to the **System** workspace. This workspace is divided into seven sections (tabs): **System**, **Appearance**, **Spectrometer**, **Imaging**, **WDS**, **Microscope** and **Stage**.

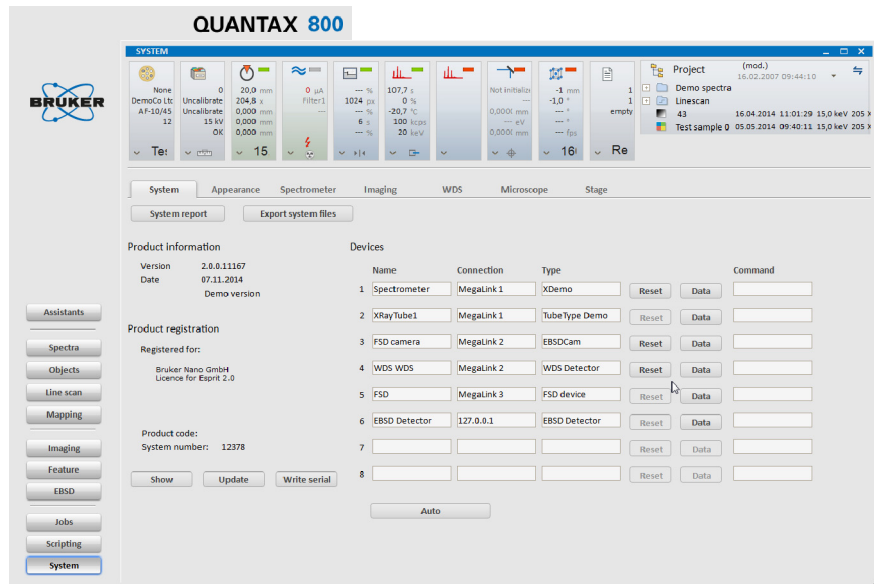
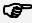


Fig. 4.7-1 Workspace System

### 4.7.1 System

The section **System** provides product information, user license administration, and control of the spectrometer settings. All items in this section are generally only of interests to Bruker service personnel. Unless explicitly requested, please do not alter any parameters within the system assignments.

**System report.** This area lists all relevant system settings. The report is valuable for troubleshooting and service. If requested by Bruker service personnel, the report can be exported via the  icon.

 The product registration is only valid for the specified system number. The system number of the current hardware is also listed in the product registration area.

**Export system files.** This dialog allows for export all system relevant setting files as a .zip file.

**Product information.** The current version of the ESPRIT software is displayed here. Please relay this information to Bruker Service upon service enquiry.

**Product registration.** The product registration area shows the owner's registration information. Further details can be accessed by clicking on the **Show** button. This will launch the PDF reader to show the current product license. If a new license file is issued, the file in the ESPRIT software can be updated by clicking the **Update** button. A copy of the license file is stored to the QUANTAX server's profile folder. The license will be checked automatically by every time the ESPRIT software is started.

**Devices.** The assignment and setup of the X-ray spectrometer(s) and other devices are normally only performed during system installation.

The assignment is completely automatic. Each detector to be assigned must be connected to either Bruker MegaLink port, a RS232 port, an USB port or via Ethernet of the QUANTAX server and powered on when the **Auto** button is clicked. All available ports are automatically scanned for valid connections. Once the spectrometer is located, the ESPRIT software must be restarted in order to register the newly found device.

After the restart of the ESPRIT software the available detectors are listed within the **Configurator bar** (refer to section 4.5).

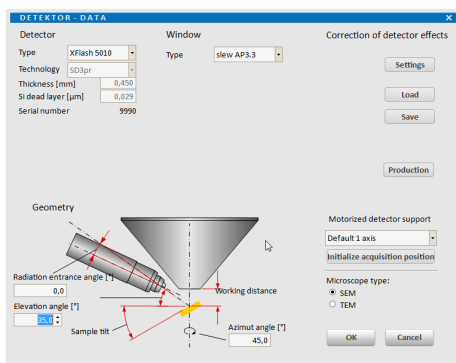


Fig. 4.7-2 Detector data dialog

Clicking the **Data** button opens a dialog for setting the type of spectrometer, the type of the radiation entrance window, and the geometry parameters. These entries are loaded from the detector description file or the detector data sheet.

The motorized detector support can be activated here. In order to record the acquisition position, click the **Initialize acquisition position** button and follow the wizard.

The **Command** field next to the **Data** button performs low-level command functions. This is for troubleshooting purposes by system engineers.

## 4.7.2 Appearance

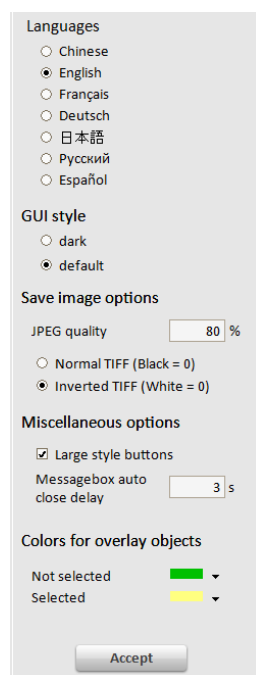


Fig. 4.7-3 Options in the Appearance tab

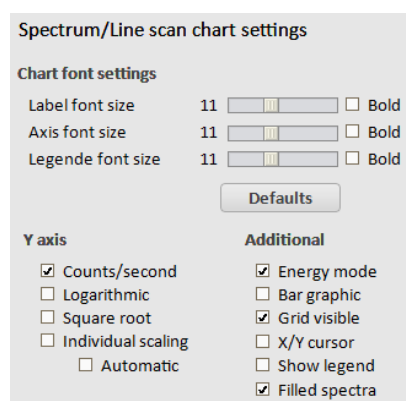



Fig. 4.7-4 Spectrum/ Line scan settings

The **Appearance** tab allows the user to configure settings related to the display of the ESPRIT software.

The subsection **Languages** allows the user to change the display language. Select a language and confirm by clicking **Accept**.

The **GUI style** allows the user to change the ESPRIT interface to the dark color scheme for use in low light conditions (e.g. in TEM rooms).

The subsection **Save image options** permits the user to reduce the quality of the images saved in jpeg format, in order to keep the file size low. For images saved in TIFF format, two types of gray palettes can be selected: **Normal** TIFF assigns black to 0, and **Inverted** TIFF assigns white to 0.

In the subsection **Miscellaneous options** the user can activate/ deactivate the **Large style buttons** mode. If the checkbox is activated, the name of the workspace or function is shown on the icon, e.g. . The **Auto close delay** option defines the time the mouse-over hint messages of the icons are displayed.

The **Colors for overlay objects** in different workspaces can be selected in this subsection. Overlay objects are objects (points, lines, rectangles, ...) to e.g. extract and display spectral information from the mapping data cube.

In the **Chart settings** tab, font sizes and properties of the spectrum chart can be modified.

To use a fixed color setting for a particular chemical element, click on the element symbol on the periodic table to change the elements properties. If no specific color is defined, a random color is assigned.

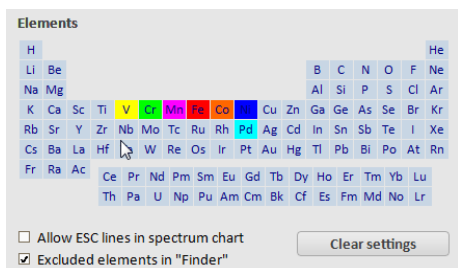


Fig. 4.7-5 Periodic table

The option **Excluded elements in 'Finder'** should be checked by default. This allows to see all possible elements when using the finder, even those that are excluded for the Auto ID (the Auto ID works best especially for noisy spectra, if the user excludes some elements that are not expected in the sample). The box should only be unchecked, if the excluded elements should not be shown in the Finder.

In the **Table settings** tab, the user can specify the default settings for the export of tables.

In the **Spectrum tables** tab, it is possible for the user to select the type of information the quantification table should display post quantification.

### 4.7.3 Spectrometer

#### Attention

The spectrometer calibration is stored in the hardware and affects all subsequent uses and users.

Click the **Spectrometer** tab to access the screen area for spectrometer calibration.

Spectrometer calibration refers to the realignment of the energy axis (abscissa) on the EDS spectrum using a calibration sample (more later) with known composition. The calibration is recommended for high precision quantitative analysis, or in instances where the line markers do not match the peak positions, or if the auto identification fails.

**Pulse throughput and Range selection.** The user can choose the different combinations of pulse throughput and energy range for spectrometer calibration. For each combination, the calibration data is maintained separately.

Do not interfere while the automatic calibration is running. If necessary, click the **Cancel** button to terminate the sequence.

**Calibration.** Follow section 5.2 to perform energy channel calibration. It is possible to return to predefined settings by clicking the **Factory defaults** button if the peaks in the spectrum deviate severely from theoretical values.

The calibration timing schedule should be adjusted to the laboratory regulations or practice and constraints defined by the application.

**Calibration report.** The calibration report lists all relevant calibration settings of the QUANTAX system. If requested, please provide this document to the Bruker service personnel.

### 4.7.4 Imaging

The **Imaging** tab gives access to image and electron microscope (EM) stage calibration.

**Image input.** The image input links the communications between ESPRIT software and the EM. The settings here are microscope dependent and should only be modified by Bruker service personnel.

**Image calibration.** The image calibration is performed to calibrate the pixel values within the ESPRIT software. It is possible to calibrate based on the electron microscope scale bars, but user should ensure that the pixel values on the microscope is up to date before performing image and stage calibration in the ESPRIT software.

A standard calibration grid or any specimen that has features of known geometry can be used for image calibration. After capturing an EM image, locate the features of the image by the green overlay line. To adjust the line, click on the line once to show the sizing handles, then drag the handles to fine tune the position of the end points of the line.

Once confirmed, enter the correct/certified values of the line length in  $\mu\text{m}$ . Click the **Accept** button to store the according calibration value.

**Stage calibration.** This option calibrates the stage movement alongside the SEM image for montage purposes. Before stage calibration, user must ensure that the line length (pixel calibration) is properly calibrated.

To start calibration, capture a SEM image, then open the acquisition settings by clicking on the  $\nabla$  icon in the Acquire button. The settings that can be changed here include the percentage of overlap (it is advised to use the same overlap value for actual montage acquisition in other modes), 2 x 2 or 3 x 3 images. Click on the  $\nabla$  icon again to confirm the settings.

For best results, it is recommended to switch off the scan rotation on the SEM to ensure that the stage movement matches the scan direction. Click the **Acquire** button to start image acquisition. A progress bar will show the progress of the image acquisition.

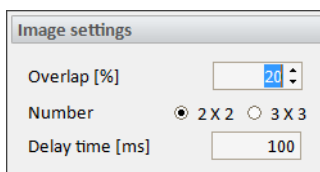


Fig. 4.7-6 Stage calibration setup

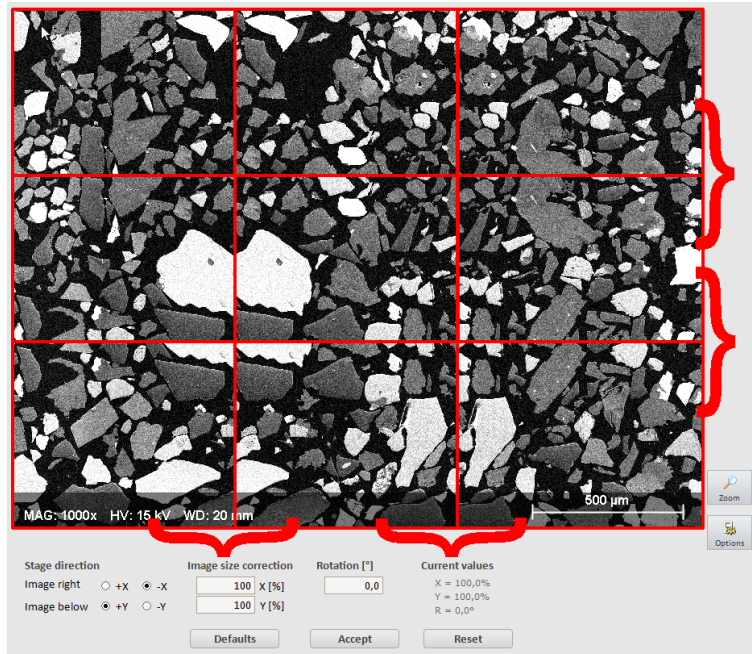


Fig. 4.7-7 Stage calibration

Once all images have been acquired, fine tune the overlap in the X/Y directions or the rotation and press ENTER on the keyboard. The correction in X will only change the overlap in the X-direction, the same applies to Y. The rotation will change the relative rotation between adjacent fields. In the image, the red curly brackets highlight the misaligned overlaps. Adjust the overlaps and rotation in small steps. Care should be taken to focus not on the corners of the image overlaps but at the straight edges (in the corners the image distortion is at its worst).

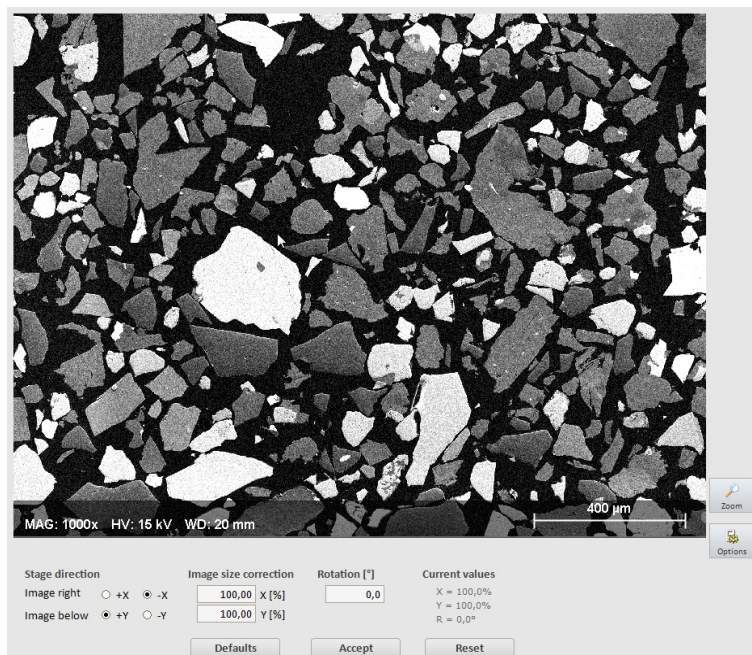




Fig. 4.7-8 Best fit of stage calibration


 It will not always be possible to get a perfect fit, especially with the first frame (top left) as the image may be affected by stage backlash.

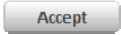
#### 4.7.5 WDS

#### 4.7.6 Microscope

 Microscope data is also controlled via **MICROSCOPE CONFIGURATION** dialog. To enable the microscope data transfer, the **Communication active** checkbox must be ticked.


#### 4.7.7 Stage

 The limits must not be changed without informing Bruker Service personnel.

Once the user has determined the best fit after adjustment, confirm it by pressing the  button. The **Current values** will then be updated.


For details refer to the QUANTAX WDS User Manual and the XSense Reference Manual.

The microscope communication settings are fixed during installation of the QUANTAX system. Image channel names are set to "Ch 1" and "Ch 2" by default. These names can be changed here to EM-specific channels (e.g., SE, BSE, HAADF,...). It is not advised for users to modify the other settings in this section unless explicitly advised by Bruker Service personnel.

To verify the stage limits, select the  tab. The stage limits for the SEM are set during installation of the QUANTAX system.

## 4.8 The Help System

**Help system.** ESPRIT features a multi-level help system. Descriptions are provided for all major control elements. Dialog boxes pop up, if manual interaction is required. Warnings to prevent performing undesired actions are displayed to the user.

**Program help.** The comprehensive user manual can be accessed from the program window using the  icon.

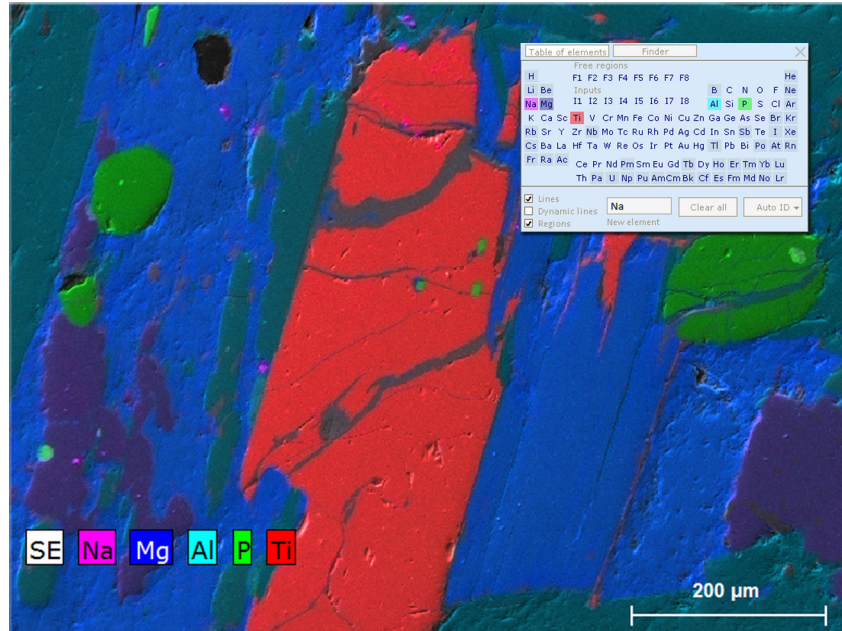
**Context sensitive help.** The mouse cursor is context sensitive. When moving the mouse over objects, e.g. buttons and labels containing short explanations are displayed.

**Assistants.** Interactive assistants provide step-by-step instructions for common analytical tasks.





# 5 Step-by-Step Guides



5.1	Preparatory Steps.....	51
5.2	Energy-Channel Calibration .....	53
5.3	Image Capture.....	55
5.4	Image Extension .....	57
5.5	Drift Correction.....	59
5.6	Image Processing.....	61
5.7	Spectrum Acquisition .....	63
5.8	Using the Spectrum Chart.....	65
5.9	Sample Coating Correction .....	68
5.10	System Factor Calibration/ Beam Current Setting .....	70
5.11	Standards Library .....	73
5.12	Quantitative EDS Analysis of Bulk Samples .....	81
5.13	Quantitative EDS of Electron Transparent Samples .....	90
5.14	Object Analysis .....	96
5.15	Line Scan.....	98
5.16	Quantitative Line Scan (QLine).....	100
5.17	Map Acquisition .....	102
5.18	Quantitative Mapping (QMap).....	104
5.19	Map Processing .....	106
5.20	Phase Analysis .....	109
5.21	Creating a Report .....	112
5.22	Managing Projects .....	114






*This chapter gives the analyst an overview of the most important tools of the ESPRIT software, permitting a quick start to perform analyses with the QUANTAX EDS system.*

*The procedures described here assume standard (default) ESPRIT settings. They also refer to the use of a scanning electron microscope (SEM) or a scanning transmission electron microscope (STEM) or similar. Differences in the analysis of bulk samples (mainly SEM and FIB) and electron transparent samples (TEM, STEM, sometimes SEM or FIB) are pointed out.*

## 5.1 Preparatory Steps

This section describes the actions needed prior to starting the analytical work with the QUANTAX EDS system.

<b>Step</b>	<b>Examples/hints</b>
<b>1</b> Prepare the sample and place it in the microscope	<i>e.g., embedding, polishing and carbon coating, TEM sample preparation</i>
<b>2</b> Adjust high voltage and working distance (SEM) or eucentric height (TEM) on the microscope. Switch on the beam. Adjust magnification.	<i>Choose high voltage (HV) based on overvoltage of elements to analyze.</i>
<b>3</b> Start ESPRIT and log in.	<i>Refer to section 4.1.</i>
<b>4</b> Use the  icon in the bottom left corner of the <b>Microscope</b> configurator to open the <b>MICROSCOPE CONFIGURATION</b> dialog and check the settings.	<i>WD, Magnification, HV values must match values set on the microscope (see section 4.5.3). Device status indicator should be green.</i>
<b>5</b> Use the  icon in the bottom left corner of the <b>EDS</b> configurator to open the <b>EDS CONFIGURATION</b> dialog and <ol style="list-style-type: none"> <li>a) Set <b>Pulse throughput</b> and <b>Maximum energy</b></li> <li>b) Check <b>Detector temperature</b>.</li> </ol>	<i><b>Normal operation mode</b> should be selected. For <b>Pulse throughput</b> and <b>Maximum energy</b> use <b>Automatic</b> settings. Change it on demand.</i>
<b>6</b> Check input count rate and dead time for the analysis.	<i>Check operation temperature. Device status indicator should be green (see section 4.5.5).</i>
<b>7</b> Use the  icon in the bottom left corner of the <b>Sample</b> configurator to open the <b>SAMPLE PROPERTIES</b> dialog and set sample name and description.	<i>The input count rate should be chosen according to the analytical task.</i>
<b>8</b> Proceed with required analysis.	<i>The sample name applies for all further measurements.</i>
	<i>Refer to section 5.7, 5.14, 5.15 or 5.17.</i>

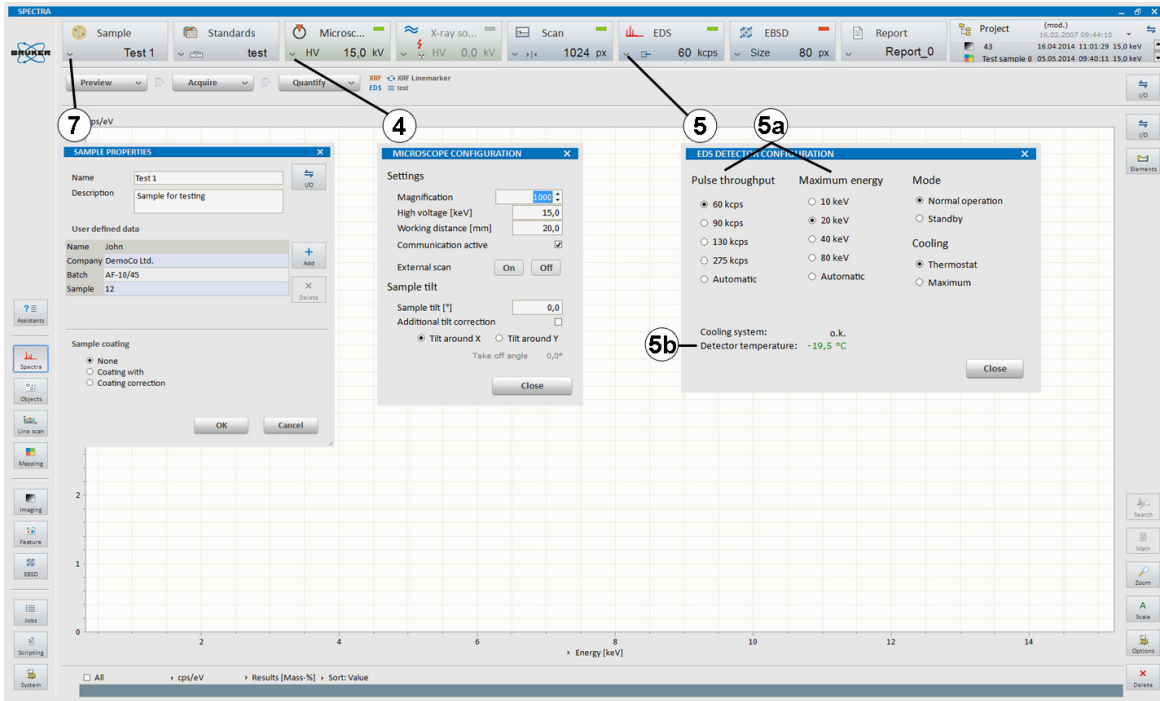
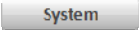
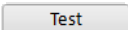





Fig. 5.1-1 Preparatory steps for measurements with the ESPRIT software

## 5.2 Energy-Channel Calibration

Correct energy-channel calibration is a prerequisite for reliable qualitative and quantitative analysis results. Calibration should be checked on a regular basis or after re-start of the signal processing unit (SVE).

Step	Examples/hints
<b>1</b> Select the  workspace and the <b>Spectrometer</b> tab.	<i>Do not confuse energy-channel calibration with system factor calibration (relevant for standard-based quantification; for details see section 5.12.3 and 5.12.4).</i>
<b>2</b> Set a calibration sample into the analysis position.	<i>Use a sample with any element in the mid energy range (Mn – Zn) without peak overlap and sufficient concentration. Preferably use single element standards (e.g. pure copper) to avoid peak misidentification, improve speed and accuracy.</i>
<b>3</b> Set the beam current to produce an intermediate count rate.	<i>Use microscope controls. Set appropriate magnification on a homogeneous sample area.</i>
<b>4</b> Select element and spectral line for calibration.	<i>Use <math>K\alpha</math> lines whenever possible.</i>
<b>5</b> Click  to check the current energy channel calibration and deviation.	<i>The test corresponds to the current <b>Pulse throughput</b> and <b>Maximum energy</b> displayed in the <b>EDS configurator</b>. The acceptable deviation value is between + 5 eV and - 5 eV. If not, proceed with step 6.</i>
<b>6</b> Select <b>Pulse throughput</b> and <b>Energy range</b> under  .	<i>Recommended value is <b>Medium</b>.</i>
<b>7</b> Choose accuracy and click  .	<i>The calibration data is not user specific. Changes in calibration will affect all users of the ESPRIT software.</i>
<b>8</b> Click  to save new calibration data.	

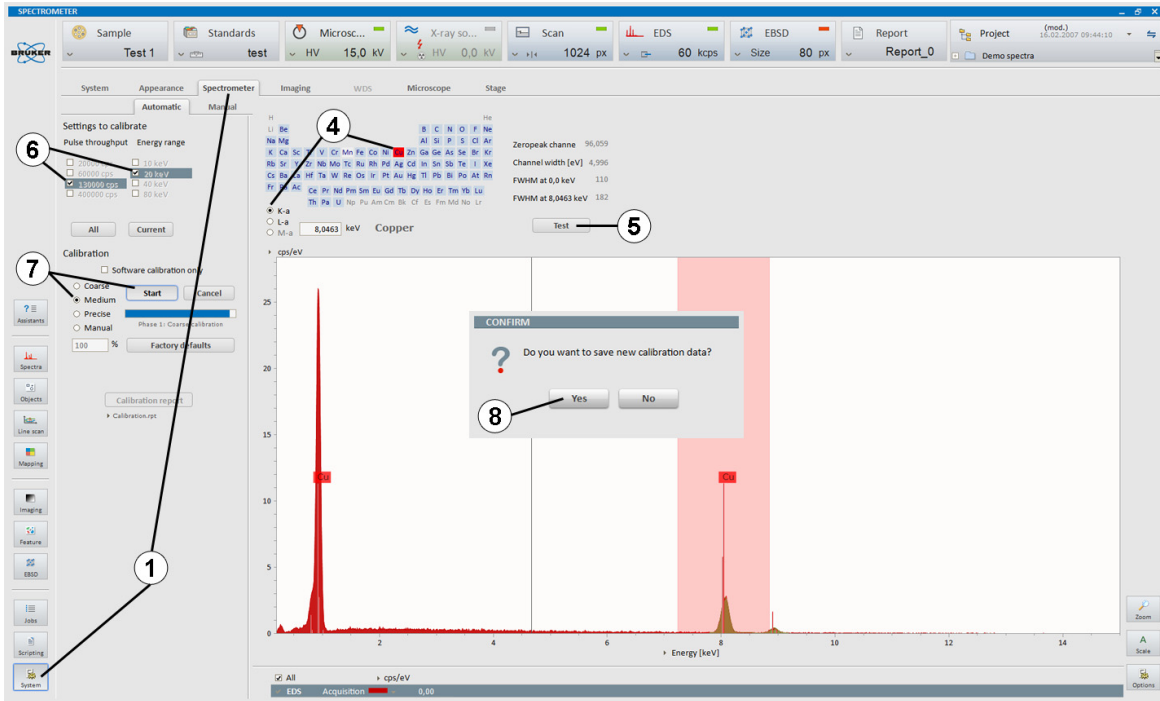

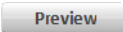

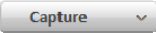
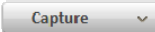
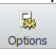
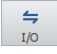


Fig. 5.2-1 Spectrometer calibration

## 5.3 Image Capture

This section describes how to capture electron microscope images with the ESPRIT software.

<b>Step</b>	<b>Example/hints</b>
<p><b>1</b> Select a workspace (<b>Objects</b>, <b>Line scan</b>, <b>Mapping</b>, or <b>Imaging</b>).</p> <p><b>2</b> Use the  icon in the bottom left corner of the <b>Scan</b> configurator to open the <b>SCAN CONFIGURATION</b> dialog.</p> <p><b>3</b> Set image resolution (pixel).</p> <p><b>4</b> Set dwell time [<math>\mu</math>s].</p> <p><b>5</b> Set line average.</p> <p><b>6</b> Click  and set image contrast and brightness on the microscope.</p> <p><b>7</b> Click  on the  button to set capture parameters and <b>Image number</b> or <b>Automatic numbering</b>.</p> <p><b>8</b>  an image.</p> <p><b>9</b> Image legend parameters can be edited using the  icon.</p> <p><b>10</b> Toggle between two image channels if available.</p> <p><b>11</b> Use the image chart  icon to</p> <ol style="list-style-type: none"> <li>a) <b>Save</b> image</li> <li>b) <b>Add image to project</b></li> <li>c) <b>Add image to report</b>.</li> </ol>	<p><i>These workspaces allow capturing an EM image.</i></p> <p><i>The parameters in steps 3, 4, and 5 affect the total <b>Image capture time</b>.</i></p> <p><i>Use histogram settings and oscilloscope overlay tools if needed.</i></p> <p><i>The image will have the name set in the <b>Sample</b> configurator. Choose <b>Single</b> to acquire one image frame, <b>Continuous</b> to update image frames or <b>Sliding average</b> to average a number of frames.</i></p> <p><i>The image name can also be changed here after capture.</i></p> <p><i>Note that the image channel names can be changed in the System workspace under Microscope to e.g. SE, BSE, HAADF... (see section 4.7.6).</i></p> <p><i>Alternatively, click with the right mouse button into the image or drag and drop it to <b>Project</b> or <b>Report</b>. To perform image processing, refer to section 5.6.</i></p>

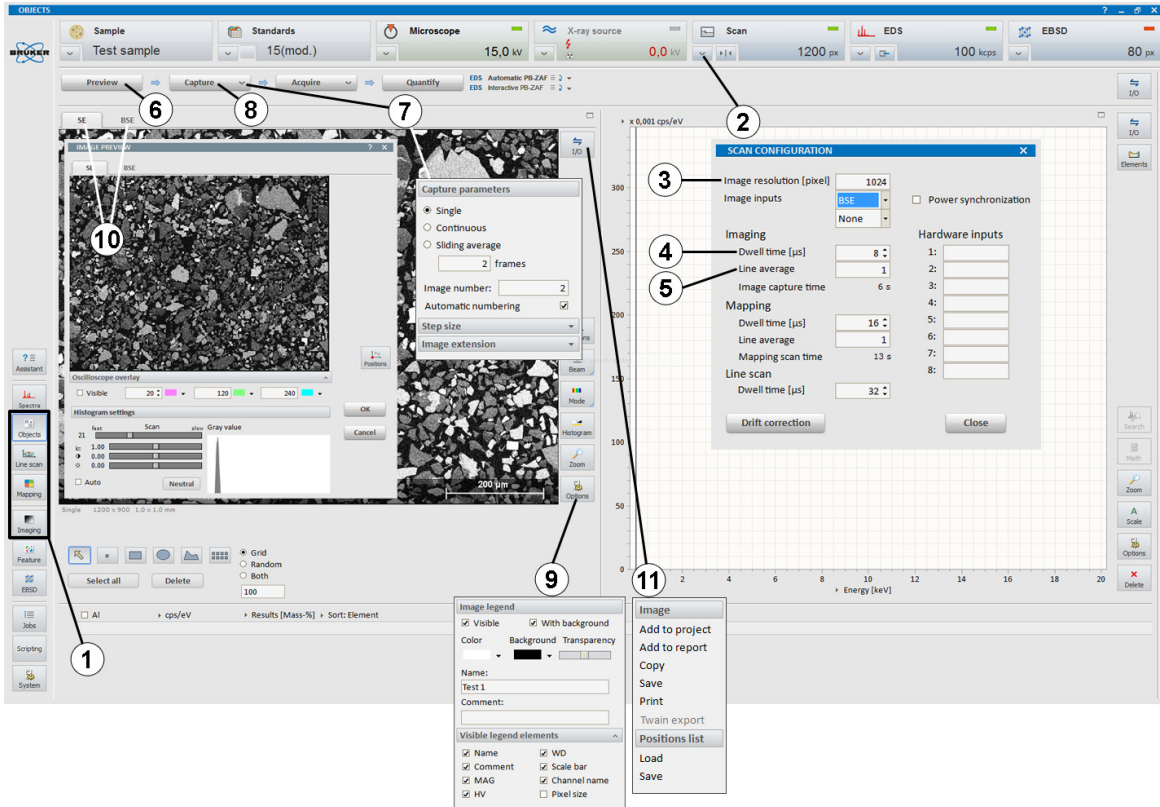


Fig. 5.3-1 Image capture





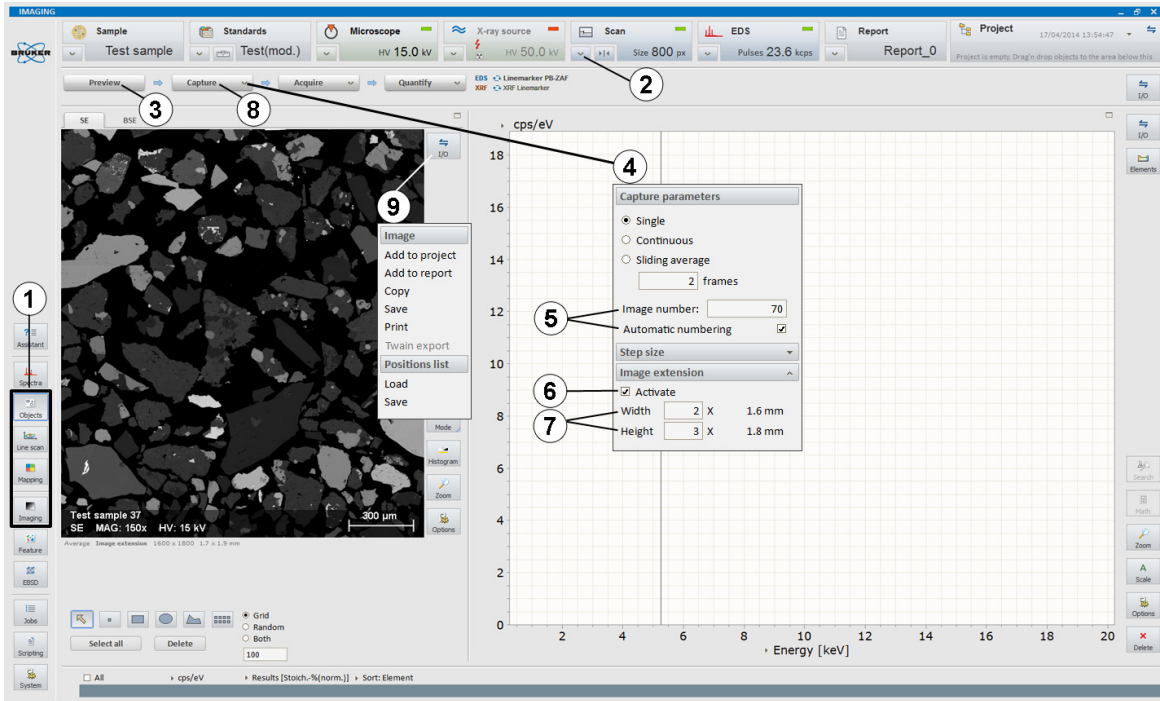


Fig. 5.4-1 Image extension



## 5.5 Drift Correction

The drift correction tool compensates sample drift when performing long and/or high-resolution acquisitions.

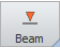
### Step

- 1 Select a workspace (**Objects**, **Line scan**, **Mapping**, or **Imaging**).

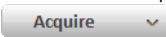
Set measurement conditions and before starting the analysis proceed as follows:


- 2 Use the  icon in the bottom left corner of the Scan configurator to open the **SCAN CONFIGURATION** dialog.
- 3 Click  to open the **IMAGE DRIFT CORRECTION SETTINGS** dialog.
- 4a When performing a Line scan or Objects analysis, select **Current input** and **Correction interval**.

- 4b When acquiring a Map, the correction time is equal to the **Mapping scan time**.

- 5 Click on  to change parking position of the beam.

- 6 Click on  to activate drift correction.

- 7 Start data acquisition by clicking the  button in the current workspace.

- 8 Right click on  to display current image shift.

### Example/hints

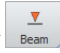
*These workspaces allow capturing an EM image and beam control.*

*Refer to the guides in section 5.3, 5.14, 5.15 or 5.17. When acquiring a Map, select a smaller map area than the captured image. Use **Fixed** or **Variable** options under **Map area** (see step 6 in 5.17).*

*Use **Current input** and 60 seconds as initial value for **Correction interval**. If the analysis stops due to extensive drift, decrease this value. Use **Current value** as **Dwell time** to capture the same quality image as during the analysis or use a lower **Special value** than used for Imaging to save time.*

*If image drift is extensive, reduce **Image capture time** by decreasing **Mapping Dwell time** and **Line average**.*

*Use a parking position outside of the analysis area to avoid beam induced contamination.*

*Right click  to change parking conditions.*

*The system starts acquisition and captures an image after each correction interval set in Step 4.*

*The latest correction image, the latest calculated drift vector (arrow) and the quality of drift correction are displayed. When the drift exceeds the limit which can be corrected by the system, the acquisition stops automatically. Acquired data remains in the workspace.*

*The starting image records the initial acquired image when the acquisition starts. The image is then compared at intervals set by the user. The quality plot records the pixel shift in X and Y directions of the image obtained at the latest interval as compared to the original image.*

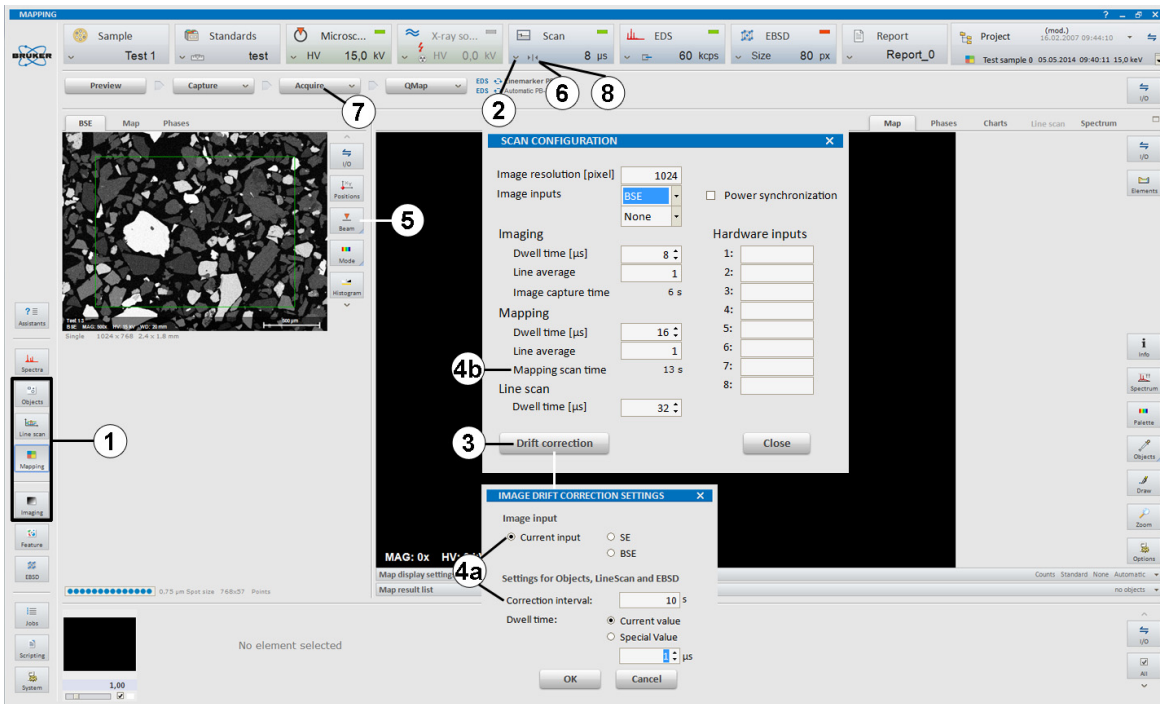
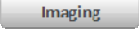
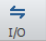
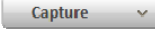
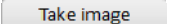

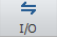
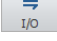
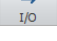


Fig. 5.5-1 Image drift correction

## 5.6 Image Processing

This section describes the workflow for the imaging workspace.

Step	Example/hints
1 Select the  workspace.	<i>Refer to section 5.3. Alternatively, load a saved image by using the workspace  icon of the <b>Imaging</b> workspace or drag an image from <b>Project</b> and drop it into the <b>Edit</b> tab in the image chart. Proceed with step 4.</i>
2  an image.	<i>The captured image appears in the <b>Edit</b> tab and is ready for processing.</i>
3 Select the <b>Filter</b> tab and click  .	<i>Perform filtering and build filter list</i> <i>Perform image masking based on grayscale intensity</i> <i>Offline compilation of mosaic images</i> <i>Perform dimension measurements on images using the  tool and display results as table.</i>
4 Select type of processing from the tab list: <b>Filter</b> <b>Segmentation</b> <b>Panorama</b> <b>Measurements.</b>	<i>Drag intermediate images to the clipboard (shown as thumbnails at the bottom of the workspace).</i>
5 Apply filters or perform other operations.	<i>Use .rti file format to save all available EM images, processed image, result tables, and drawn items into one file.</i>
6 Use the workspace  icon to a) <b>Save</b> processed image b) <b>Add</b> processed image <b>to project</b> c) <b>Add</b> processed image <b>to report.</b>	<i>Alternatively, drag and drop the processed image to <b>Project</b> or <b>Report</b>.</i>  <i>Use .bmp, .jpg, .png, or .tif file format.</i>
7 Use the image chart  icon of the <b>Edit</b> tab to a) <b>Save</b> processed image b) <b>Add</b> processed image <b>to project</b> c) <b>Add</b> processed image <b>to report.</b>	<i>Alternatively, drag and drop the processed image to <b>Project</b> or <b>Report</b>.</i>
8 Use the thumbnail bar  icon to a) <b>Save</b> processed image b) <b>Add</b> processed image <b>to project</b> c) <b>Add</b> processed image <b>to report.</b>	<i>Alternatively, drag and drop the processed image to <b>Project</b> or <b>Report</b>.</i>

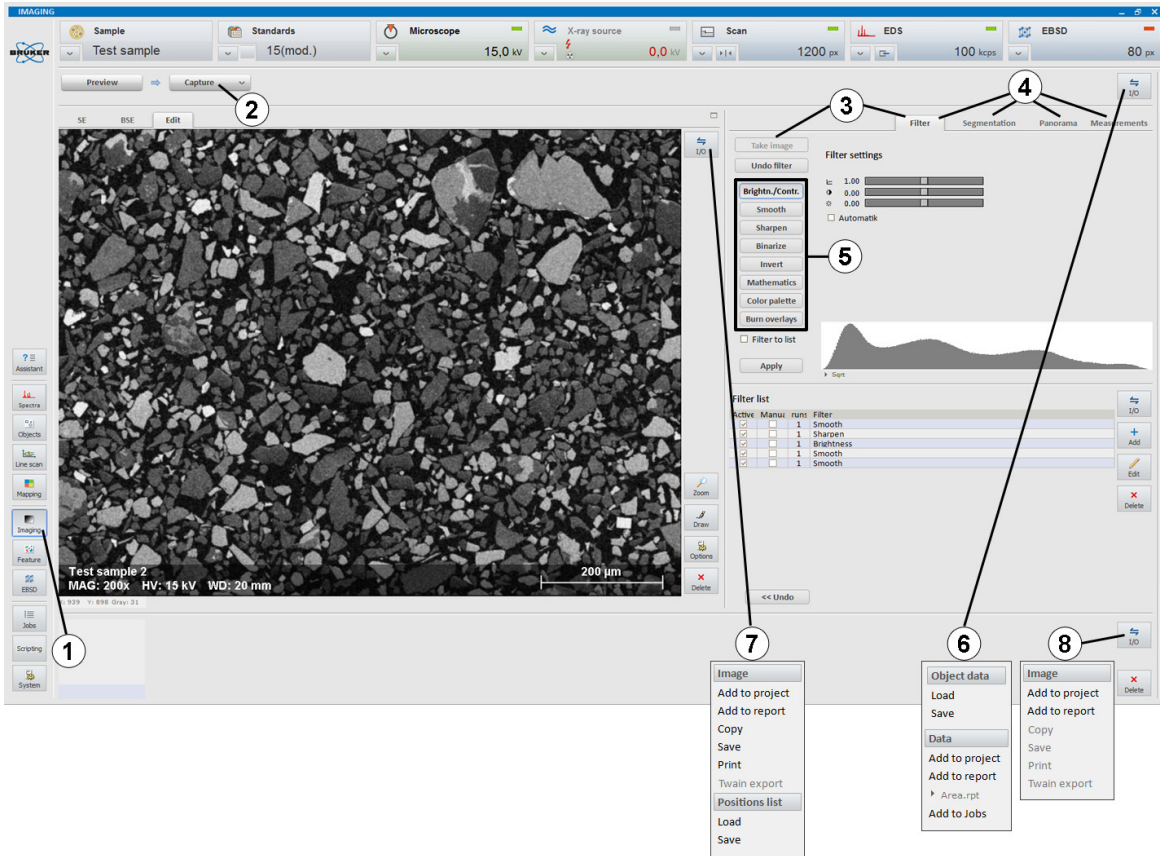


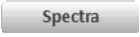
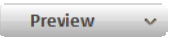

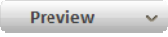

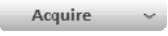
Fig. 5.6-1 Image processing

## 5.7 Spectrum Acquisition

This section describes the workflow for spectrum acquisition within the Spectra workspace. It is recommended to acquire a preview spectrum before proceeding with analysis in various workspaces.

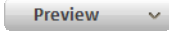
### Step

Choose appropriate measurement conditions on the electron microscope.

- 1 Select the  workspace.
- 2 Click the  button to acquire a live spectrum.
- 3 Click  on the  button to set timing parameters in the **Preview timing** submenu.
- 4 Click  on the  button
  - a) Set **Acquisition parameters**
    - Automatic**
    - Manual**
    - Real Time**
    - Live Time**
    - Counts**
  - b) Set **Automatic quantification**
  - c) Activate **Cyclic acquisition**
  - d) **Spectrum numbering**.


### Examples/hints

Refer to the previous sections 5.1, 5.2, and 5.3.

The live spectrum appears in the spectrum window. Click the  button again to stop acquisition.

Set acquisition time.



**Fast** (50,000 counts) for major elements,  
**Precise** (250,000 counts) for minor elements,  
**Exhaustive** (1,000,000 counts) for elements close to the detection limit.

If **Manual** is selected, the acquisition has to be stopped manually by clicking the  button.

The acquisition will stop after the time entered in the dialog box has elapsed.

The acquisition will stop after the dead time corrected acquisition time has elapsed.

The acquisition will stop after the predefined number of counts is recorded.

If **Continuous/After acquisition** is selected, the spectrum is quantified during/after acquisition. Click  to access the method editor (see description in 5.12.1). Load a quantification method using the  button.

This option acquires several spectra with identical settings from the same area. Set Cycle count and Pause [s]. Use **Add to project** to automatically add acquired spectra to project.

**Automatic numbering** can be activated. The numbering will start with the entered **Spectrum number**.


5  a spectrum.

The acquired spectrum appears in the spectrum chart and the quantification results in the spectrum list (refer to section 5.8). The spectrum can be further processed (element identification, quantification).

6 Use the workspace  icon to

- a) **Add data to project**
- b) **Add data to report.**

Alternatively, drag and drop the spectrum from the spectrum list to **Project** or **Report**. The project has to be saved manually (see section 5.22).

7 Use the spectrum chart  icon to

- a) **Save** spectrum (or result list, element selection, graphic)
- b) **Add item to project**
- c) **Add item to report.**

Highlight spectra in the spectrum list to save several spectra. Alternatively, right click into the spectrum. Various file formats are possible\*. To include all meta data use the .spx format. Refer to section 5.8 and 5.12.1.

Alternatively, drag and drop the spectrum from the spectrum list to **Project** or **Report**. The project has to be saved manually (see section 5.22).

\*optional, license-based

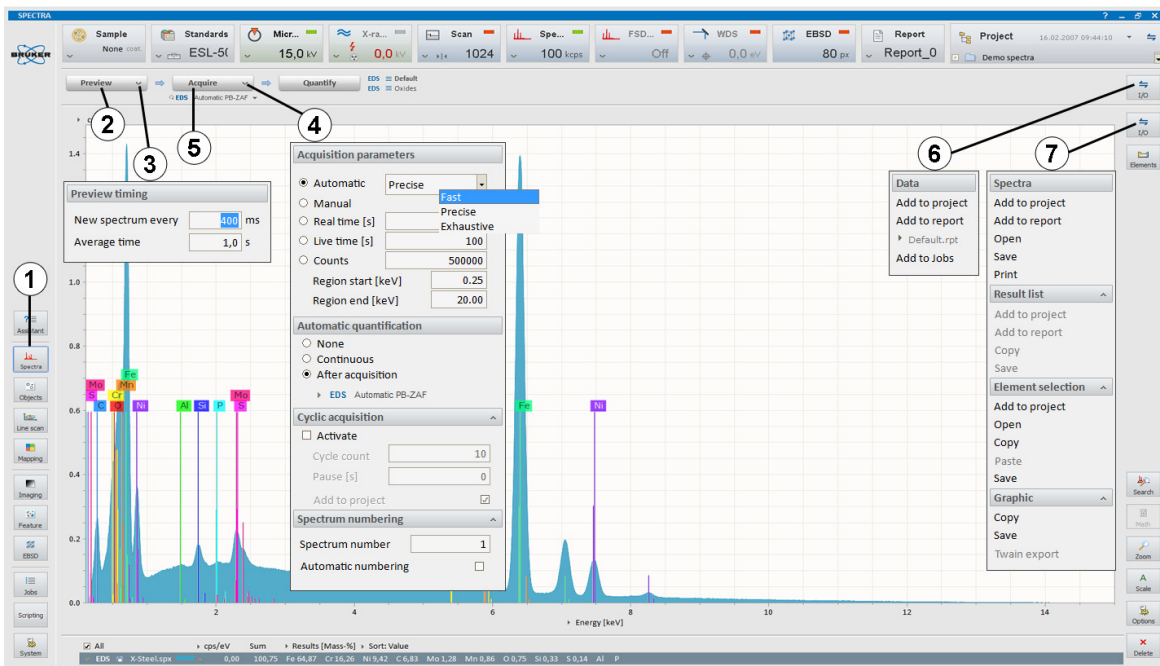


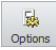






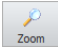
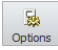


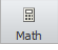
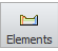


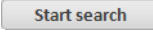
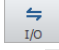
Fig. 5.7-1 Spectrum acquisition



## 5.8 Using the Spectrum Chart

This section describes the features of the Spectrum chart, available in the Spectra, Objects, Line scan, and Mapping workspaces. The assigned numbers do not indicate that a certain order needs to be followed.

Step	Example/hints
The spectrum list shows:	
<b>1a</b> Spectrum type: excitation and/or analysis method	
<b>1b</b> Spectrum name	<i>The spectrum name has to be unique in the spectrum list. It can be changed after clicking on the name.</i>
<b>1c</b> Spectrum color	<i>Click on the color bar to toggle between filled and unfilled display. Click on  to select the color.</i>
<b>1d</b> Options	<i>Click  to select Pulses, cps, Net counts, energy resolution of peak selected by the cursor (FWHM fit).</i>
<b>1e</b> Factor: scaling factor of y-axis	<i>Available when Individual scaling is checked in the  menu.</i>
<b>1f</b> Results.	<i>Click  to select Spectrum information, identification or Display of quantification results (Mass-%, Mass-% (norm.), Atom-%, Stoich.-%, Stoich.-% (norm.)).</i>
<b>2</b> Multiple spectra can be selected (checkbox) or highlighted by clicking on the spectrum line (gray outline).	<i>Checkbox selects spectra for display, highlighted spectra can be processed and exported. To highlight several spectra use SHIFT or CTRL keys and left click on the spectra. When exporting multiple spectrum results, the result table contains methodical error values (Sigma).</i>
<b>3</b> To delete a spectrum press the DEL key on the keyboard or  on the spectrum chart.	<i>Alternatively, drag and drop spectrum from the spectrum list onto the  icon.</i>
To scale spectrum diagram or zoom, use either:	
<b>4a</b> Mouse scroll wheel to change x-scale	
<b>4b</b> Click and hold mouse scroll wheel to move spectrum area	
<b>4c</b> Use CTRL key + left mouse click into the spectrum diagram + drag the mouse to scale x- and y-axis	
<b>4d</b> Right click on x- or y-axis: scale values can be entered manually	

- 4e** Use the  icon for Automatic scaling (spectrum fills display area)
- 4f** Use the  icon to zoom in the spectrum window.
- 5** Use the  icon or right click on the spectrum chart and select **Properties...** to change Spectrum display properties.
- 6** Click  next to the x-axis dimension.
- 7** Click  next to the y-axis dimension.
- 8** Right click on the x or y scale bar.
- 9** Double click on a spectrum in the spectrum list to access Spectrum properties (Results, Sample info, Detector, Parameter, and Spectrum).
- 10** Click the  icon to select the desired operator.
- 11** Click the  icon to select or de-select an element.
- 12** Select the tab  in the periodic table.
- 13** Highlight a region in the spectrum (by dragging the spectrum cursor with the right mouse button pressed) and select a free region (F1-F8).
- 14** Click the  icon on the right side of the spectrum chart and choose the location where the reference spectra are located. Then click .
- 15** Use the spectrum chart  icon to select Add to project in the **Element selection** section.

*Set a rectangle by holding the left mouse key to zoom into the area of interest.*

*To normalize multiple spectra for comparison, select **Individual scaling** and **Automatic** and select an energy region with right mouse button.*

*Toggle between energy (keV) and channel.*

*Toggle between cps/eV or channel and pulse/eV or channel.*

*Adjusting the x and y scaling.*

*Quantification results, sample info, detector (detector parameters), parameter (Acquisition parameters), spectrum (Energy calibration data) can be here retrieved.*

*The **SPECTRA ARITHMETIC** dialog opens. Available from the dropdown list are sum, absolute difference, relative difference, quotient, maximum and minimum. The resulting spectrum will be added to the spectra list.*

*Any element can be selected or de-selected just by clicking the according symbol.*

*The **Finder** option supports the identification of unknown peaks in the spectrum. Place the spectrum cursor over the center of a peak or highlight the peak range by dragging the cursor with pressed right mouse button to display a list of all possible elements. The first element of this list is that one with the highest probability of actually being present in the sample.*

*Up to 8 spectral regions can be assigned to monitor non-analytical peaks or background levels.*

**Free regions**  
F1 F2 F3 F4 F5 F6 F7 F8

***SEARCHING FOR SIMILAR SPECTRA** dialog opens. Found spectra will be added to the spectrum list in the dialog. Use the sensitivity slider (cross correlation factor) to optimize search results.*

*Add the element selection to the project.*

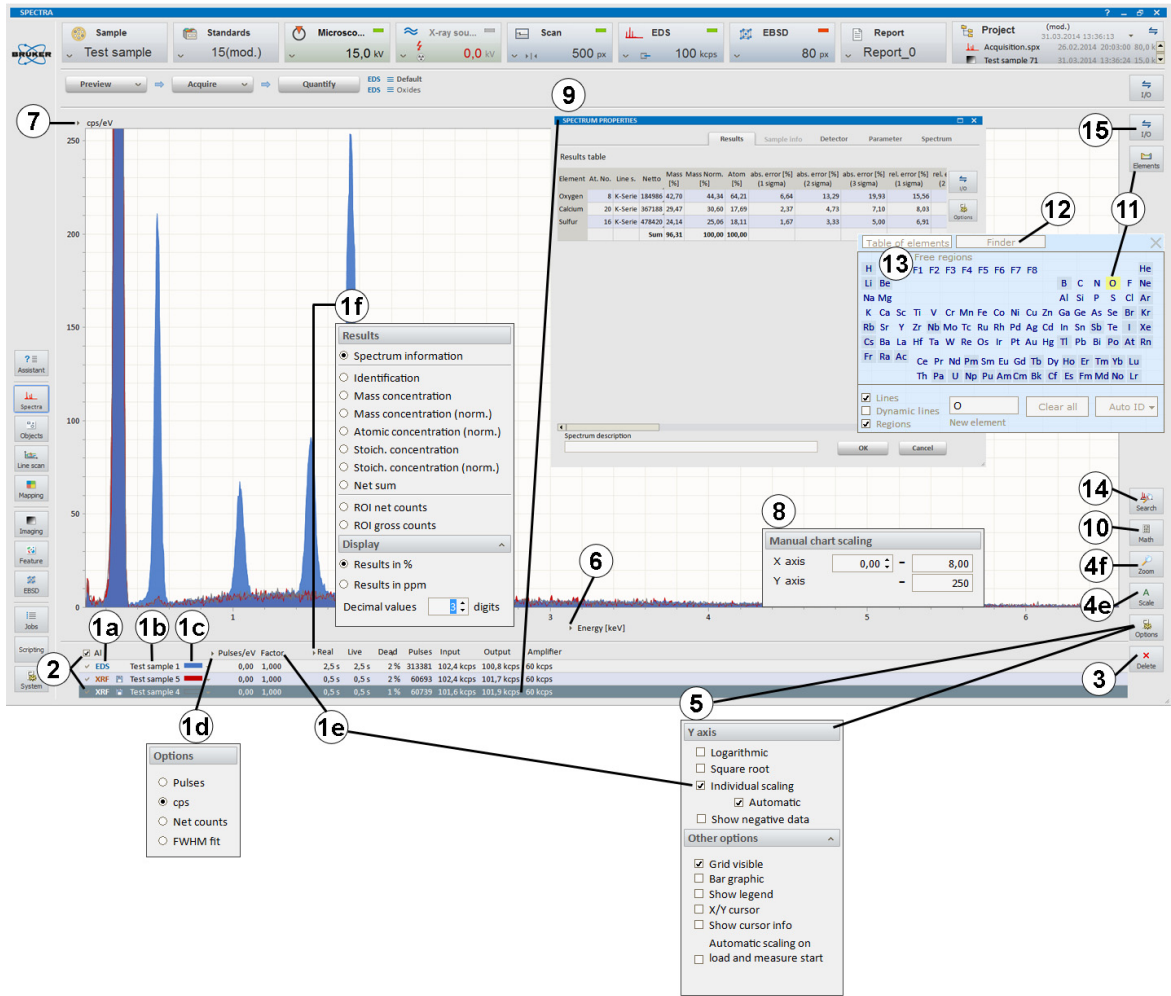


Fig. 5.8-1 Features in the spectrum chart

## 5.9 Sample Coating Correction


This section describes how to use sample coating correction in two ways:

1. Set an element (e.g. carbon) as coating material and omit the net counts originating from it.
2. Measure a reference coating spectrum and subtract this from measured spectra.

### Step

### Example/hints

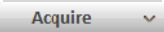
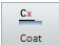
**Set an element (e.g. carbon) as coating material and omit the net counts of this element in the quantification process:**

- 1 Click the  icon in the bottom left corner of the **Sample configurator** to open the **SAMPLE PROPERTIES** dialog.
- 2 Select option **Coating with**.
- 3 Click on an element.

The **Select element** dialog pops up.

The selected element will be deconvolved during the quantification, but not quantified during automatic quantification in subsequently acquired spectra, line scans, and maps.

**Acquire a coating correction file of a coated sample and use this to subtract from all subsequent spectrum acquisitions:**

- 4  a spectrum of a coated sample (as coating reference).
- 5 Perform quantification of the spectrum.
- 6 Click on an element in the result list.
- 7 Click on the  icon.
- 8 Save the .ccc file.


Make sure that the sample does not contain the same element as the coating material.

Refer to section 5.12.2.

The **Save coating correction data** dialog appears.

This file contains the net counts from the coating reference sample.

**Apply the saved .ccc file for coated samples**

- 9 Click the  icon in the bottom left corner of the **Sample configurator** to open the **SAMPLE PROPERTIES** dialog.
- 10 Select option **Coating correction**.
- 11 Load the previously saved .ccc file.

The **LOAD COATING CORRECTION DATA** dialog opens.

Make sure that the coating reference sample and the unknown sample are coated identically.

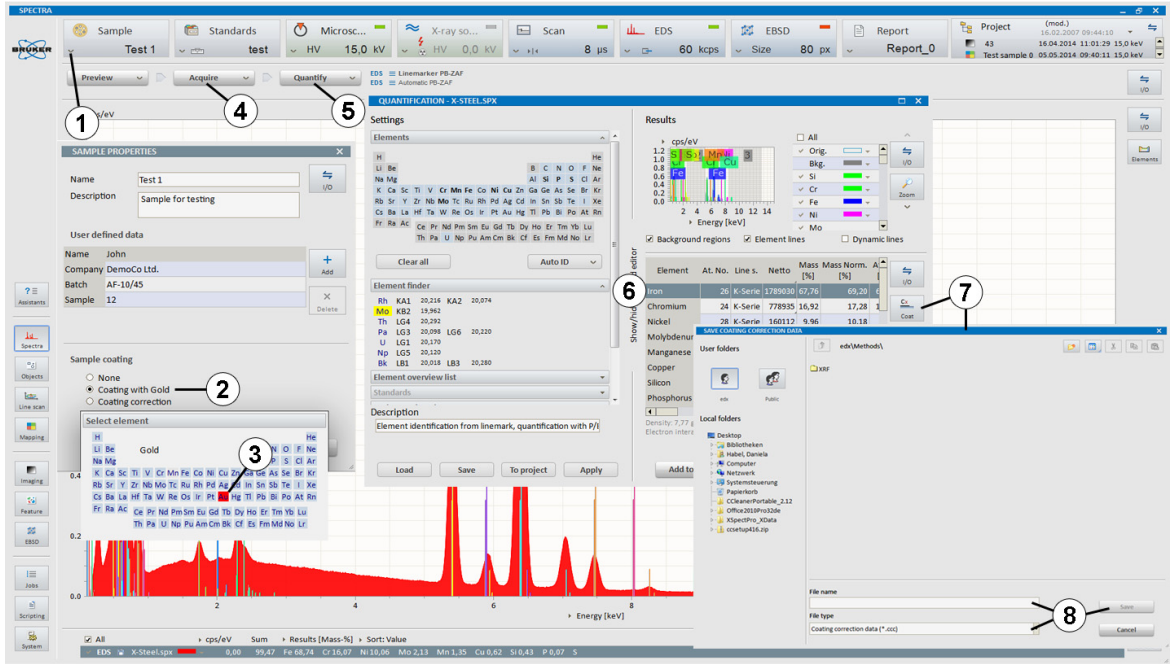


Fig. 5.9-1 Sample coating correction – Step 1 to 8

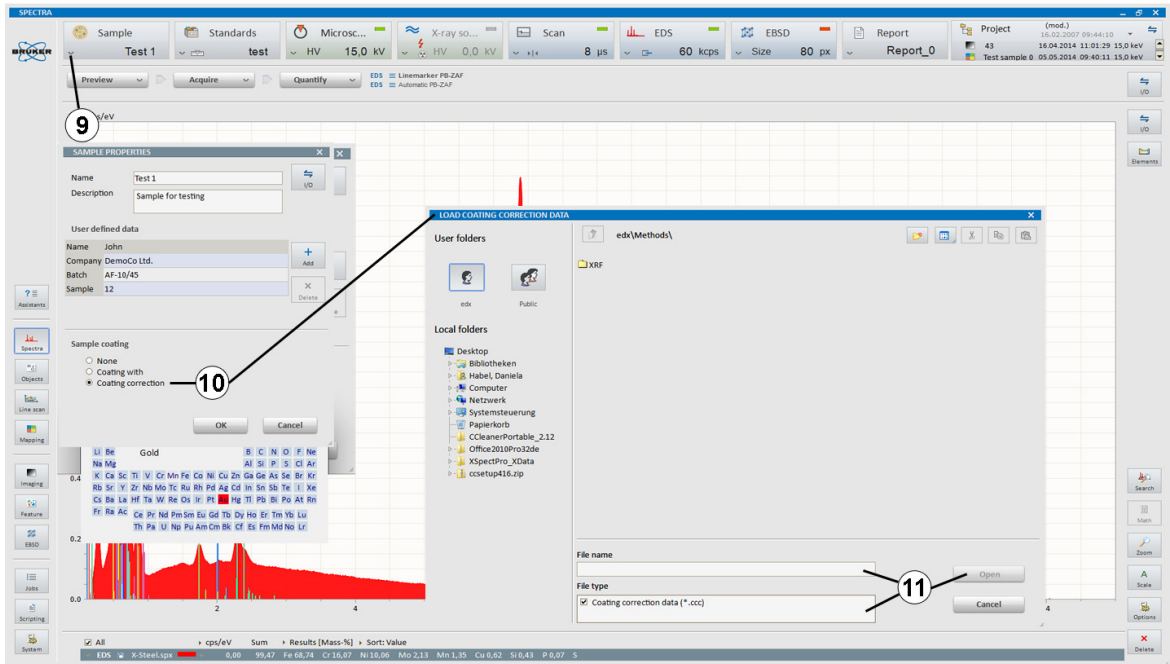


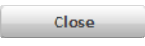


Fig. 5.9-2 Sample coating correction – Step 9 to 11

## 5.10 System Factor Calibration/ Beam Current Setting

For standard-based quantification, a probe current-dependent system factor has to be determined. There are two ways to proceed depending on sample type and whether a picoamperemeter / beam current monitor is available or not. The system factor value is stored in the ESPRIT software and saved along with the following measurements until the software is closed. The quantification of electron transparent samples with the Zeta-factor method requires the knowledge of beam/probe current.

### System Factor Calibration

Step	Example/hints
1 Move the calibration sample into the analytical position and adjust microscope parameters.	<i>Use a single element calibration sample (Co, Mn, Cu). Select a homogeneous area for spectrum acquisition.</i>
2 Load or create a standards library.	<i>Follow section 5.11.1.</i>
3 Click the <b>Standards configurator's</b> calibration icon  to open the <b>SYSTEM FACTOR CALIBRATION</b> dialog.	
4 Select the material of the <b>Calibration sample</b> from the <b>Select the sample to use for calibration</b> drop-down list.	
5 Set <b>Acquisition time</b> parameters.	
6 Click  to calibrate the system factor.	<i>After the calibration is finished the message <b>System is now calibrated</b> appears. Do not change microscope parameters afterwards.</i>
7  the <b>SYSTEM FACTOR CALIBRATION</b> dialog.	<i>Spectra that are obtained after this calibration will contain the measured system factor.</i>
8 Proceed with measurement of standard samples.	<i>Continue with 5.12.3 for standard-based measurement using a temporary reference. Continue with 5.12.4 for standard-based measurement using a standards library.</i>

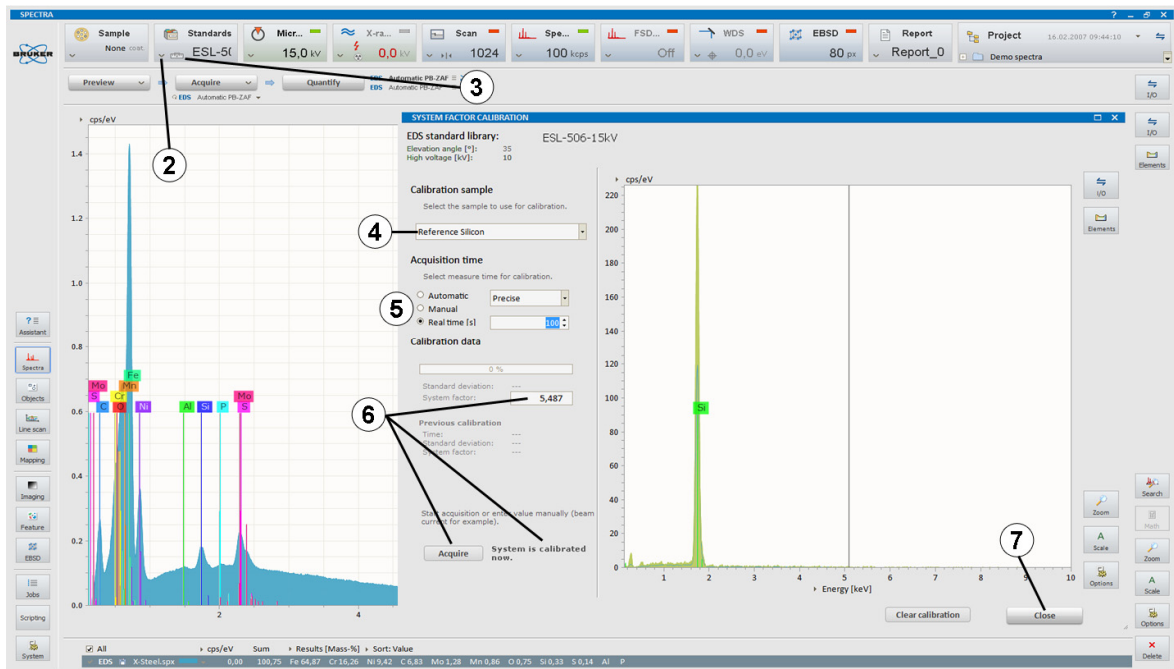

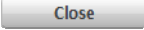


Fig. 5.10-1 System factor calibration

### Beam Current Setting

Use this guide for electron transparent samples and for bulk samples when a picoamperemeter / beam current monitor is available to measure the beam current.

Step	Example/hints
1 Measure beam/ probe current.	Use a Faraday cup or a calibrated electron screen.
2 Load or create a standards library.	Follow section 5.11.1 for bulk samples or sections 5.11.2 and 5.11.3 for electron transparent samples.
3 Click the <b>Standards configurator's</b> calibration icon  to open the <b>SYSTEM FACTOR CALIBRATION</b> dialog.	
4 Enter probe/beam current value manually in the System factor field and press <Enter>.	For Zeta quantification of electron transparent samples, the probe current needs to be known and entered in pA.
5  the <b>SYSTEM FACTOR CALIBRATION</b> dialog.	The system factor is now updated with the beam/probe current and measured spectra will contain this information.

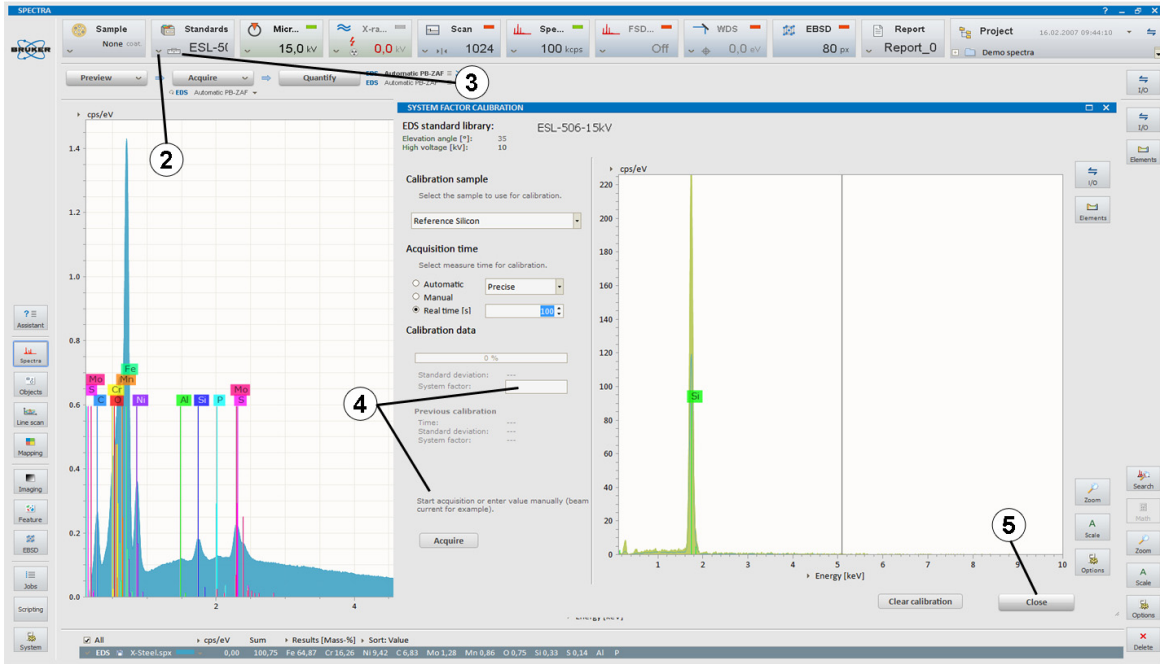




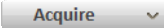
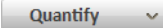


Fig. 5.10-2 Beam current setting



## 5.11 Standards Library

### 5.11.1 Maintain Standards Library for Bulk Samples

This guide describes how to create and maintain a standards library for the standard-based quantification of bulk samples.


Step	Example/hints
<p><b>Create/open a standards library</b></p>	<p>See Fig. 5.11-1.</p>
<p>1 Click the  icon in the bottom left corner of the <b>Standards configurator</b> to open the <b>ADMINISTRATION OF STANDARD LIBRARY</b> dialog.</p>	
<p>2 Use the <b>standards library</b>  icon to</p>	
<p>a) <b>Load</b> a standards library or</p>	<p><i>Load a previously saved *.esl file which corresponds to the current detector geometry parameters. Proceed with step 5.</i></p>
<p>b) Select <b>New</b> to create an empty standards library</p>	<p><i>The question “Do you want to load a spectrum measured with corresponding system settings?” pops up.</i></p>
<p>i. Select <b>Yes</b></p>	<p><i>If you are using the software as Data station (not connected to a microscope). The <b>LOAD SPECTRUM</b> dialog pops up. Load a spectrum .spx which was acquired with the corresponding detector geometry.</i></p>
<p>ii. Select <b>No</b>.</p>	<p><i>If you are using the ESPRIT software connected to the microscope.</i></p>
<p><b>Update standards library with standard spectra</b></p>	
<p>3 Perform system factor calibration.</p>	<p><i>Follow section 5.10.</i></p>
<p>4 Move the reference sample into analysis position and  a reference spectrum (within the <b>Spectra</b> workspace).</p>	<p><i>Follow section 5.7 to acquire a spectrum.</i> <b>Important:</b> Use automatic spectrum numbering or rename spectra after acquisition.</p>
<p>5  the reference sample with a standardless quantification method (within the <b>Spectra</b> workspace).</p>	<p><i>Load a standardless quantification method or deselect <b>Use standards</b> under <b>Quantification model</b> in the method editor.</i> <b>Important:</b> Set the quantification method to  (interactive) to access the quantification dialog. Follow section 5.12.1.</p>
<p>6 Click  in the quantification dialog.</p>	<p>See Fig. 5.11-2. <b>EDIT STANDARD PROPERTIES</b> dialog pops up. Confirm the assignment for all chemical elements to be referenced to that standard. If there is an element missing, it needs to be added in the <b>QUANTIFICATION</b> dialog under <b>Settings/Elements</b>.</p>

7 Add reference concentrations and click **Add to standard library**.

8 Close quantification dialog with **OK**.  
Repeat step 4 to 8 until all standard samples are processed.

**Edit existing standard values**

See Fig. 5.11-1.

9 Click the  icon in the bottom left corner of the **Standards configurator** to open the **ADMINISTRATION OF STANDARD LIBRARY** dialog.

10 Click on an element to select it.

11 Select the standard spectrum assigned to the selected element.

The assigned spectrum appears in spectrum chart.

12 Click **Edit** to change standard concentrations.

The **EDIT STANDARD** dialog pops up, where the assigned concentrations can be changed.

13 Click **Validate** to re-quantify the standard spectrum.

The **QUANTIFICATION** dialog pops up, where the re-quantification of standard spectrum can be performed. Follow section 5.12.1.

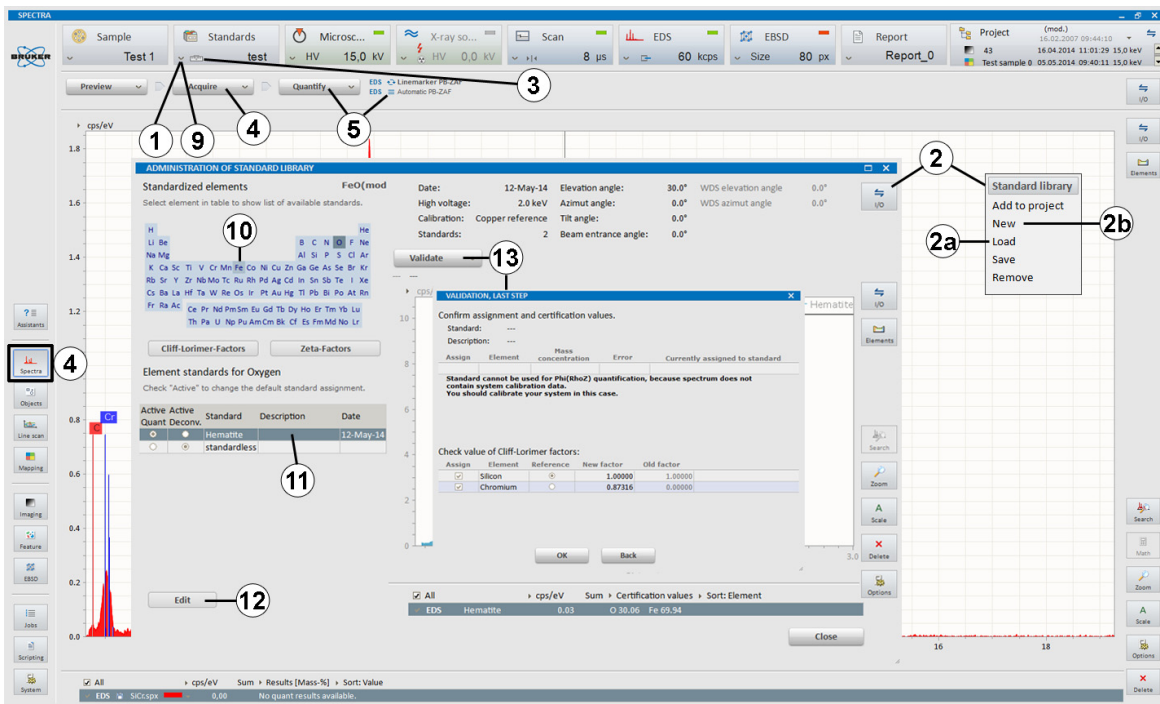


Fig. 5.11-1 Creating and maintaining standards library

**Settings**

Elements: [List of elements]

Element finder: I KBr1 72.294 KBr3 72.218

Element overview list: Standards, Background settings, Deconvolution settings, Quantification model, Additional settings

High voltage [kV]: 200.0

Sample tilt angle [°]: 0

Fast quantification:

**Results**

cps/eV vs Energy [eV]

Peaks: 1 (Cr), 2 (Si), 3 (Cr)

**EDIT STANDARD PROPERTIES**

Name: SiCr  
 Description:   
 Real time [s]: 35.517  
 Life time [s]: 33.790  
 Specification in:  Mass-%  Atomic-%  Stoich-%

Element	Concentration [%]	Error [%]
Silicon	50	
Chromium	50	

Sum of concentrations [%]: 100.00%

Element	At. No.	Line s.	Netto [%]	Mass Norm. [%]	Norm. At. [%]
Silicon	14	K-Series	4340	30.03	34.54
Chromium	24	K-Series	10015	56.92	65.46
<b>Sum</b>			<b>86.94</b>		<b>100.00</b>

Description: Automatic element identification, quantification with P/B-ZAF


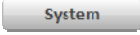
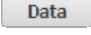

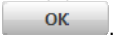
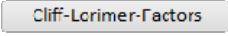
Density: 4.18 g/cm<sup>3</sup>  
 Electron interaction depth: 95.5 µm Radius: 53.5 µm

Buttons: Add to standard library (7), Add as temporary reference, Add to standards (6), OK (8), Cancel

Fig. 5.11-2 Quantification dialog


### 5.11.2 Maintain Standards Library for Electron Transparent Samples – Cliff-Lorimer Factors

This guide describes how to check and calculate theoretical Cliff-Lorimer factors for standardless quantification of electron transparent samples using the standards library dialog.

Step	Example/hints
<p><b>1</b> Click the  icon in the bottom left corner of the <b>Standards configurator</b> to open the <b>ADMINISTRATION OF STANDARD LIBRARY</b> dialog.</p> <p><b>2</b> Check the detector geometry parameters.</p>	<p><i>These parameters have to match the parameters of measured data and/or settings of your EDS configuration on the microscope. A spectrum measured with the correct system settings will also contain the respective values. These can be checked in the file or by double click on the text line below the spectrum chart. The geometry parameters of a measurement system can be checked in the  workspace by clicking on  and should only be changed by Bruker service. If the geometry parameters do not match, proceed with step 3. If they match, proceed with quantification as described in section 5.13.1 or 5.13.2.</i></p>
<p><b>3</b> Use the <b>standards library</b>  icon to</p> <p>a) <b>Load</b> a standards library or</p> <p>b) Select <b>New</b> to create an empty standards library</p> <p>i. Select <b>Yes</b></p> <p>ii. Select <b>No</b>.</p>	<p><i>Load a previously saved *.esl file which corresponds to detector geometry parameters. Proceed with step 5.</i></p> <p><i>The question <b>“Do you want to load a spectrum measured with corresponding system settings?”</b> pops up.</i></p> <p><i>If you are using the software as Data station (not connected to a microscope). The <b>LOAD SPECTRUM</b> dialog pops up. Load a spectrum *.spx which was acquired with corresponding detector geometry.</i></p> <p><i>If you are using the ESPRIT software connected to the microscope.</i></p>
<p><b>4</b> Check <b>NEW STANDARD LIBRARY</b> geometry parameters, enter a name and click .</p>	<p><i>Check or update values of Energy, Elevation angle, Azimuth angle, Tilt angle, Beam entrance angle. They should match the values of your microscope/detector configuration (see step 2).</i></p>
<p><b>5*</b> Click on  to review.</p>	<p><i>The <b>EDIT CLIFF-LORIMER FACTORS</b> dialog pops up: Cliff-Lorimer factors for K-, L- and M-lines are listed. Zoom the graph: left click into it and use the arrow keys on keyboard.</i></p>

6\* Click on a Cliff-Lorimer value to edit it manually.

*It is advised to calculate theoretical factors or set standards as described in section 5.13.2.*

7 Use the **standards library**  icon to save the standards library.

*Save the standards library (containing Cliff-Lorimer factors) as .esl file.*

\* steps can be omitted

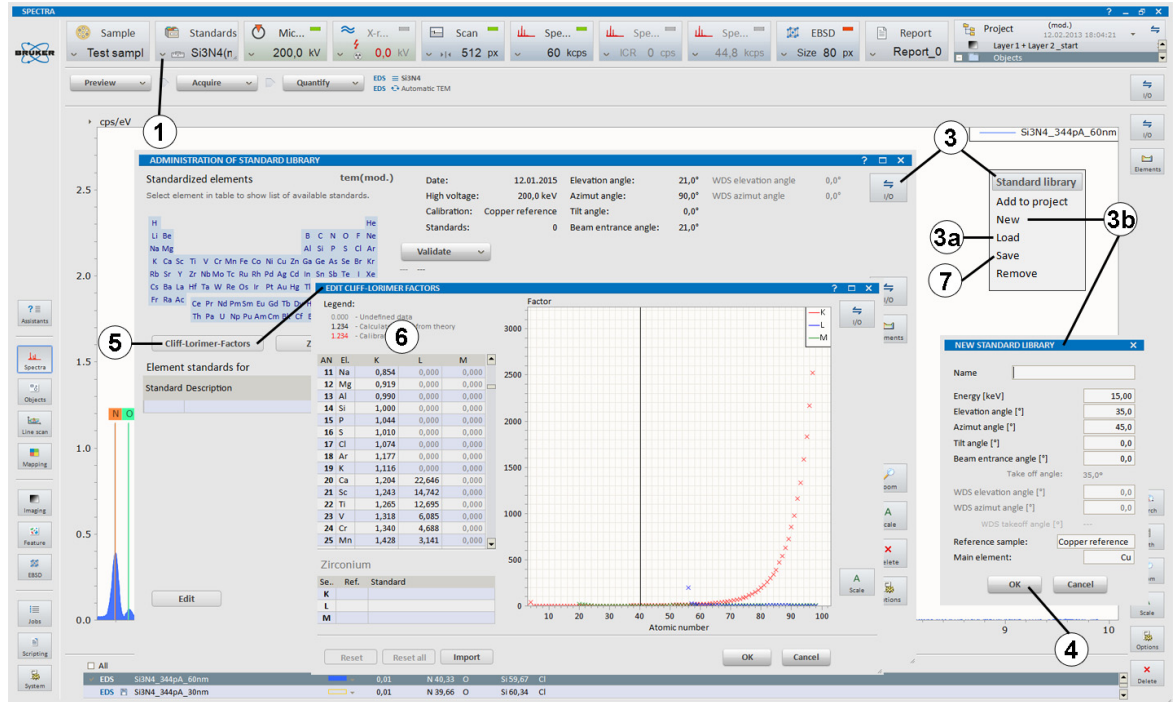


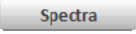

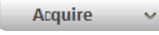
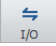
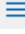

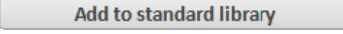
Fig. 5.11-3 Maintain standards library for Cliff-Lorimer factors

### 5.11.3 Maintain Standards Library for Electron Transparent Samples – Zeta Factors

This guide describes how to set up the Zeta factor standards library. Quantification of electron transparent samples with the Zeta factor method requires knowledge of the beam current in [pA] for standard and “unknown” sample and sample thickness of the standard sample.

#### Experimental Zeta Factors

Follow this guide to update Zeta factors for the elements contained in the standard sample.

Step	Example/hints
1 Set up a Cliff-Lorimer standards library.	<i>Follow step 1-4 in section 5.11.2 since the geometry setup for Cliff-Lorimer and Zeta factors is the same.</i>
2 Select the  or the  workspace and	<i>Acquire or open a spectrum of the reference/standard sample. The composition and thickness of the sample has to be known. The beam current during measurement also has to be known and has to be entered as described in 5.13.3 or, if offline, as spectrum parameter: double click on the highlighted line of the active spectrum in the spectrum list (below the spectrum) to open the <b>SPECTRUM PROPERTIES</b> dialog and enter the beam current in pA as system factor in the <b>Parameter</b> tab.</i>
a)  a spectrum	
b) Use the Spectrum chart  icon to open a spectrum.	
3 Follow the standardless quantification steps 6 to 13 to identify elements, set background regions and deconvolution.	<i>Refer to section 5.12.2. Set quantification method mode to  (interactive) to access the <b>QUANTIFICATION</b> dialog.</i>
4 Select Zeta factor method under Quantification model.	
5 Click  .	<i>The <b>EDIT STANDARD PROPERTIES</b> dialog pops up.</i>
6 Enter reference element concentration values and sample thickness for standard sample.	<i>Review Beam current. If missing or incorrect, enter the correct value and also enter the correct value in the spectrum <b>Parameter</b> tab (see step 2). The density will be calculated internally using the sample composition, if changed manually, the thickness will change, respectively.</i>
7 Click  .	<i>The <b>VALIDATION</b> dialog pops up with new and old Zeta factors. Initial Zeta factors have the same values as theoretical Cliff-Lorimer factors (and are considerably smaller).</i>

8 Click **OK** to close the **VALIDATION** dialog.

The updated Zeta factors for elements contained in the reference sample will be marked red in the Zeta factor list.  
NOTE: Save the updated standards library.

9 Click **OK** to close the **QUANTIFICATION** dialog.

If the "unknown" sample contains other elements than the standard sample, proceed with the "Fit Zeta Factors to Standard" procedure below.

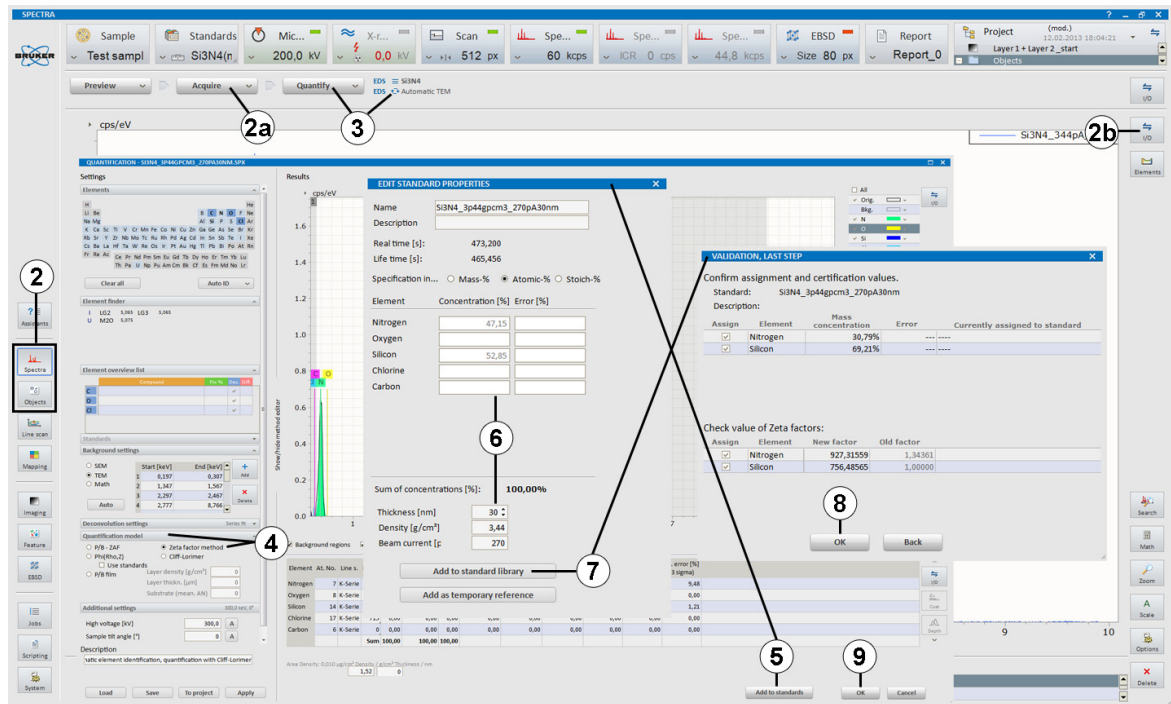



Fig. 5.11-4 Maintain standards library for Zeta factors

### Fit Zeta Factors to Standard

Follow this guide to update Zeta factors for all elements by performing a polynomial fit based on Zeta factors of elements set using standard sample(s).

#### Step

- 1 Set up Zeta factors for the standard sample.
- 2 Click the  icon in the bottom left corner of the **Standards configurator** to open the **ADMINISTRATION OF STANDARD LIBRARY** dialog.
- 3 Click **Zeta-Factors**.

#### Example/hints

Follow section 5.11.3 "Experimental Zeta factors" and see Fig. 5.11-4.

The **EDIT ZETA FACTORS** dialog pops up. The list shows Zeta factors for standard sample in red (default initial values are theoretical Cliff-Lorimer factors). To view the graph, left click into it and zoom by using the arrow keys.

4\* Click on a Zeta-factor value to edit it manually.

\* It is advised to proceed with step 5.

5 Click **Fit to Stds** to perform a polynomial fit based on the already existing standard composition values.

A polynomial is fitted on the standard elements Zeta factors. Zeta factors for elements not contained in the standard sample(s) are interpolated.

6 Press **OK** to confirm the new Zeta factors.

7 Use the updated Zeta factors for quantification.

Follow section 5.13.3.

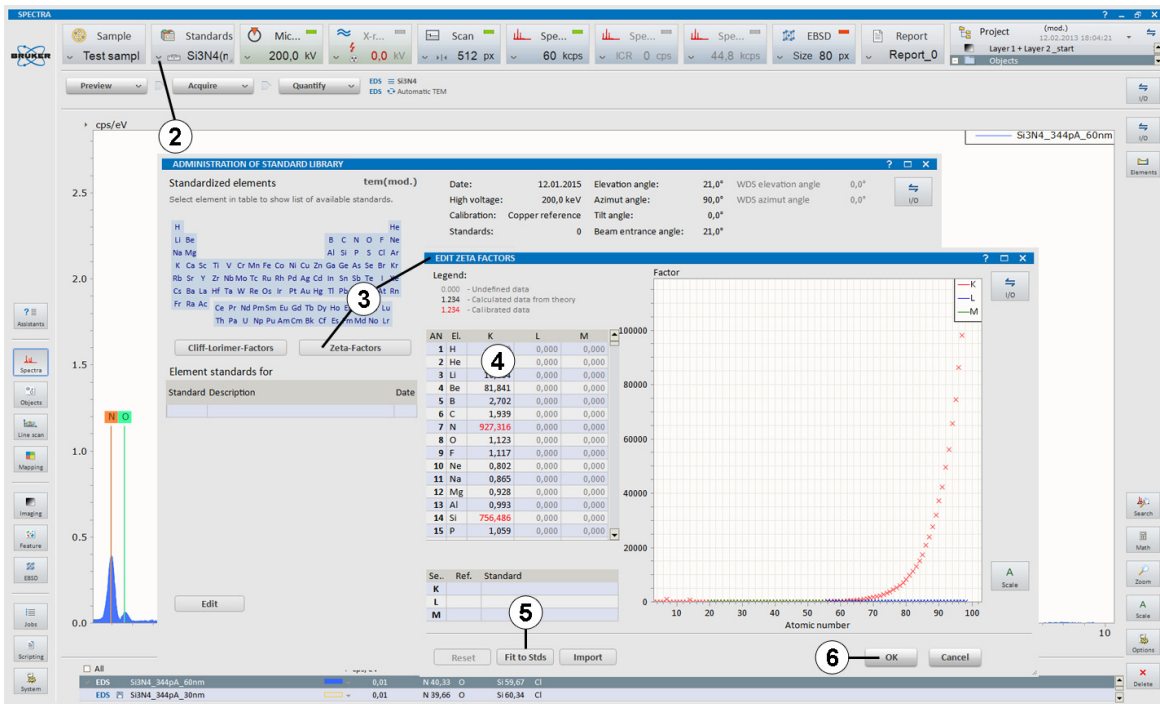







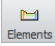
Fig. 5.11-5 Fitting the Zeta factors to standard



## 5.12 Quantitative EDS Analysis of Bulk Samples

### 5.12.1 Using the Method Editor

This section describes how to use the Method editor and how to set up a quantification method for the automatic quantification of multiple spectra. Setting up a quantification method is also important to turn a HyperMap into a QMap (see section 5.18).

<b>Step</b>	<b>Example/hints</b>
<p><b>1</b> Select the <b>Spectra, Objects, Line scan</b> or <b>Mapping</b> workspace.</p>	<p>In the <b>Line scan</b> and <b>Mapping</b> workspaces, spectra are accessible under the <b>Spectrum</b> tab.</p>
<p><b>2</b> Click  on the  button and select <b>Load...</b> to load a quantification method.</p>	<p>Select a quantification method (eg. <b>Default, Oxides, Standards</b> or <b>TEM</b>). Alternatively drag and drop a quantification method from the project to the spectrum chart.</p>
<p><b>3</b> Click  on the  button and</p> <p>a) Select <b>Edit...</b></p> <p>b) Double click on the method name next to the  button.</p>	<p>This opens the <b>METHOD EDITOR</b> to set up a method before performing the quantification.</p>
<p><b>4</b> Use the <b>Elements</b> menu to</p> <p>a) choose how peaks will be labelled</p> <p>b) access element properties</p> <p>c) change settings for the AutoID (when using the <b>Search additional elements</b> option).</p>	<p><b>Use spectrum elements</b> will use element markers that you have previously set in the spectrum via the  icon.</p> <p><b>Use list elements</b> will only use the elements that you have selected (bold) in the periodic table within the method editor.</p> <p><b>Search additional elements</b> will perform an AutoID (see also step 4c).</p> <p>Double click or right click on an element to access the element properties dialog. Use this dialog to set <b>Deconvolution only</b> or to change spectral line for quantification.</p> <p>Set <b>Minimum concentration</b> to an appropriate value to avoid that the AutoID finds too many peaks (especially when spectra are noisy). Elements with a gray background in the periodic table are excluded from the AutoID. To include them click on the specific element.</p>
<p><b>5</b> Use the <b>Element overview list</b> menu to display non-default settings for selected elements.</p>	<p>Use this list to change or review element parameters.</p>
<p><b>6</b> Use <b>Background settings</b> menu to select background fit areas and fit model</p>	<p>The background fit area can be set to <b>Automatic</b> or <b>Preset</b>.</p>

a) Use **Auto** to generate an automatic background fit

b) Activate **+** and drag with right mouse key in the spectrum diagram to set a background fit area manually.

Deactivate the **+** button to use spectrum diagram for Element finder (see step 8)

7 Select the **Deconvolution settings** menu.

Use **Series fit** as default.

8 Select the **Quantification model** menu.

Use **PB/ZAF** as default. Use (Rho,Z) for light elements. See section 5.13.1, 5.13.2, and 5.13.3 for further reference.

9 Use the **Additional settings** menu for

a) Setting high voltage correction

b) Setting sample tilt angle correction

c) Selecting fast quantification.

Click **A** to perform automatic high voltage or sample tilt angle correction.

10 **Save** and rename the modified quantification method for further use.

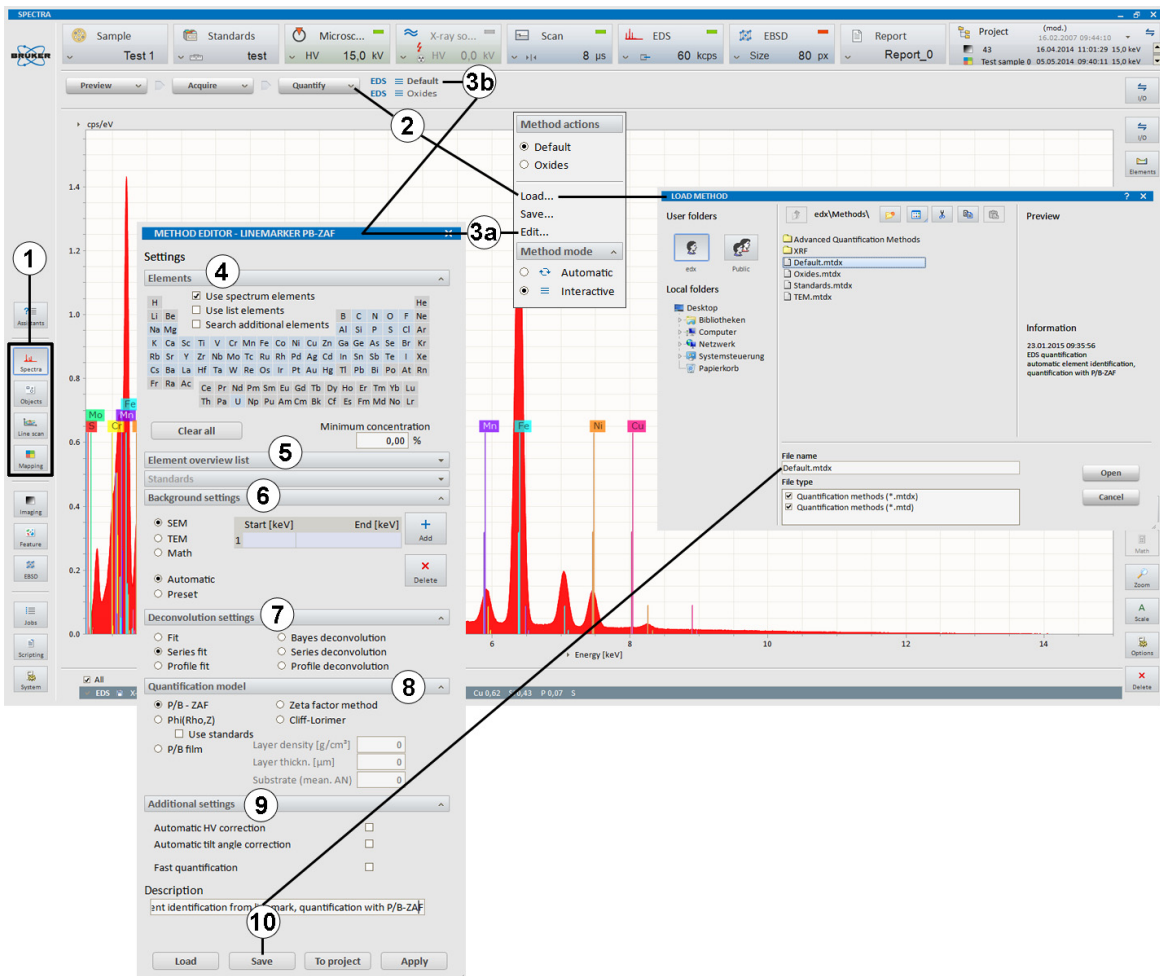
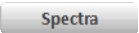


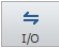








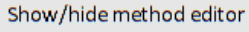






Fig. 5.12-1 Using the method editor

### 5.12.2 Standardless Quantification


This section describes the steps of automatic standardless quantification and quantification using the quantification dialog.

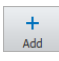
<b>Step</b>	<b>Examples/hints</b>
<p><b>1</b> Select the  or  workspace and</p> <p>a)  a spectrum or</p> <p>b) Use the Spectrum chart  icon to <b>Open</b> a spectrum.</p> <p><b>2</b> Click  on the  button and select <b>Load...</b> to load a quantification method.</p> <p><b>3</b> Click  on the  button and select <b>Method mode</b>.</p> <p><b>4</b> Click  on the  button and select <b>Edit...</b></p> <p><b>5</b> Start quantification by clicking the  button.</p> <p><b>6</b> Edit quantification parameters using the quantification dialog</p> <p>a) Use  to drop down the different settings menus</p> <p>b) Click on  to access/hide the settings of the quantification method.</p> <p><b>7</b> Use the <b>Elements</b> tab to identify elements.</p> <p><b>8</b> The <b>Element finder</b> displays elements with spectral lines at the cursor position in the spectrum diagram.</p>	<p><i>This workflow also applies to any spectrum chart available in the ESPRIT software.</i></p> <p><i>Refer to section 5.7.</i></p> <p><i>Optionally, drag and drop a spectrum from the <b>Project</b> into the spectrum workspace</i></p> <p><i>Select a quantification method (eg. Default, Oxides, Standards or TEM). Alternatively drag and drop a quantification method from the project to the spectrum chart.</i></p> <p> <i>Interactive method</i></p> <p> <i>Automatic method</i></p> <p><i>Use the method editor to set up or edit a quantification method before performing the quantification. This step can be omitted.</i></p> <p><i>If quantification is set to automatic (  ), the quantification results appear in the spectrum list (see section 5.7).</i></p> <p><i>If quantification is set to interactive (  ), the <b>QUANTIFICATION</b> dialog pops up. In this case proceed with step 6.</i></p> <p><i>The left dialog part has the same functionality as the method editor. All changes have an immediate effect on the quantification results displayed in the result table.</i></p> <p><i>Select an element by clicking on the ID in the periodic table. Double click or right click on an element to access the element properties dialog. Use this dialog to set <b>Deconvolution only</b> or to change spectral line for quantification.</i></p> <p><i>Use right click and drag mouse to select a spectral region: all elements are listed in the Finder which have spectral lines in the selected spectrum region.</i></p>

**9** The **Element overview list** displays non-default settings for selected elements.

*Use this list to change or review element parameters.*

**10** Use **Background settings** to set background fit areas and fit model

a) Use  to generate an automatic background fit

b) Activate  and drag with right mouse key in the spectrum diagram to set a background fit area manually.

*Deactivate Add button to use spectrum diagram for Element finder (see point 8)*

**11** Select **Deconvolution settings**.

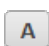
*Use **Series fit** as default.*

**12** Select **Quantification model**.

*Use PB/ZAF as default. Use Phi(Rho,Z) for light elements.*

**13** Use **Additional settings** for

- High voltage correction
- Sample tilt angle correction
- Selecting fast quantification.


*Click  to perform automatic high voltage or sample tilt angle correction, if needed.*

**14** Review spectrum chart.

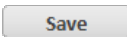
*Compare deconvolution results with measured spectrum.*

**15** Review quantification result table.


**16** Click

a)  to load a quantification method


*Use this to load a previously saved or a default quantification (\*.mtdx) file.*

b)  to save current settings as a quantification method

*Use this to save the current settings as quantification method (\*.mtdx) file.*


c)  to add a method with current settings to project

*Alternatively, use  on the  button.*


d)  to update the active quantification method with the current quantification settings.

*Use this to add current settings to project as a quantification method file.*

*Use this to transfer current quantification settings to method editor.*

**17** a) Use the Spectrum chart  icon to export or save spectrum diagram

*Use this to export deconvolution results or background-subtracted spectra.*

b) Use the Results  icon to **Copy**, **Save** or **Add to Project** or **Add to report**.

**18** Click  to close the **QUANTIFICATION** dialog.

*The results will be displayed in the spectrum list. The display options of the spectrum diagram are described in section 5.8.*

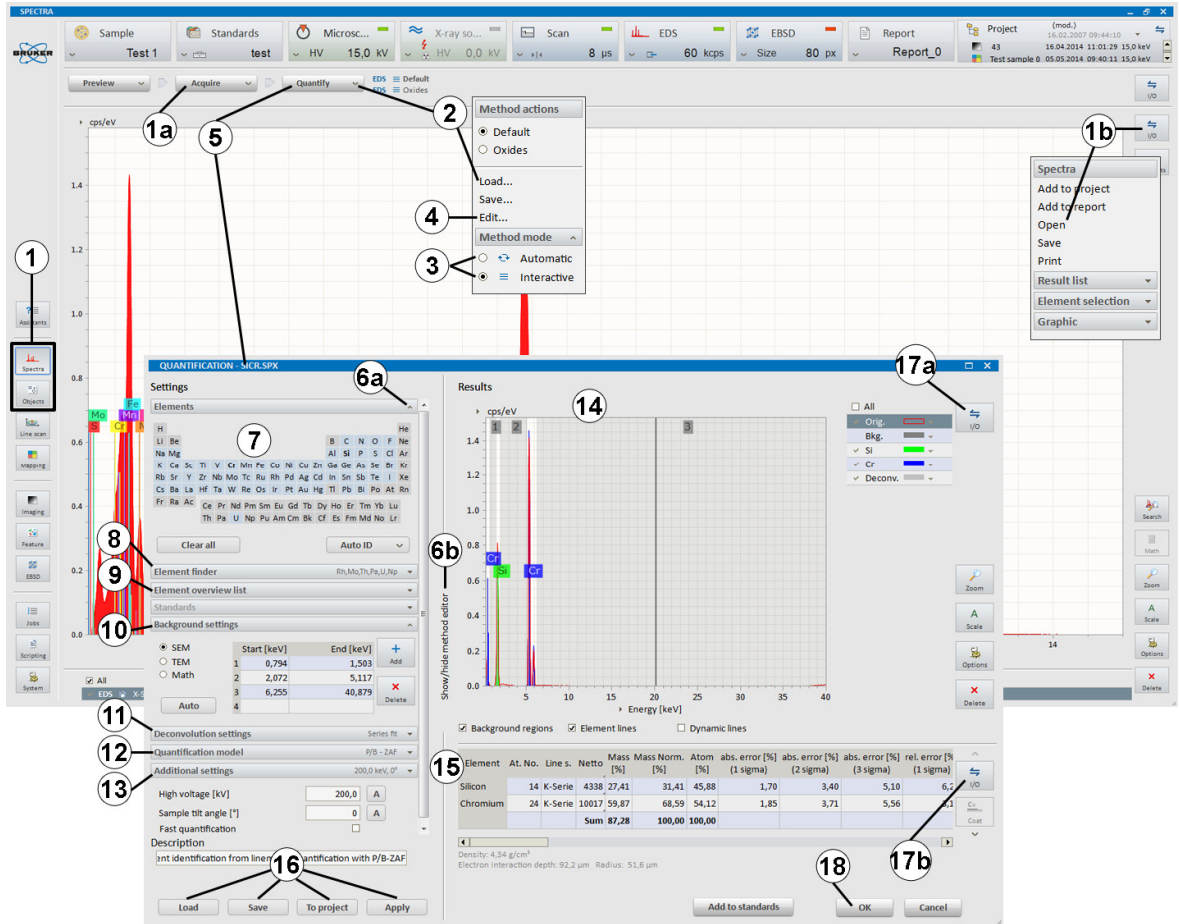
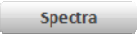






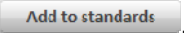
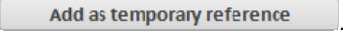
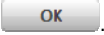
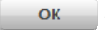
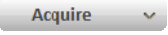


Fig. 5.12-2 Automatic standardless quantification

### 5.12.3 Standard-based Quantification with a Temporary Reference

This guide describes how to perform standard-based quantification without calibration of system factor, measuring reference and unknown samples with same microscope/beam conditions.

<b>Step</b>	<b>Example/hints</b>
<p><b>1</b> Prepare an “unknown” sample, reference sample(s) and a sample for system factor calibration. Select workspace.</p>	<p><i>“Unknown” sample, reference sample(s) and sample for system factor calibration must be available at the same time in the SEM chamber.</i></p>
<p><b>2</b> Calibrate the system factor.</p>	<p><i>For details see section 5.10. Do not change microscope parameters afterwards.</i></p>
<p><b>3</b> Select the  or  workspace.</p>	
<p><b>4</b> Move the reference sample into analytical position and  a spectrum.</p>	<p><i>Follow 5.7 to acquire a spectrum. Use automatic spectrum numbering or rename spectrum after acquisition.</i></p>
<p><b>5</b> Click  at the  button and select <b>Load...</b></p>	<p><i>Load a quantification method.</i></p>
<p><b>6</b> Set <b>Method mode</b> to Interactive ().</p>	
<p><b>7</b> Click .</p>	<p><i>The <b>QUANTIFICATION</b> dialog opens. Set quantification parameters (Element ID, EDS background, Deconvolution) as needed.</i></p>
<p><b>8</b> Click .</p>	<p><i><b>EDIT STANDARD PROPERTIES</b> dialog opens</i></p>
<p><b>9</b> Enter reference concentrations and click .</p>	<p><b>NOTE:</b> <i>The reference values are TEMPORARILY stored until the ESPRIT software is closed. They will not be added into the standards library. For advanced standard-based quantification options see section 5.12.4.</i></p>
<p><b>10</b> The <b>VALIDATION</b> dialog pops up. Close it with .</p>	
<p><b>11</b> Click  to close the <b>QUANTIFICATION</b> dialog.</p>	
<p><b>12</b> Repeat steps 4-11 for other reference samples.</p>	
<p><b>13</b> Move the “unknown” sample into the analytical position.</p>	<p><i>Do not change any relevant microscope settings (beam current, focus, high voltage...)</i></p>
<p><b>14</b>  a spectrum.</p>	<p><i>If automatic quantification during or after acquisition is enabled (see 5.7), the spectrum is quantified automatically and the results appear in the result list.</i></p>

15 Click on the **Quantify** button and select **Load...**

Load a standard-based quantification method e.g. **Linemarker Standards.mtdx**.

Alternatively, open the Method editor and select **Use standards** under **Quantification model**.

16 Click

...unless automatic quantification during or after acquisition is enabled.

17 Use the button of the result table to **Copy**, **Save** or **Add to project** or **Add to report**.

Save the modified quantification method for further use. Use on the button or **Save** in the quantification dialog.

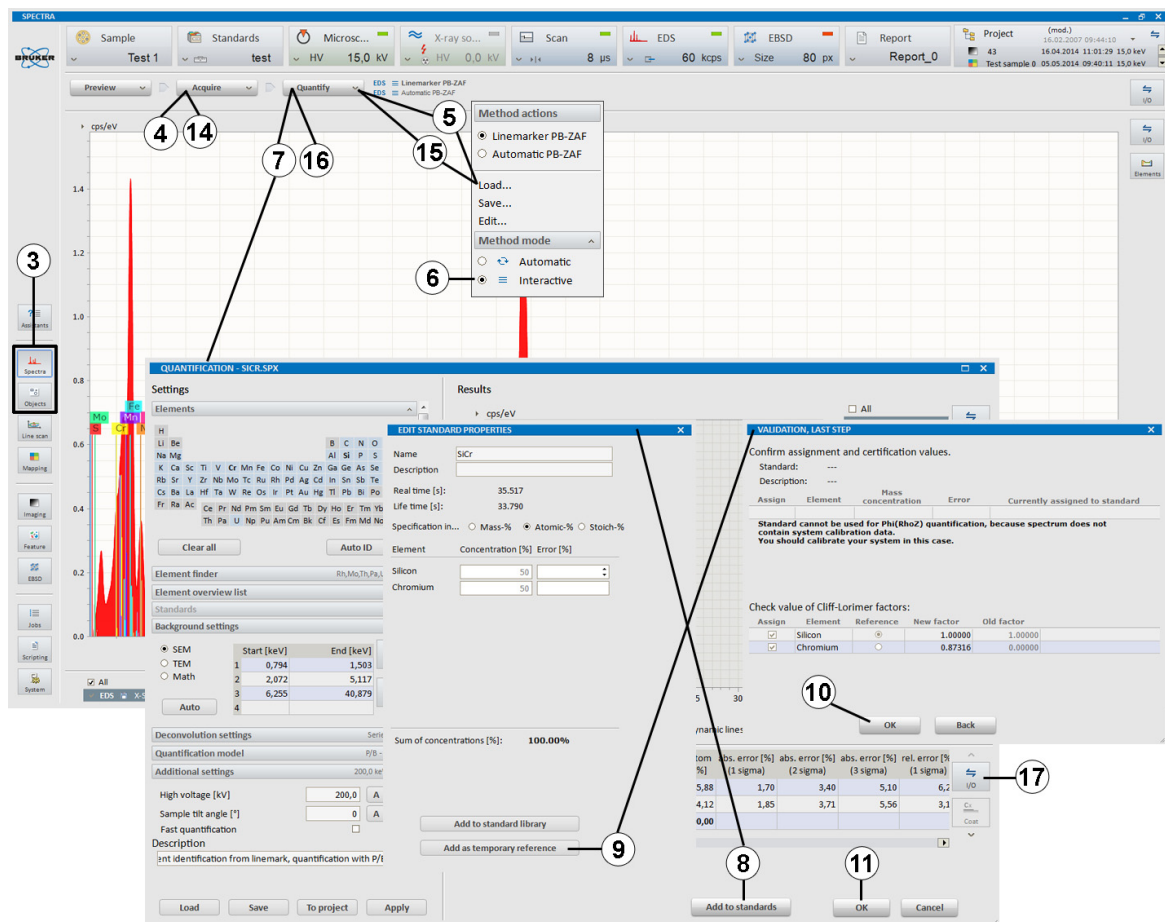

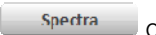

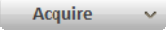
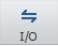



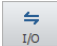


Fig. 5.12-3 Standard-based quantification with temporary reference

### 5.12.4 Standard-based Quantification with Standards Library

Standard-based quantification using a standards library allows the measurement of standard and unknown samples at different times using beam current dependent system factor.

#### Step

- 1 Click the  icon in the bottom left corner of the **Standards configurator** to open the **ADMINISTRATION OF STANDARD LIBRARY** dialog and
  - a) Load a standards library
  - b) Create a New standards library.
- 2 Perform system factor calibration.
- 3 Select the  or  workspace and
  - a)  a spectrum
  - b) Use the Spectrum chart  icon to open a spectrum.
- 4 Click  on the  button and select **Load...** to load a quantification method.
- 5 Click .
- 6 Use the  button of the result table to **Copy, Save** or **Add to Project** or **Add to report**.

#### Example/hints

*Follow section 5.11.1 to measure standard samples and update the standards library.*

*Refer to section 5.10.*

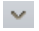

*This workflow also applies to any spectrum chart available in the ESPRIT software.*

*Follow section 5.7.*

*Optionally, drag and drop a spectrum from the **Project** into the spectrum workspace.*

*Select a standard-based quantification method, e.g.: **Linemarker Standard.mtdx**.*

*Follow section 5.12.1. Make sure, that **Use standards** is selected under **Quantification model** in the method editor panel.*

*Save the modified quantification method for further use. Use  on the  button or **Save** in the quantification dialog.*



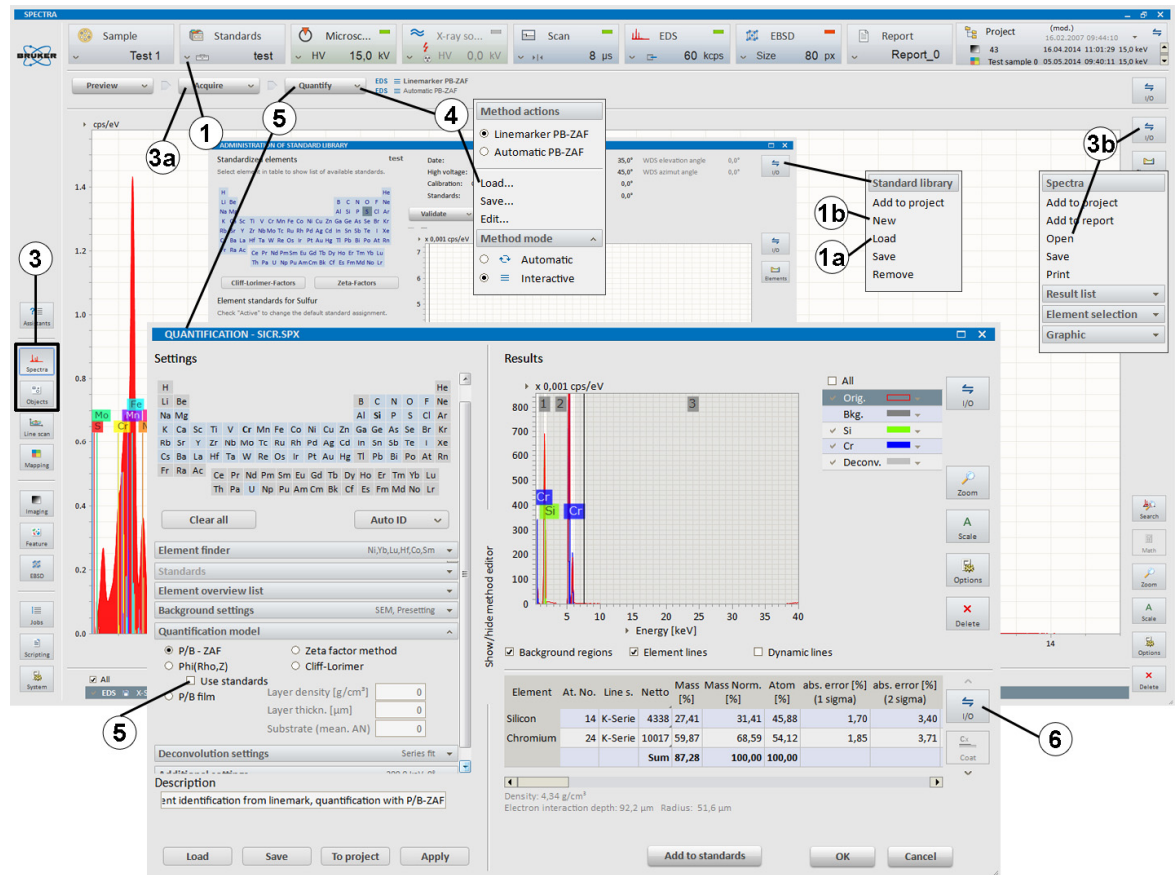
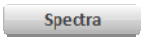



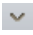





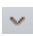







Fig. 5.12-4 Standard-based quantification with a standards library



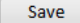
## 5.13 Quantitative EDS of Electron Transparent Samples

### 5.13.1 Standardless Cliff-Lorimer Quantification

This guide describes the steps of standardless quantification of electron transparent samples.

<b>Step</b>	<b>Example/hints</b>
<p><b>1</b> Load or create a standards library containing theoretical Cliff-Lorimer factors.</p>	<p>Follow section 5.11.2.</p>
<p><b>2</b> Select the  or  workspace and</p> <p>a)  a spectrum</p> <p>b) Use the Spectrum chart  icon to open a spectrum.</p>	<p>The following workflow also applies to any spectrum chart available in the ESPRIT software, e.g. also within in the <b>Mapping</b> workspace to quantify complete maps.</p> <p>Follow section 5.7.</p> <p>Optionally, drag and drop a spectrum from the <b>Project</b> into the spectrum chart. The spectrum chart in the <b>Mapping</b> workspace can also be used.</p>
<p><b>3</b> Click  on the  button and select <b>Load...</b> to load a quantification method.</p>	<p>Select a TEM quantification method based on Cliff-Lorimer. Alternatively drag and drop a quantification method from the project to the spectrum chart. Two different quantification methods can be loaded, the active one is displayed in bold.</p>
<p><b>4</b> Make sure that geometrical EDS detector and HV parameters of spectrum and Cliff-Lorimer factor database (standards library) match.</p>	<p>If the information message "<b>Cliff-Lorimer / Zeta factor: Wrong primary energy in standards library</b>" pops up, update the standards library as described in 5.11.2.</p>
<p><b>5</b> Click  on the  button and select <b>Method mode</b>.</p>	<p> Interactive method</p> <p> Automatic method.</p>
<p><b>6*</b> Click  on the  button and select <b>Edit...</b></p>	<p>Use the method editor to set up or edit a quantification method before performing the quantification. For details see section 5.12.1. * step can be omitted</p>
<p><b>7</b> Start quantification by clicking the  button.</p>	<p>If quantification is set to automatic (  ), the quantification results appear in the spectrum list (see section 5.8 step 1f).</p> <p>If quantification is set to interactive (  ), proceed with step 8.</p>
<p><b>8</b> Edit quantification parameters in the quantification dialog.</p>	<p>The left dialog part has the same functionality as the method editor. All changes have an immediate effect on the quantification results displayed in the result table. Refer to section 5.12.1.</p>

9 Use the  icon of the result table to **Copy, Save or Add to Project or Add to report.**

Save the modified quantification method for further use. Use  on the **Quantify**  button or the **Save**  button in the quantification dialog.

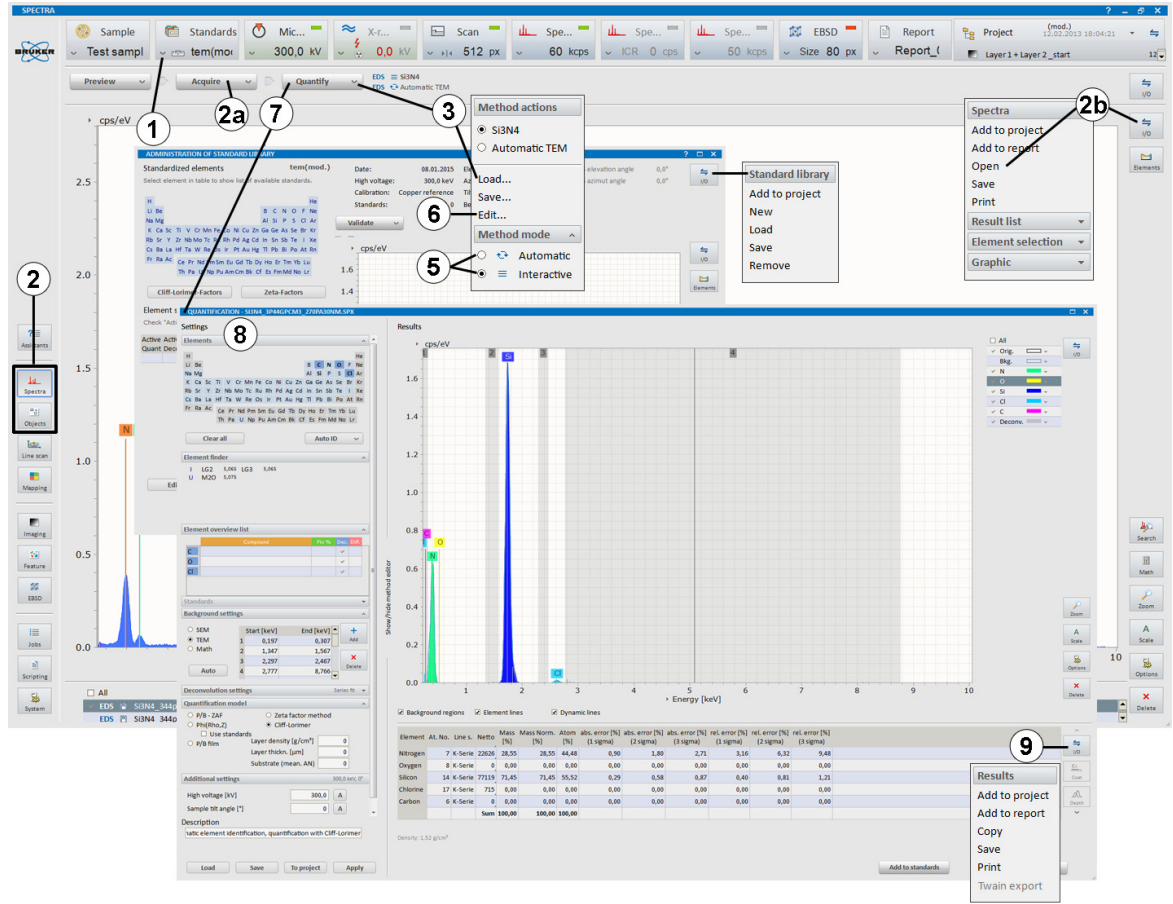
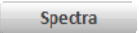
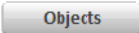
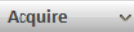




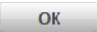
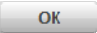


Fig. 5.13-1 Standardless Cliff-Lorimer quantification

### 5.13.2 Standard-based Cliff-Lorimer Quantification


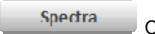

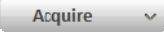




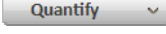
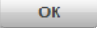
This guide describes the Cliff-Lorimer quantification of electron-transparent samples using standard samples.

<b>Step</b>	<b>Example/hints</b>
<b>1</b> Load or create a standards library containing theoretical Cliff-Lorimer factors.	<i>Follow section 5.11.2.</i>
<b>2</b> Select the  or  workspace and a)  a spectrum b) Use the Spectrum chart  icon to open a spectrum.	<i>Acquire or open a spectrum of the reference/standard sample. The sample should contain elements which are present in the "unknown" sample.</i>
<b>3</b> Follow step 3 to 7 of standardless Cliff-Lorimer quantification to identify elements, set background regions and ensure correct deconvolution.	<i>Refer to section 5.13.1. Set quantification Method mode to <b>Interactive</b> () to access the <b>QUANTIFICATION</b> dialog.</i>
<b>4</b> Click  .	<i>The <b>EDIT STANDARD PROPERTIES</b> dialog pops up.</i>
<b>5</b> Enter reference element concentration values for standard sample.	<i>Use Mass% or Atomic-%.</i>
<b>6</b> Click  .	<i>The <b>VALIDATION</b> dialog pops up with new and old Cliff-Lorimer factors.</i>
<b>7</b> Select an element as reference.	<i>The CL-factor of this element will not be changed. Advice: Keep a higher atomic number value as reference.</i>
<b>8</b> Click  to close the <b>VALIDATION</b> dialog.	<i>The modified Cliff-Lorimer factors will be marked red in the Cliff-Lorimer factor list in the standards library (see section 5.11.2). <b>NOTE:</b> Save the updated standards library.</i>
<b>9</b> Click  to close the <b>QUANTIFICATION</b> dialog.	
<b>10</b> Perform quantification of "unknown" sample.	<i>Follow section 5.13.1. The process is the same as for standardless Cliff-Lorimer factors. The updated standard-based Cliff-Lorimer factors will be used for the quantification.</i>



### 5.13.3 Standard-based Zeta Quantification

This section describes how to perform standard-based quantification of electron transparent samples using beam current and thickness of standard sample as reference values.

Step	Example/hints
<p><b>1</b> Click the <b>Standards configurator's</b> calibration icon  to open the <b>SYSTEM FACTOR CALIBRATION</b> dialog.</p>	
<p><b>2</b> Measure and enter the beam current value as system factor.</p>	<p>Follow section 5.10 for electron transparent samples. Enter the beam current in [pA] units.</p>
<p><b>3</b> Check or Load the appropriate Zeta factors.</p>	<p>Follow section 5.11.3.</p>
<p><b>4</b> Select the  or  workspace and</p> <p>a)  a spectrum</p> <p>b) use the Spectrum chart  icon to open a spectrum</p> <p>c) drag and drop a spectrum from project to Spectrum chart.</p>	<p>Enter the beam current as spectrum parameter: double click on the highlighted line of the active spectrum in the spectrum list (below the spectrum) to open the <b>SPECTRUM PROPERTIES</b> dialog and enter the beam current in pA as system factor in the <b>Parameter</b> tab.</p>
<p><b>5</b> Click  on the  button and <b>Load...</b> a quantification method.</p>	<p>Select a TEM quantification method. Alternatively drag and drop a quantification method from the project to spectrum.</p>
<p><b>6</b> Set the quantification method to <b>interactive</b> ().</p>	
<p><b>7</b> Click on the  button to access the <b>QUANTIFICATION</b> dialog.</p>	
<p><b>8</b> Set <b>Quantification model</b> to Zeta factor method and perform quantification of "unknown" sample.</p>	<p>See section 5.12.1 to optimize quantification and to save or export quantification method and results. Calculated thickness is shown below the result table. Density is calculated based on element composition. If density is known, it can be entered and sample thickness will change correspondingly.</p>
<p><b>9</b> Click  to close the <b>QUANTIFICATION</b> dialog.</p>	

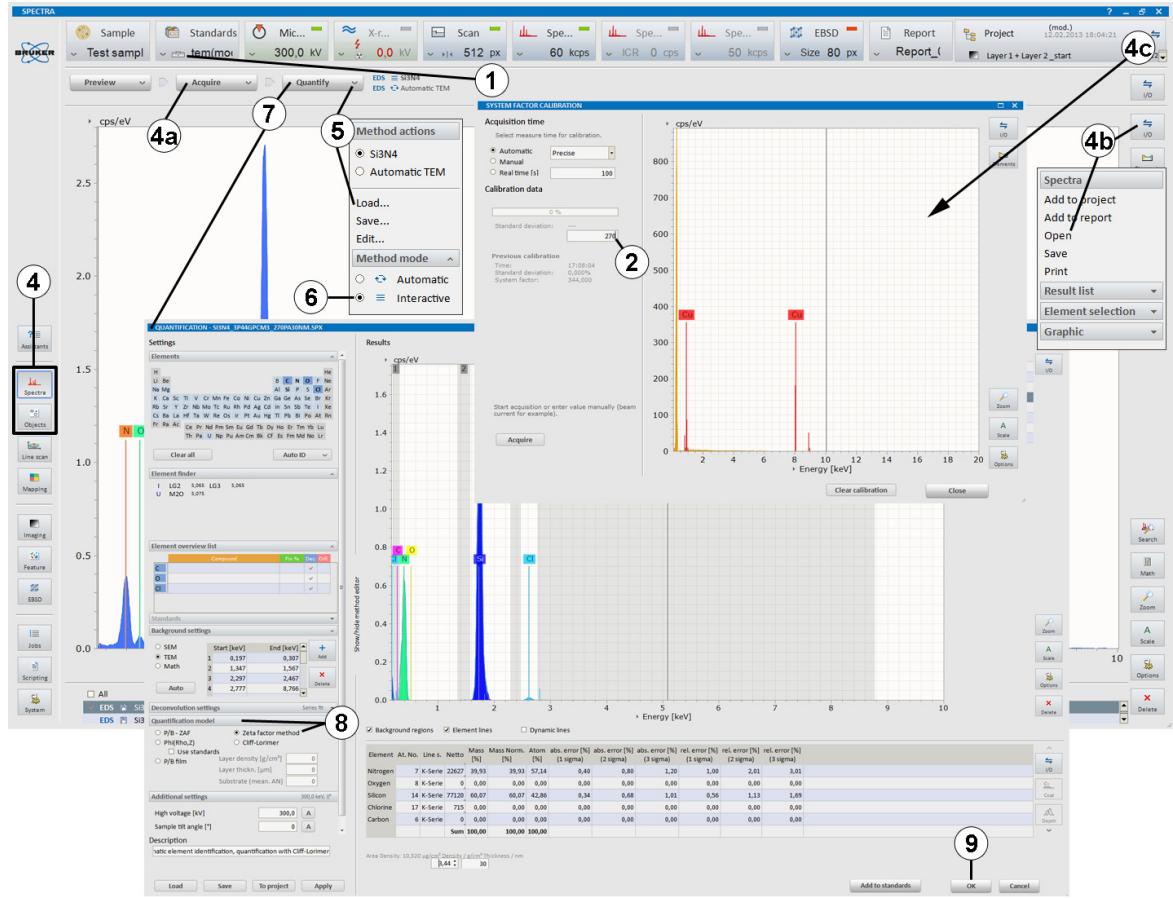

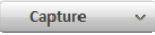




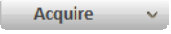





Fig. 5.13-3 Standard-based Zeta quantification

## 5.14 Object Analysis

This section describes the steps of user-defined area analysis with image capture using the Objects workspace.

Step	Example/hints
1 Select the  workspace.	
2  an image.	<i>Refer to section 5.3.</i>
3 Select the desired object type.	<i>Objects types are point, rectangle, ellipse, and polygon or a grid of points (optional grid, random, both).</i>
4 Draw an object in the captured image.	<i>The object can be edited (moved or resized) when the  icon is activated. Click  to delete selected objects.</i>
5 Use  to highlight all objects.	
6 Click  at the  button to set acquisition parameters.	<i>If automatic quantification is desired, check if automatic analysis <b>After acquisition</b> or <b>Continuous</b> is selected.</i>
7  a spectrum.	<i>Objects are automatically acquired in sequence. Progress is indicated by the acquisition bar. If <b>Manual</b> is selected, the acquisition has to be stopped manually.</i>
8 The acquired spectra will appear in the spectrum list.	<i>Refer to section 5.8 for further reference.</i>
9 Use the workspace  icon to a) <b>Save</b> object data	<i>Use .rto file format. This file type contains all available EM images including objects and corresponding spectra.</i>
b) <b>Add data to project</b>	<i>Alternatively, drag and drop data to <b>Project</b> or <b>Report</b>.</i>
c) <b>Add data to report.</b>	
10 Use the spectrum chart  icon to	
a) <b>Save</b> spectrum (or Result list, Element selection, Graphic)	<i>Alternatively, click with the right mouse button into the image or spectrum.</i>
b) <b>Add item to project</b>	
c) <b>Add item to report.</b>	



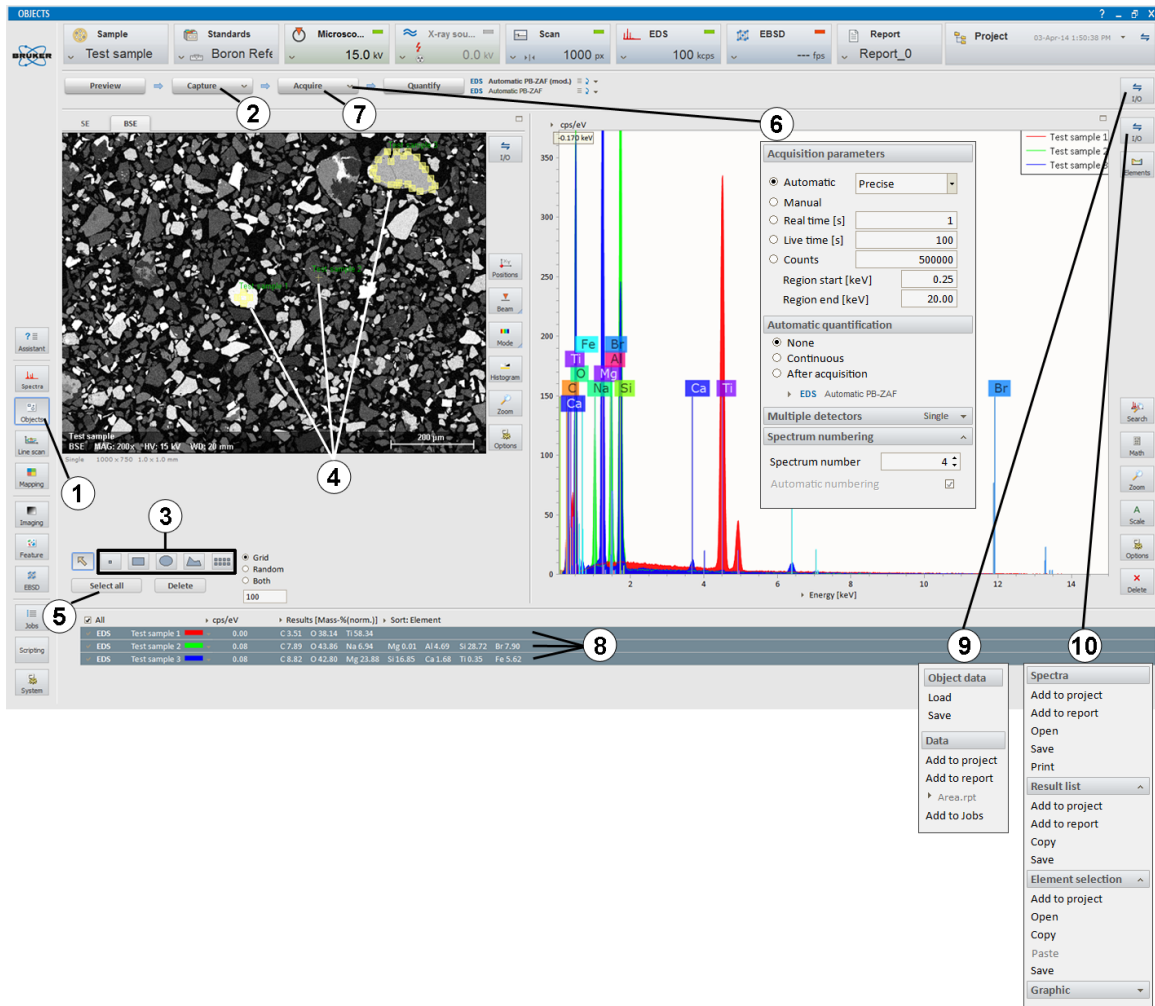
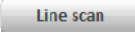
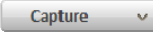

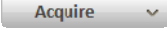



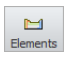
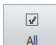
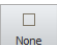
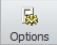

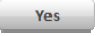


Fig. 5.14-1 Object analysis

## 5.15 Line Scan

This section describes the steps of line scan analysis using the Line scan workspace.

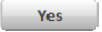
Step	Example/hints
1 Select the  workspace.	
2  an image.	Refer to section 5.3.
3 Highlight the line and drag and adjust the endpoints to the desired position.	
4 Set <b>Point count</b> of the line scan.	Alternatively select <b>Distance [<math>\mu\text{m}</math>]</b> between measurement points.
5 Click  at the  button to set <b>Scan time</b> .	Choose <b>Automatic</b> , <b>Manual</b> or <b>Measurement time [s]</b>
6  a line scan.	The line scan is automatically terminated if <b>Automatic</b> or <b>Measurement time [s]</b> is selected. When <b>Manual</b> is selected, acquisition can be stopped manually by clicking the  button. One click terminates scan after finishing the last scan (the button changes into  , a second click terminates the measurement immediately.
7 Use the  icon to identify elements.	Elements are automatically identified by <b>Auto ID</b> . Add and delete elements by clicking on the <b>Element ID</b> or use the <b>Finder</b> in the <b>Spectrum</b> tab. Identified elements are displayed in the line profile thumbnail bar. The element selection can be changed at any time during or after the line scan acquisition.
8 Select elements in the thumbnail bar by ticking the boxes below the individual element images to display their profiles in the <b>Profiles</b> tab.	Optionally use  or  on the right of the thumbnail bar. Use the  icon to edit the line profile display settings, e.g. scaling of the x-axis, selection of result types, filter strength, etc. (see also Fig. 5.16-1).
9 Use right mouse key in the scan image to extract region of interest spectrum from the line scan.	Extracted spectrum (named as <b>Range</b> ) appears in the <b>Spectrum</b> tab. (optional step). The <b>Scan</b> spectrum in the spectrum tab presents the sum spectrum of the line.
10 Use the workspace  icon to	
a) <b>Save</b> line scan	Use <b>.rtl</b> file format to save line scan data including EM and scan images and point spectra.
b) <b>Add data to project</b>	Image(s), composite element profiles and scan image will be transferred.
c) <b>Add data to report.</b>	Confirm the pop-up window <b>Do you want to save point spectra too?</b> by clicking  to transfer spectral data.


11 Use the Profiles tab  icon to

- a) Save data
- b) Add data to project
- c) Add data to report.

Save the composite element profile as .bmp, .jpg, .png, .tif file format including **Composite element profile, Scan image or both.**

Alternatively, right click into the profile and use the local **Line scan** menu.

Confirm the pop-up window **Do you want to save point spectra too?** by clicking  to transfer spectral data.

12 Use the thumbnail bar  icon to **Add Line scan to report.**

Highlight elements in profile thumbnails. Select **Individual element profiles** to add separate profiles for each element. Select **Composite element profile** to add all selected elements in one diagram and the scan image.

13 To quantify a line scan proceed with section 5.16.

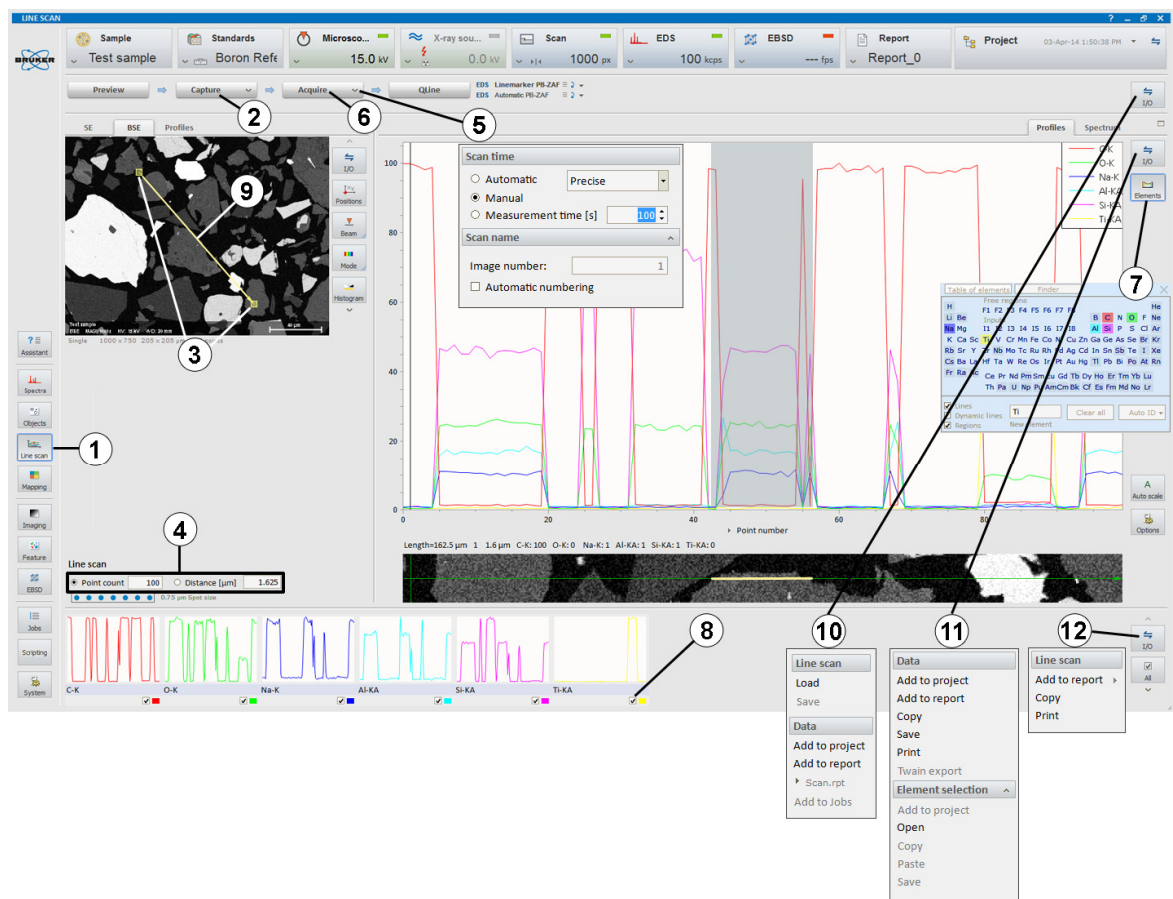
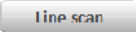




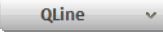
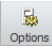

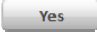

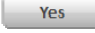
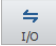


Fig. 5.15-1 Line scan acquisition

## 5.16 Quantitative Line Scan (QLine)

This section describes how to quantify a line scan.

Step	Example/hints
1 Select the  workspace.	
2  a line scan or <b>Load</b> a stored line scan file (.rtl) from the workspace  menu, or retrieve it from the <b>Project</b> folder.	See section 5.15.
3 Click  on the  button and a) <b>Load...</b> an optimized quantification method	Use the Scan spectrum of the Line scan to optimize a quantification method. Save this method to disk or add to project and drag and drop from the project into the Profiles tab. See section 5.12.1 for details. For electron transparent samples use a suitable Cliff-Lorimer or Zeta method.
b) Set <b>Quant range</b> .	To enhance spectra statistics, use values 3, 5, 7, 9, 11, 13 to bin neighboring point spectra.
4 Click  to start the quantification of the line scan database.	
5 Use the  icon in the <b>Profiles</b> tab to change profile display options.	Select from: <b>Counts, Net sum, Mass percent, Mass percent (norm.) and Atomic percent (norm.)</b>
6 Use the workspace  icon to  a) <b>Save</b> the line scan  b) <b>Add data to project</b>  c) <b>Add data to report.</b>	When adding the data into the project, or saving the line scan .rtl file format, the point spectra for each point in the line scan can be added to the <b>Project</b> by confirming the pop-up window <b>Do you want to save point spectra too?</b>
Use .rtl file format to save quantified line scan data including scan image and point spectra.	
Image(s), composite element profiles and scan image will be transferred.	
Confirm the pop-up window <b>Do you want to save point spectra too?</b> by clicking  to transfer spectral data.	
7 Use the <b>Profiles</b> tab  icon to  a) <b>Save</b> data  b) <b>Add data to project</b>  c) <b>Add data to report.</b>	Save as .bmp, .jpg, .png, .tif file format including <b>Composite element profile, Scan image</b> or <b>both</b> .
Alternatively, right click into the profile and use the local <b>Line scan</b> menu. Confirm the pop-up window <b>Do you want to save point spectra too?</b> by clicking  to transfer spectral data.	

- 8 Use the thumbnail bar  icon to **Add** Line scan **to report**.

Highlight elements in profile thumbnails. Select **Individual element profiles** to add separate profiles for each element. Select **Composite element profile** to show all selected elements in one diagram.

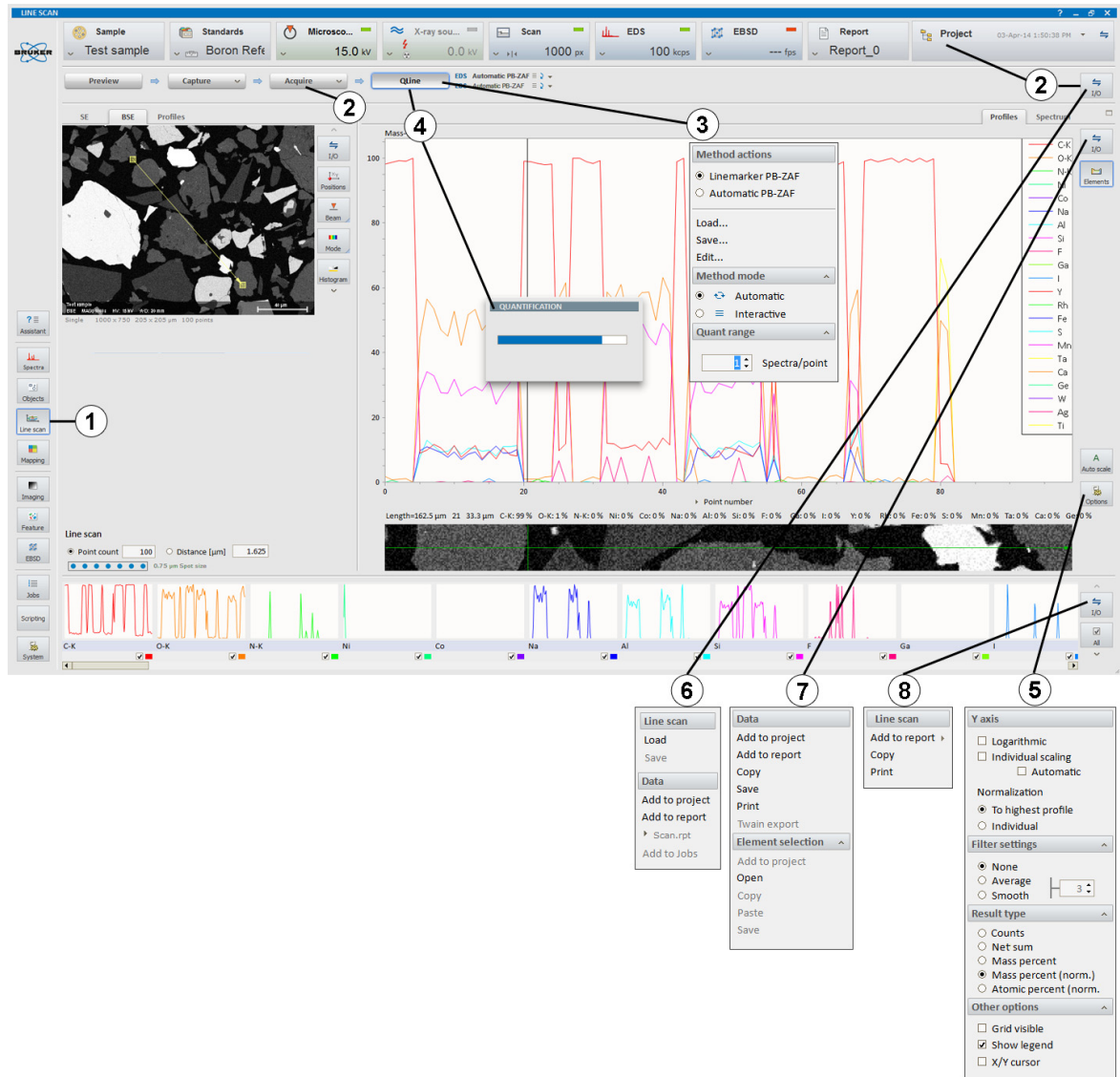
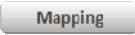

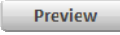
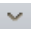

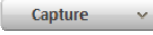

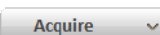
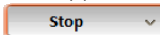



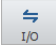


Fig. 5.16-1 Quantitative line scan

## 5.17 Map Acquisition

This section describes the acquisition of maps using the Mapping workspace. Maps can be saved as element distribution images or datacube (HyperMap data)\*.

<b>Step</b>	<b>Example/hints</b>
1 Select the  workspace.	<i>Use high count rate for good count statistics.</i>
2 Click  on the scan configurator to set <b>Image resolution, Mapping dwell time</b> and <b>Line average</b> .	<i>Image resolution defines the pixel resolution.</i>
3 Click  and adjust image.	<i>Adjust brightness, contrast and magnification on the microscope.</i>
4 Click  on the  button to open the <b>Capture parameters</b> menu.	<i>Adjust <b>Capture parameters</b> for the image: set single or continuous (see section 5.3 for further details).</i>
5  an image.	
6 Click  on the  button and set parameters for <b>Map time, After measurement</b> and <b>Map area</b> .	<i>Mapping is automatically terminated if <b>Measurement time [s]</b> or <b>Cycles</b> is set. When <b>Manual</b> is selected acquisition has to be stopped manually by clicking the  button. One click terminates scan after finishing the last frame, a second click terminates the measurement immediately.</i>
7  a map.	<i>The processing of map data is described in section 5.19.</i>
8 Use the workspace  icon to <ol style="list-style-type: none"> <li>a) <b>Save</b> Map data</li> <li>b) <b>Add data to project</b></li> <li>c) <b>Add data to report.</b></li> </ol>	<i>Use file format*:  <b>.bcf</b>: to save HyperMaps (datacube, hyperspectral data set). Spectra for each pixel are saved. Further processing is possible only, when data is saved as a .bcf file.  <b>.rtm</b>: To save element distribution images without point spectra.  <b>.raw</b>: To save hyperspectral datacube for further processing with third party softwares (NIST Lispix format).  <i>Note: Only the EM image, composite element images and Map spectrum will be added to a <b>Project</b> and <b>Report</b>, not the whole hyperspectral database (HyperMap).</i></i>
9 Use the chart  icon of <b>Map</b> tab to <ol style="list-style-type: none"> <li>a) <b>Save</b> map image (Element selection, Settings, Result table)</li> <li>b) <b>Add item to project</b></li> <li>c) <b>Add item to report</b></li> </ol>	<i>Use image file formats (.bmp, .png, .jpg, .tif) to save the composite element image. Alternatively, click with the right mouse button into the map.          Composite element image will be added to <b>Project</b> or <b>Report</b>.</i>

10 Use the thumbnail bar  icon to

a) **Save** Images

*Individual element images of selected thumbnails will be saved. To select them mark the individual element images with left mouse-click + SHIFT or CTRL key.*

b) **Add images to project**

*Composite and individual element images will be added to **Project**.*

c) **Add images to report.**

*Individual element images will be added to **Report**. Alternatively, click with the right mouse button into the thumbnail bar.*

11 Process Map data as desired.

*Refer to section 5.19. For quantitative mapping see section 5.18. For phase analysis see section 5.20.*

\*optional, license-based

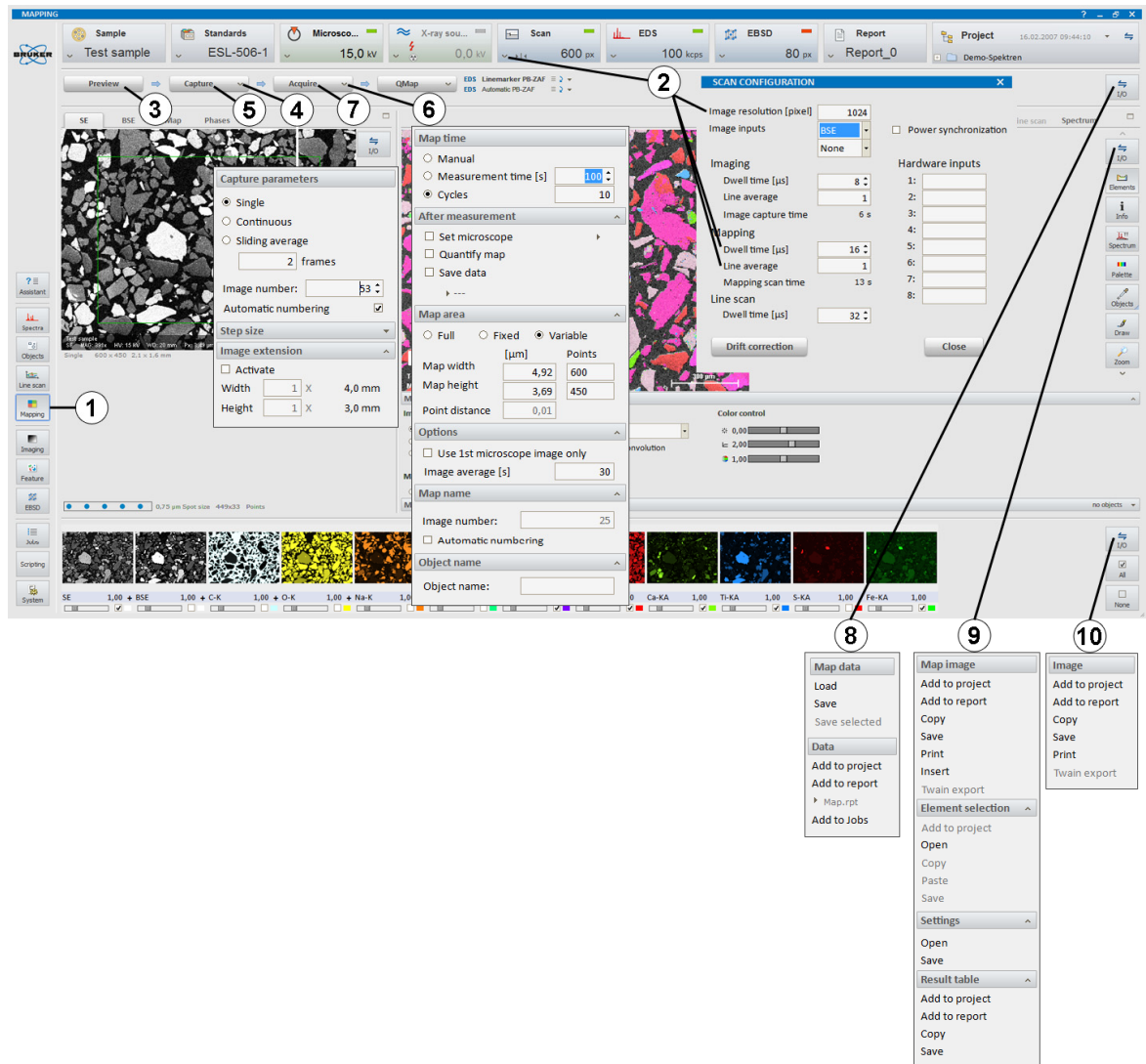
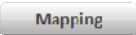
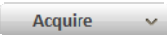

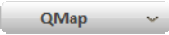
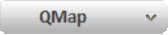
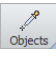


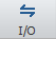


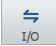
Fig. 5.17-1 Map acquisition

## 5.18 Quantitative Mapping (QMap)

This section describes the quantification of mappings.

Step	Example/hints
1 Select the  workspace.	
2 a)  a map or b) open a .bcf file.	<i>Good spectrum statistics may improve the quantification accuracy.</i>
3 Click  on the  button and <b>Load...</b> an appropriate quantification method.	<i>Use a previously optimized method (presetting element IDs, background fit regions and quantification algorithm).</i>
4 Set <b>QMap options</b> .	<i>Use tile size of 3x3, 5x5 or 9x9 pixels to reduce processing time and enhance count statistics. The <b>Estimated time</b> to complete the QMap will be displayed.</i>
5 Click  to start the quantification of the map.	<i><b>QMap</b> will terminate automatically after the last pixel has been quantified.</i>
6 Adjust colors and image controls.	<i>See <b>Map Display settings</b> and refer to section 5.19.</i>
7 Use the  tool to extract quantification results of regions of interest.	<i>Quantification results appear in the Map result list.</i>
8 Click on  to switch to pseudo colors display.	<i>The pseudo colors display presents numerical quantification values. Choose one element only for pseudo color display. If several elements are selected, the map displays summarized concentration values.</i>
9 Set <b>Result type</b> in <b>Map display settings</b> submenu.	<i>Select Net sum, Mass %, normalized Mass % or Atomic %. Select <b>Counts</b> to display non-quantified intensity maps.</i>
10 Use the workspace  icon to a) <b>Save</b> Map data b) <b>Add data to project</b> c) <b>Add data to report.</b>	<i>Save the quantified datacube (Hypermap) as .bcf file. Alternatively, drag and drop composite element image to project or right click into the map and use the local Mapping menu.</i>
11 Use the chart  icon in <b>Map</b> tab to a) <b>Save</b> Map image (Element selection, Settings, Result table) b) <b>Add item to project</b> c) <b>Add item to report.</b>	<i>Alternatively, click with the right mouse button into the map image.</i>



12 Use the thumbnail bar  icon to

a) **Save Images**

*Individual element images of selected thumbnails will be saved.*

b) **Add images to project**

*Composite and individual element images will be added to the **Project**.*

c) **Add images to report.**

*Individual element images will be added to the **Report**. Alternatively, click with the right mouse button into the thumbnail bar.*

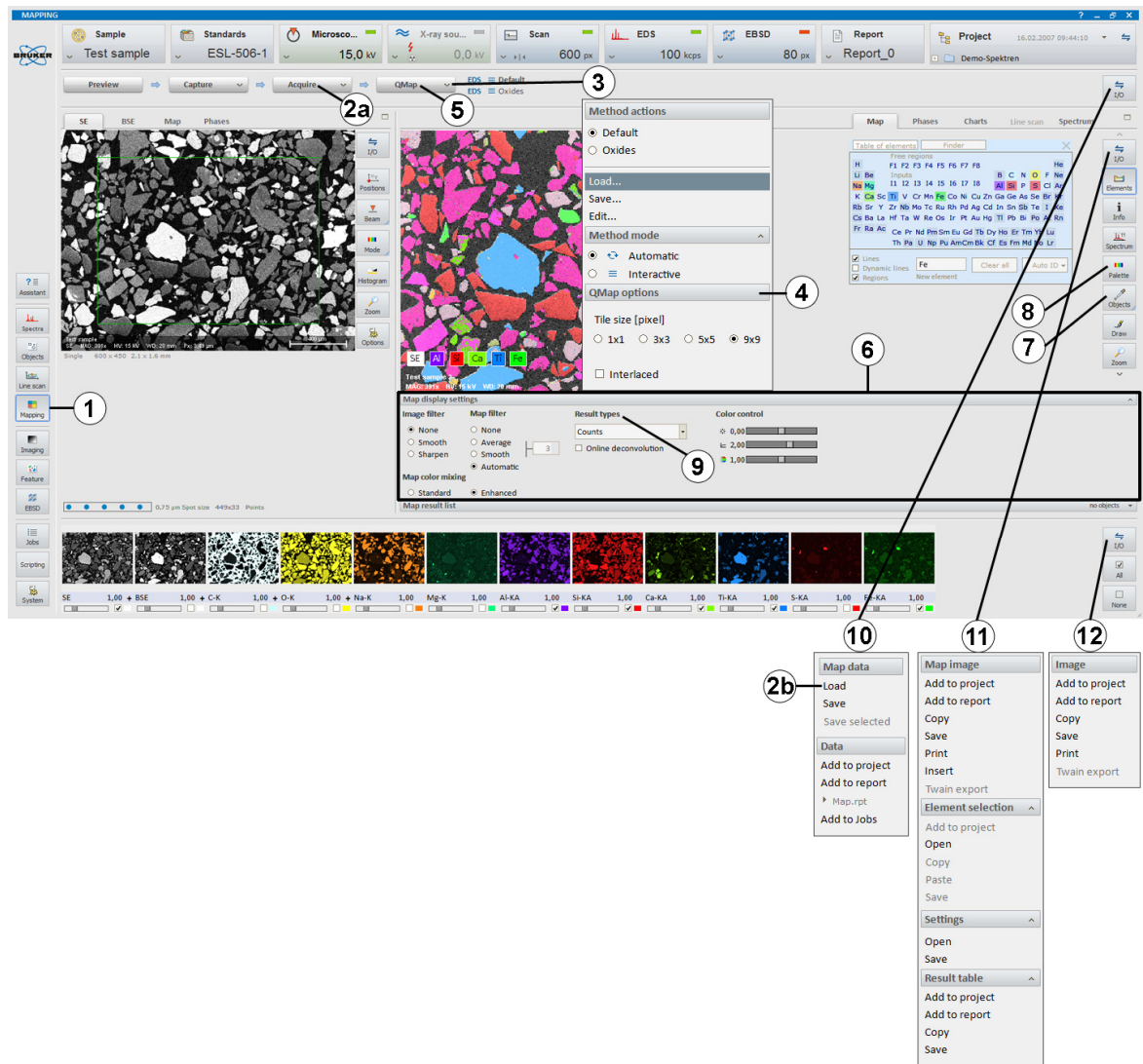
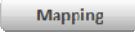


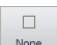
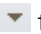
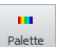
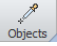
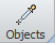



Fig. 5.18-1 Quantitative mapping

## 5.19 Map Processing

This section describes the online and offline processing of HyperMap data (datacube/.bcf files\*).

Step	Example/hints
<p>1 Select the  workspace.</p>	
<p>2 Use the  icon to open the periodic table and identify elements.</p>	<p>Select elements by clicking on the <b>Auto ID</b> or use the <b>Finder</b> in the <b>Spectrum</b> tab. Identified elements are displayed in the map thumbnail bar.</p>
<p>3 Select/Deselect element(s) in the thumbnail bar by ticking/unticking the boxes below the individual element images to display their distribution in the large/ mixed map image.</p>	<p>Optionally use  or  on the right of the thumbnail bar.</p>
<p>4 Click  to open the <b>Map display settings</b> submenu and adjust</p> <p>a) <b>Image filter</b></p> <p>b) <b>Map filter</b></p> <p>c) <b>Result types</b></p> <p>d) <b>Color control</b></p> <p>e) <b>Map color mixing.</b></p>	<p>Alternatively, click with the left mouse button into the <b>Map display</b> bar.</p> <p>Applies filters on the overlay EM image.</p> <p>Applies filters on the element distribution map.</p> <p>Use <b>Online deconvolution</b> to display deconvolved elements distribution maps.</p> <p>Use sliders to apply image brightness, gamma correction and color saturation.</p> <p>Use <b>Enhanced</b> to improve the display of EM image and element map overlay.</p>
<p>5 Click on  to switch to pseudo colors display.</p>	<p>Choose one element only for pseudo color display. Right click on the icon to adjust <b>Palette settings</b>. Note: If the map is not quantified, the values are normalized to the highest displayed intensity value.</p>
<p>6 Extract point/area spectra from the map:</p> <p>i. Right click on  to access object tool list</p> <p>ii. Select the preferred tool</p> <p>iii. Click on  to activate the object tool</p> <p>iv. Move the cursor to the area of interest, press left mouse key and drag cursor to expand object.</p>	<p>View and quantify the extracted point/area spectra in the <b>Spectrum</b> tab.</p> <p>Select point, line, rectangle, ellipse or polygon.</p> <p>Object spectra appear in <b>Spectrum</b> tab. Line profile appears in the <b>Line scan</b> tab. Enhance line profile statistics: use + and – keys to set line width. Line scan can be transferred to the Line Scan workspace via the Project. To quantify a map or a line scan, see section 5.18 or 5.16.</p>
<p>7 Click on  to generate a MaximumPixelSpectrum</p>	<p>Identification of locally enriched elements with low total concentration. The MaximumPixelSpectrum is displayed in the <b>Spectrum</b> tab.</p>


- 8 Select the **Chart** tab to see element correlations or element histogram.

- 9 Use the workspace  icon to

a) **Save** Map data


b) **Add data to project**

c) **Add data to report.**

- 10 a) Use the chart  icon in the **Map** tab to **Save** the map image (Element selection, Settings, Result table)

b) Use the chart  icon in the **Charts** tab to **Save** the charts image

c) Use the chart  icon in the **Spectrum** tab to **Save** the spectrum.

- 11 Use the thumbnail bar  icon to

a) **Save** images

b) **Add images to project**

c) **Add images to report.**

\*optional, licence-based

Choose **Color settings**, **Mode** and displayed **Elements** in the **Chart settings** below the Chart diagram. Make sure to have good count statistics in the map.

Use file format\*:

**.bcf**: to save HyperMaps (datacube, hyperspectral data set). Spectra for each pixel are saved. Later offline (re-) processing is only possible when data is saved as a .bcf file.

**.rtm**: To save element distribution images without point spectra.

**.raw**: To save hyperspectral datacube for further processing with third party softwares (NIST Lispix format).

Note: Only the EM image, composite element images and Map spectrum will be added to a **Project** and **Report**, not the whole hyperspectral database (HyperMap).

Use image file formats (.bmp, .png, .jpg, .tif) to save composite element image. Alternatively, click with the right mouse button into the map. The options **Add to project** and **Add to report** are also available.

The options **Add to project** and **Add to report** are also available.

Alternatively, click with the right mouse button into the spectrum. The options **Add to project** and **Add to report** are also available.

Individual element images of selected thumbnails will be saved. To select them mark the individual element images with left mouse-click + SHIFT or CTRL key.

Composite and individual element images will be added to **Project**.

Individual element images will be added to **Report**. Alternatively, click with the right mouse button into the thumbnail bar.

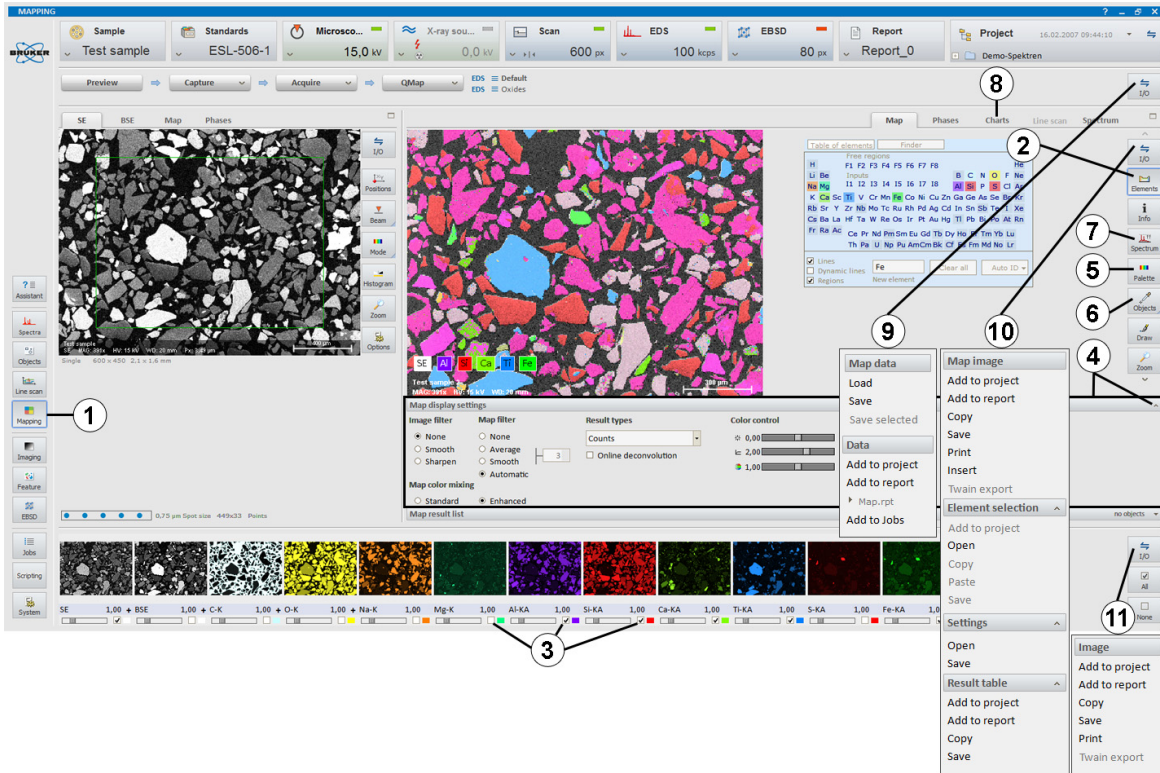
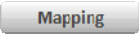

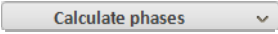
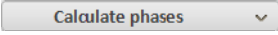

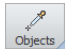


Fig. 5.19-1 Map processing

## 5.20 Phase Analysis

The AutoPhase tool calculates images that reveal the local distribution of different phases in a sample.

<b>Step</b>	<b>Example/hints</b>
<p><b>1</b> Select the  workspace and acquire a map or open a bcf. or rtm. file.</p> <p><b>2</b> Select the <b>Map</b> tab and the elements of interest for phase calculation in the thumbnail bar.</p> <p><b>3</b> Select the <b>Phases</b> tab.</p> <p><b>4</b> Use  on the  button to open the corresponding menu and activate <b>Automatic update</b> option.</p> <p><b>5</b> Click .</p> <p><b>6</b> Use  on the <b>AutoPhase settings</b> sub-menu to open it.</p> <p><b>7</b> Select the AutoPhase <b>Method</b></p> <p style="margin-left: 20px;"><b>i. PCA</b></p> <p style="margin-left: 20px;"><b>ii. Clusters</b></p> <p style="margin-left: 20px;"><b>iii. Objects.</b></p> <p><b>8</b> Select and adjust the <b>Options</b> for the phase calculation</p>	<p><i>Refer to section 5.17. Phase analyses can be performed on qualitative and quantitative maps.</i></p> <p><i>Select elements by ticking the boxes below the individual element images in the thumbnail bar. Only selected elements will be used for phase calculation.</i></p> <p><i><b>Automatic update</b> is useful when experimenting with methods and parameters. The phase image will update after every change.</i></p> <p><i>Phases will be calculated and appear in the thumbnail bar.</i></p> <p><i>Set parameters for phase determination.</i></p> <p><i>All three AutoPhase methods work based on the number of element maps chosen. Neighboring pixels containing the same chemical information represent a particular phase. Depending on the chosen sensitivity, pixels with similar composition get merged into one phase.</i></p> <p><i><b>PCA</b> (Principal Component Analysis) is a specific type of vector analysis reducing the dataset to the most relevant chemical information. It starts its operation with an n-dimensional "element" space. The intensity of every element selected in the map thumbnail bar (below the left-hand side <b>Map</b> tab) spans one axis of this space. Map pixels that contain the same chemical information form data clouds which are separated by the PCA dimension reduction algorithm.</i></p> <p><i>The <b>Cluster</b> method compares the composition of neighboring pixels (by running through the map pixel by pixel).</i></p> <p><i>The phases are user-defined by manually defining objects using drawing tools accessible through the  icon.</i></p> <p><i>Use the sliders to change the values.</i></p>

i. **Sensitivity**

Represents how similar (high sensitivity) or different (low sensitivity) the chemical compositions of the distinguished phases area.

ii. **Area**

Represents the minimum size of a phase, ranging from 0 to 1 area % of the map image.

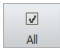
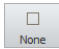
iii. **Edges.**

Represents the sensitivity used to connect small phases.

**9** Phase results will appear in the **Phase result list**.

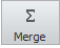
Choose area %, area pixels and area  $\mu\text{m}/\text{mm}$ . You can rename P1, P2, ... by selecting a phase from the list and clicking on its name. Rename and press <ENTER>.


**10** Select or deselect the phases in the thumbnail bar.

Optionally use  or  on the right of the thumbnail bar.

**11** Click on a phase thumbnail to highlight it and

a) Merge phases using the  icon

Several phases can be merged into a single phase by highlighting the particular phases in the clipboard (press CTRL or SHIFT key to highlight multiple thumbnails) and pressing the  icon.

b) Extract a spectrum of a highlighted phase using the  icon.

The spectrum will be displayed in the **Spectrum** tab.

**12** Use the workspace  icon to

a) **Save** Map data

Save datacube (Hypermap) as .bcf file or map data as .rtm file.

b) **Add** data **to project**

Alternatively, drag and drop data to the **Project** or **Report**.

c) **Add** data **to report**.

**13** Use the **Phases** tab  icon to

a) **Save** Phase data

Alternatively, right click into composite image and use the **Phases** menu.

b) **Add** Phase data **to project**

c) **Add** Phase data **to report**.

**14** Use the thumbnail bar  icon to

a) **Save** Phase images

Individual phase images of selected thumbnails will be saved.

b) **Add** Phase images **to project**

Composite and individual phase images will be added to **Project**.

c) **Add** Phase images **to report**.

Individual phase images will be added to **Report**.

The screenshot displays the Quantax EDS software interface for phase analysis. The main window shows a color-coded phase map of a sample. The 'Phases' panel on the right contains 'AutoPhase options' and 'AutoPhase settings'. The 'Phase result list' table provides quantitative data for four phases (P1, P2, P3, P4). A context menu is open over the map, showing options for 'Map data', 'Phase data', and 'Image'.

Counts/Pixel	Area	BSE	C-K	O-K	Na-K	Mg-K	Al-KA	Si-KA	Ca-KA	Ti-KA	S-KA	Fe-KA
P1	46.4 %	3798	8337.48	31.80	57.49	69.15	76.49	67.00	35.53	26.68	57.72	17.78
P2	21.7 %	22107	19693	113.97	2023.40	822.23	79.45	1372.08	3597.66	51.53	37.75	55.45
P3	10.9 %	17662	14406	111.74	1952.20	59.82	78.25	116.63	5747.47	44.67	35.76	52.29
P4	8.4 %	39026	50063	160.13	1296.72	91.90	91.54	125.53	2330.24	96.94	1339.26	108.33

Fig. 5.20-1 Performing phase analysis

## 5.21 Creating a Report

This section describes how to create a report based on the data obtained or processed in all workspaces of QUANTAX ESPRIT.

<b>Step</b>	<b>Example/hints</b>
<b>1</b> Select a workspace and perform an analysis or process data.	<i>Reports can also be composed from files or project entries.</i>
<b>Add data to report:</b>	
<b>2a</b> Use the workspace  icon and select <b>Add to report</b>	<i>All results contained in the workspace will be added to the report.</i>
<b>2b</b> Use the chart  icon and select <b>Add to report</b>	<i>Alternatively, right click into the chart and use the local menu.</i>
<b>2c</b> Use the thumbnail bar  icon and select <b>Add to report</b> .	<i>Available in the <b>Mapping</b> or <b>Line scan</b> workspace. Alternatively, a right mouse click into the thumbnail bar opens the corresponding local menu.</i>
<b>3</b> Use the  icon in the bottom left corner of the report configurator to open the <b>Report</b> .	<i>The <b>Report preview/Report</b> is either docked to the right side of the screen or to the configurator bar. The display can be changed using the  icon in the <b>Project</b> tool.</i>
<b>4</b> Edit/ review the report	
a) Use <b>Zoom</b> to set display	
b) Use <b>Properties</b> to set object layer order	<i>Objects are displayed as layers. Uncheck <b>Fixed</b> when moving an object. Check <b>Pass on</b> to show object on every page.</i>
c) Use <b>Objects</b> to add text	
d) Use <b>Tools</b> for drawing.	<i>Draw <b>Point, Arrow, Line, Rectangle, Circle/Ellipse</b>. Set fill/color/size options for selected element.</i>
e) Use the  icon in the <b>Report</b> editor to import data.	<i>Open image, spectrum or text.</i>
<b>5</b> Rearrange order of items.	<i>Drag objects (maps, images, spectra...) and move them to the desired position.</i>
<b>6</b> Use the <b>Report</b>  icon to <b>Save as</b> report.	<i>Alternatively, use the <b>Report Preview</b>  icon. The report is saved as .rpt.</i>
<b>7</b> Use the <b>Report</b>  icon to <b>Export to WORD</b> .	<i>If Microsoft Word® is available.</i>
<b>8</b> Click the  button to print the report.	<i>Check page <b>Setup</b> before printing. Use printer or generate a .pdf file.</i>



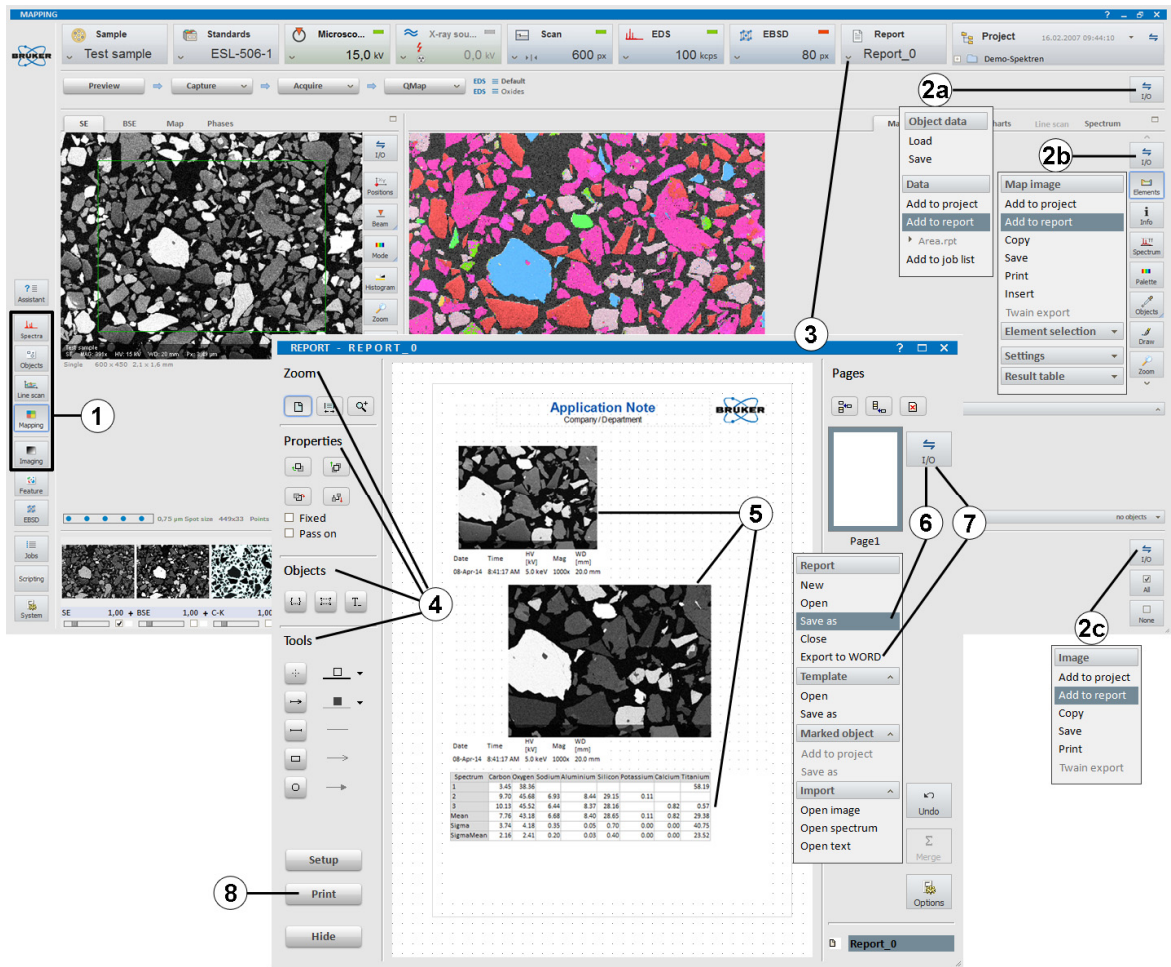


Fig. 5.21-1 Creating a report

## 5.22 Managing Projects

This section describes how to use Project to manage and archive data obtained or processed by ESPRIT. Use Project as a data clipboard to transfer data between the workspaces or save all data added to Project as one file to disk. Add images, spectra, element selection, quantification methods, result tables, map/phase images, line scan data (including point spectra).

Step	Example/hints
<p><b>1</b> Use the <b>Project</b>  icon to</p> <p>a) Generate a <b>New</b> Project</p> <p>b) <b>Open...</b>a Project file from disk.</p> <p><b>2</b> Select a workspace and perform analysis or process data.</p>	<p><i>An empty project list is generated.</i></p>
<b>Add data to project:</b>	
<p><b>3a</b> Use the workspace  icon and select <b>Add to project</b></p>	<p><i>All results contained in the workspace will be added to the project.</i></p>
<p><b>3b</b> Use the chart  icons and select <b>Add to project</b></p>	<p><i>Alternatively, right click into the image chart or result chart and use the local menu.</i></p>
<p><b>3c</b> Use the thumbnail bar  icon and select <b>Add to project</b></p>	<p><i>Available in the <b>Mapping</b> workspace.</i></p>
<p><b>3d</b> Drag and drop an item from the current workspace/chart to the <b>Project</b></p>	<p><i>The item appears in the <b>Project</b> list.</i></p>
<p><b>3e</b> Use the <b>Project</b>  icon to <b>Import</b> an item from the hard disk.</p>	<p><i>Use the  icon to toggle the position of <b>Project</b>. It is either docked to the right side of the screen or to the configurator bar. Double click an item in the project to review.</i></p>
<b>Retrieve an item from Project to a workspace:</b>	
<p><b>4</b> Drag and drop an item from <b>Project</b> to the current workspace/chart.</p>	<p><i>Use this function as a data clipboard for data transfer between workspaces of ESPRIT.</i></p>
<b>Save project to hard disk:</b>	
<p><b>5</b> Use the <b>Project</b>  icon to <b>Save</b> the project to hard disk.</p>	<p><i>The project is saved as .rtx file. Use <b>Save</b> to overwrite the current project and use <b>Save as...</b> to add a new name.</i></p> <p><i>Note: (mod.) is displayed after the project name if an opened project file has been modified and the current changes are not saved.</i></p>

Please note:

Items in the project are not automatically saved to the hard disk! Save the project file manually to archive your data.

Do not use Project to store HyperMap data. The current map display can be added to the project, but not the HyperMap data (point spectra). HyperMaps should be saved as described in 5.17.

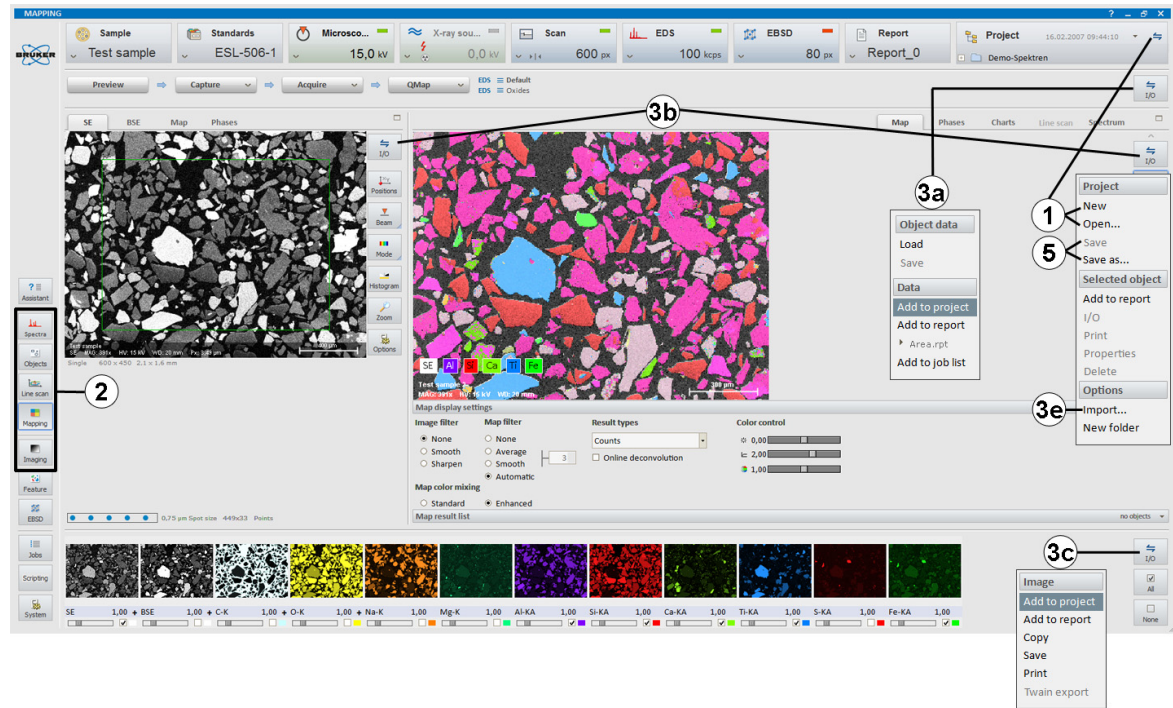


Fig. 5.22-1 Creating a project



# Index

## A

Access rights .....	21
Administrator rights .....	18
Analysis	
Line scan .....	98
Map .....	106
Objects .....	96
Assistants .....	28, 47
Automation .....	29
AutoPhase .....	109

## B

Backing up data .....	21
Beam current setting .....	71
Beryllium	
Safety issues .....	10
Blind flange, original .....	9
Broken window .....	20
BSE detector .....	30

## C

Calibration	
Energy-Channel .....	53
Spectrometer .....	44, 53
Capturing images .....	55
Client-server architecture .....	17, 25
Cliff-Lorimer factors	
Experimental .....	76
Theoretical .....	76
Coating correction .....	68
Coating sample .....	32
Command field .....	42
Communication active, control .....	33
Communication server .....	26
Compatibility, electromagnetic .....	11
Configurator	
Microscope .....	33
Project .....	38
Report .....	38
Sample .....	32
Scan .....	34
Spectrometer .....	36
Cooling mode .....	20, 36
Copy to clipboard .....	39

## D

Data	
Backing up .....	21
File formats .....	39
Load file .....	39
Loading .....	39
Save file .....	39
Saving .....	39

Transfer of SEM- .....	17
Datcube .....	102, 106
Dead time .....	16
Delete	
User .....	27
Detector	
Broken window .....	10, 20
Window .....	20
Device settings .....	42
Display settings .....	43
Drift correction .....	59
Dwell time .....	34

## E

EBSD .....	28
Electromagnetic interference .....	11
Electron beam	
Deflection .....	17
Electron transparent samples	
Cliff-Lorimer quantification .....	90, 92
Zeta quantification .....	94
Electronic noise .....	16
Electrostatic damage .....	20
Element distribution images .....	102, 106
Element identification	
Exchanging data .....	39
Saving data .....	39
Element selection	
Storing .....	39
Energy	
Range .....	36
Resolution .....	16
Energy dispersive spectrometry .....	15
Energy-Channel calibration .....	53
Error recovery .....	21
Export	
Graphic .....	39
Extension, Image .....	57

## F

Feature .....	28
File	
Data .....	39
Format .....	39
License .....	18
Project .....	39
Filter time .....	16
Folder .....	26, 39
Font size .....	43

## G

Grounding	
mains supply .....	10

**H**

Hardware key .....	18
Help	
Assistants.....	47
Context sensitive .....	47
Interactive .....	47
Online user manual .....	47
HyperMap .....	28, 102, 106
Hyperspectral data set .....	102, 106

**I**

Image	
Calibration .....	45
Capture.....	55
Detector selection.....	30, 34
Drift correction .....	59
Extension .....	57
Processing.....	61
Resolution .....	34
Image capture	
Area.....	30
Timing .....	34
Imaging .....	28
Import/Export menu .....	39
IO-Scan, Bruker-.....	18

**J**

Jobs .....	29
------------	----

**L**

Language settings.....	43
License.....	18
Light element window .....	20
Line average.....	34
Line scan.....	28
Acquisition.....	98
Qualitative .....	98
Quantitative.....	100
Load	
Image .....	39
Spectra, multiple .....	39
Spectrum.....	39

**M**

Magnification calibration.....	45
Mains voltage.....	10
Managing Project .....	114
Map	
Acquisition.....	102
Mapping .....	28, 102, 106
Quantitative.....	104
Maximum	
Energy (setting) .....	36
Maximum cooling (detector setting) .....	20, 36
Method editor .....	81
Microscope configuration.....	33

Multi channel analyzer .....	15
Multi-user systems .....	25

**O**

Objects .....	28
Online manual.....	47
Option panels.....	30

**P**

Phase analysis .....	109
Photon, counting.....	15
Position tagged spectrometry .....	17
Preamplifier.....	15
Product license .....	42
Project	
File .....	39
Managing .....	114
Project tool.....	38
Projects.....	39
Pulse throughput.....	16

**Q**

Quantification	
Bulk samples- .....	83, 86, 88
Electron transparent samples- .....	90, 92, 94
Standard-based- .....	86, 88
Standard-based Cliff-Lorimer- .....	92
Standard-based Zeta- .....	94
Standardless- .....	83
Standardless Cliff-Lorimer-.....	90
With library.....	88
With temporary reference.....	86
Quantum energy.....	15

**R**

Radiation safety .....	9
Real time.....	15
Registration.....	18
Remote access .....	25
Report.....	112
Report editor .....	38, 112
Resolution	
Image.....	34
Result	
Chart .....	30
Table .....	39
RS232	
Microscope interface .....	17

**S**

Safety	
Beryllium.....	10
Electrical- .....	10
Electromagnetic compatibility- .....	11
Radiation-.....	9
Sample	

Coating .....	32	System Factor Calibration .....	70
Coating correction .....	68	<b>T</b>	
Configuration .....	32	Thermostat mode (detector setting) .....	20, 36
Stress .....	34	Twain export .....	39
Tilt .....	33	<b>U</b>	
Save		USB Dongle .....	19
Image .....	39	User	
Spectrum .....	39	Deleting of- .....	27
Scan		Name .....	25, 26
Configuration .....	34	New- .....	26
Switch box .....	17	Password .....	25, 26
System .....	17	Profile .....	25
Scripting .....	29	Using Spectrum chart .....	65
SDD detector .....	15	<b>V</b>	
SE detector .....	30	Volumes	
Semiconductor detector, X-ray .....	15	Private .....	26
Server workstation .....	17	<b>W</b>	
Signal processing unit .....	15	Windows 7 .....	18
Software options .....	18	Working distance .....	51
Software version .....	42	Workspace	
Spectra .....	28	EBSD .....	28
Spectrometer		Feature .....	28
Calibration .....	53	Imaging .....	28, 55, 57, 59, 61
Configuration .....	36	Jobs .....	29
Electronic .....	15	Line scan .....	28, 98, 100
Energy dispersive- .....	15	Mapping .....	28, 102, 104, 106
Standby .....	36	Object .....	96
Spectroscopy signal processor .....	16	Objects .....	28
Spectrum .....	15	Scripting .....	29
Acquisition .....	63	Spectra .....	28, 63, 65
Chart .....	30	System .....	29, 41
Graphic output .....	39	Workstation	
Saving .....	39	Server .....	17
Spectrum acquisition		<b>X</b>	
Maximum speed .....	16	XFlash® detector .....	15, 19
Spectrum analysis		X-ray radiation .....	9
<i>Automatic</i> .....	96	<b>Z</b>	
Spot .....	30	Zeta factors	
Stage calibration .....	45	Experimental .....	78
Standard-based quantification .....	88	Reset .....	78
Standardless quantification .....	83		
Standards library .....	88		
Bulk samples .....	73		
Cliff-Lorimer factors .....	76		
Zeta factors .....	78		
Standby .....	19, 36		
Stress, sample- .....	34		
Supply voltage .....	10		
Switch Box, scan- .....	17		
System			
Assignments .....	41		
Control panels .....	29		
Report .....	41		
Settings .....	41		
Workspace .....	41		